

---

# HJ-150 Compact Scale

## Operating Instructions


---

### Cautions

- Avoid installing / leaving the scale in direct sunlight / high humidity.
- Place the scale on a flat and solid surface in a place free from wind.
- Avoid vibration, strong electric or magnetic fields and rapid changes in temperature or humidity.
- Avoid using the scale near other electronic equipment.
- Do not mix the battery types. Replace all of the batteries at same time.
- If the scale is not to be used for a long period of time, remove the batteries.
- Avoid over / shock load that could cause damage to the scale.
- Keep the protective cover attached when you carry the scale.
- Wait for more than 6 seconds to let the scale warm up after turning on the power.
- When the location to be used or the ambient environment has changed, calibrate the scale before use.

### Specifications

Capacity	150 x 0.1 g / 5.29 x 0.01 oz
Display	7 segment LCD with backlight (Character height 8 mm)
Power	2 x "AAA" size batteries
Pan size	60 (W)x 64 (D) mm / 2.36 (W) x 2.52 (D) in.
Accessories	Instruction manual, 100g weight, 2 monitor batteries

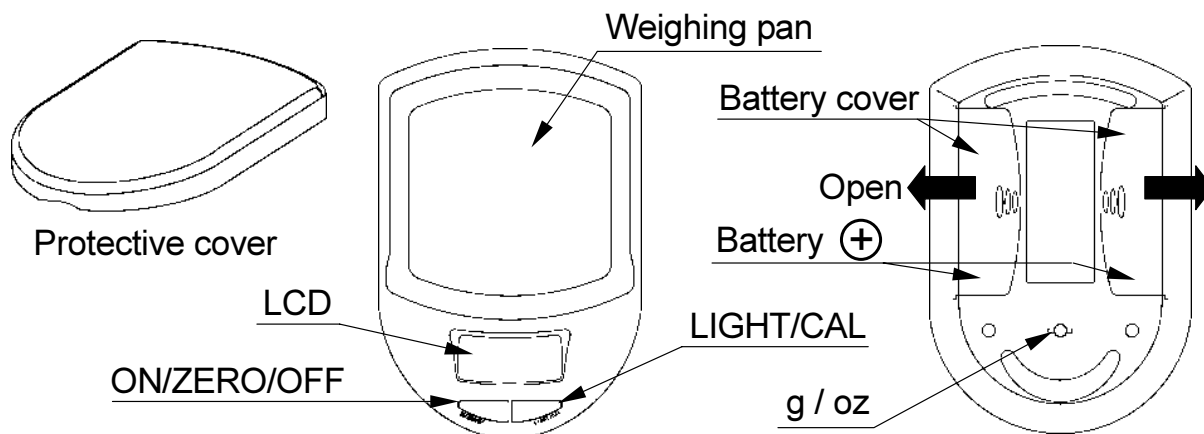

## Switches

**ON/ZERO/OFF** Switch:

Press to turn the scale on. In the operation mode, press briefly to set the display to zero. Press and hold approximately 3 seconds to turn the scale off.

**LIGHT/CAL** Switch:

In the operation mode, press briefly to turn the LCD backlight on. Press and hold approximately 3 seconds to enter the calibration mode.

**g / oz** Switch:

In the operation mode, press the switch located on the bottom of the scale to change the weighing unit between “g” and “oz”.

## Installing / Replacing batteries

Remove the battery covers and insert 2 AAA size batteries into each housing, taking care with the polarities. Attach the battery covers again.

- Keep the protective cover attached when installing and/or replacing batteries, to prevent excessive force from being added to the weighing pan.

## Weighing procedures

### 1. Turn on / off the scale

Press the **ON/ZERO/OFF** key to turn on the scale. The scale shows all segments for a few seconds and displays zero. The backlight will turn on for about one minute. Press and hold approximately 3 seconds to turn the scale off.

### 2. Select the weighing unit

Press the **g / oz** switch to select a weighing unit from “g” and “oz”.

### 3. Weighing

Verify the reading is zero. If not, press the **ON/ZERO/OFF** key to zero the display. If you use a container for weighing, place an empty container on the weighing pan and press the **ON/ZERO/OFF** key to zero the display.

Place the objects to be weighed on the pan or in the container. Wait for the display to become stable and read the weight.

- E** means overload. Remove the objects from the pan immediately.


## Automatic power-off function

The automatic power-off function turns power off after approximately 2 minutes of stable display during the object(s) applied. When there is no object applied, it will turn off after 30 seconds.

## Calibration

1. Make sure that there is nothing on the weighing pan, press and hold the **LIGHT/CAL** key approximately 3 seconds. The display shows **CAL**. Release the key.
  2. The scale memorizes the zero point and automatically displays **1000 g**.
  3. Place the calibration weight of 100g on the center of the pan within 10 seconds.
  4. After approximately 8 seconds, the scale will display **F** for success and **E** for failure. Then, the scale automatically turns off.
  5. Turn the scale on again to start the operation.
- If calibration fails, check if anything is touching the weighing pan, there is vibration or wind, and the calibration weight is correct.

## Dimension

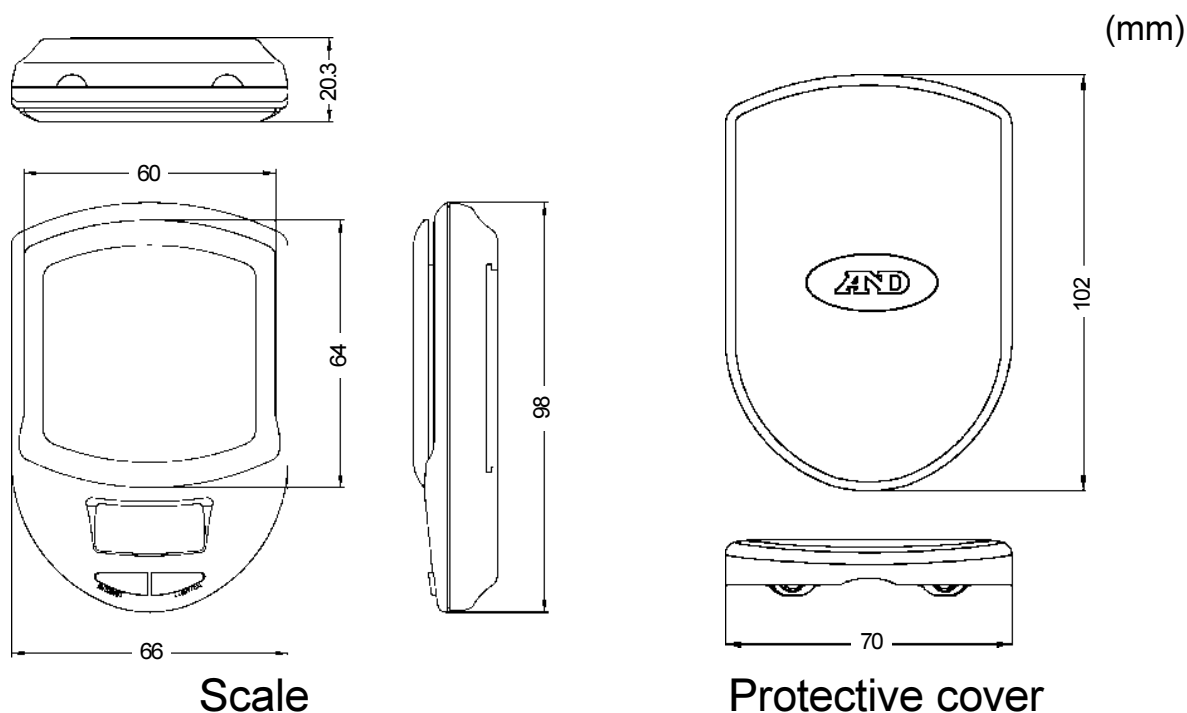



**A&D Company, Limited**

3-23-14 Higashi-Ikebukuro, Toshima-ku, Tokyo 170-0013 JAPAN  
Telephone: [81] (3) 5391-6132 Fax: [81] (3) 5391-6148

**A&D ENGINEERING, INC.**

1555 McCandless Drive, Milpitas, CA. 95035 U.S.A.  
Telephone: [1] (408) 263-5333 Fax: [1] (408) 263-0119

**A&D INSTRUMENTS LTD.**

Unit 24/26 Blacklands Way Abingdon Business Park, Abingdon, Oxon OX14 1DY U.K.  
Telephone: [44] (1235) 550420 Fax: [44] (1235) 550485

**<German Sales Office>**

Große Straße 13b 22926 Ahrensburg GERMANY  
Telephone: [49] (0) 4102 459230 Fax: [49] (0) 4102 459231

**A&D MERCURY PTY. LTD.**

32 Dew Street, Thebarton, South Australia 5031 AUSTRALIA  
Telephone: [61] (8) 8301-8100 Fax: [61] (8) 8352-7409

**A&D KOREA Limited**

Manhattan Bldg. 8F, 36-2 Yoido-dong, Youngdeungpo-gu, Seoul, KOREA  
Telephone: [82] (2) 780-4101 Fax: [82] (2) 782-4280