# BM series

# INSTRUCTION MANUAL

# **Analytical Balance**

BM-5

BM-5D

BM-20

BM-22

BM-252

BM-200

BM-300

BM-500



© 2019 A&D Company, Limited. All rights reserved.

No part of this publication may be reproduced, transmitted, transcribed, or translated into any language in any form by any means without the written permission of A&D Company, Limited.

The contents of this manual and the specifications of the instrument covered by this manual are subject to change for improvement without notice.

Windows, Word and Excel are registered trademarks of Microsoft Corporation.

# Contents

<b>Basic Operation</b>
------------------------

1. 1.1. 1.2.	IntroductionFeaturesCompliance	4
2.	Unpacking the Balance	7
3. 3.1. 3.2. 3.3. 3.4. 3.5.	Installing the Balance and Precautions Installing the Balance Precautions Before Use Precautions During Use Precautions After Use Precautions for Power Supply	10 13 15 16
4.	Display Symbols and Key Operation	17
5. 5.1. 5.2. 5.3.	Weighing Units Units Storing Active Units Selecting Units and Weighing Modes	18 22
6. 6.1. 6.2. 6.3. 6.4. 6.5. 6.6. 6.7.	Weighing Basic Operation (Gram Mode) Smart Range Counting Mode (PC) Percent Mode (%) Ionizer (Neutralization Device) Monitoring the Clock and Environmental Sensors Repeatability Test	
Adaptir	ng to the Environment	
7. 7.1. 7.2.	Response Adjustment	35
8. 8.1. 8.2. 8.3. 8.4. 8.5. 8.6. 8.7.	Calibration Calibration Group Automatic Self Calibration Calibration Using the Internal Mass Calibration Test Using the Internal Mass Calibration Using an External Weight Calibration Test Using an External Weight Correcting the Internal Mass Value	
Selectin	ng Functions	
9. 9.1. 9.2.	Function Switch and Initialization Permit or Inhibit Initializing the Balance	44
10. 10.1. 10.2. 10.3. 10.4.	Function Table	46 48 50

10.5.	Description of the Item "Data Format"	53
10.6.	Data Format Examples	
10.7.	Clock and Calendar Function	58
11.	ID Number and GLP Report	59
11.1.	Setting the ID Number	59
11.2.	GLP Report	60
12.	Data Memory	66
12.1.	Notes on Using Data Memory	66
12.2.	Data Memory for Weighing Data	
12.3.	Data Memory for Calibration and Calibration Test Data	70
13.	Underhook	72
14.	Density Measurement (Specific gravity)	73
15.	Password Function	76
15.1.	Balance Software Version 1.50 to 1.64	76
15.2.	Balance Software Version 1.660 or later	
15.3.	Using the Password Function	
15.4.	Entering the Password Before Weighing	
15.5.	Logging Out	
15.6. 15.7.	Storing or Changing the PasswordChanging the Password	
15.7.	Deleting the Password (USER [] I~I[])	
15.0.	Forgotten Password	
	ce and Communication	
16.	Standard Input and Output Interface	
16.1.		
16.2.	Connecting to Peripheral Equipment	
16.3.	Commands	
17.	Key Lock Function	
17.1.	J	
17.2.	Locking the Specified Keys	96
18.	Checking the Balance Software Version	96
Mainte	nance	
19.	Maintenance	97
19.1.		
19.2.	Error Codes	97
19.3.	Other Display	
19.4.	5	
19.5.	Asking for Repair	
20.	Specifications	
20.1.		
20.2.	Options and Peripheral Equipment	104
21.	Terms/Index	108
21.1.	Terms	
21.2.	Index	109

2

# 1. Introduction

This manual describes how the **BM** series balances work and how to get the most out of them in terms of performance. Read this manual thoroughly before using the balance and keep it at hand for future reference.

#### About this manual

This manual consists of the following five parts:

operation.

Adapting to the environment......... Describes response (and stability) adjustment to adapt to the environment where there is vibration or drafts, the way to maintain weighing precision in a variation of ambient temperature, calibration and calibration test.

Basic operation ...... Describes precautions, the balance's construction and basic

Selecting functions ...... Describes functions of the balance.

Interface and communication...... Describes the serial interface used for communicating with a

computer that requests weighing data and controls the

balance, and for use with a printer

Maintenance ...... Describes maintenance, error codes, troubleshooting,

specifications and options.

#### **Balance software version**

Some balance operations vary depending on the balance software version. For details on checking the software version, refer to "18. Checking the Balance Software Version".

3

# 1.1. Features

- A built-in ionizer can eliminate static electricity from the weighing sample before weighing, reducing weighing errors.
- Each electrode unit of the ionizer is designed to be removed, cleaned and replaced.
- Automatic self calibration, using the internal mass, adapting to changes in temperature.
- Response adjustment adapting to drafts and/or vibration in the environment automatically.
- Memory function to store weighing data and calibration data.
   When only weighing data is stored, a maximum of 200 data can be stored.
   Interval mode to store the weighing data periodically.
- Good laboratory practice (GLP) / Good manufacturing practice (GMP) data can be output using the RS-232C serial interface.
- A built-in clock and calendar that can add the time and date to the output data. (Changing the clock setting can be limited to the administrator. Refer to "15. Password Function".)
- The password function can be used to restrict balance users or changes to the function settings.
- The key lock function disables the balance key operation and enables the balance to be operated only by commands from external devices.
- Underhook, for measuring density and weighing magnetic materials.
- Multiple weighing units with most of the common units used around the world.
   Gram, Milligram, Counting mode, Percent mode, Ounce (Avoir), Troy Ounce, Metric carat, Momme,
   Pennyweight, Grain (UK), Tael (preset at the factory), Tola (India), Messghal and Density mode
- The BM-5D and BM-22 are equipped with a smart range function to weigh in the precision range (readability: 1 μg) after zeroing a tare weight, if within the weighing capacity.
- □ The micro tube holder is included in the accessories of the BM-5, BM-5D, BM-20 and BM-22.
- Density mode for calculating the density of a solid.
- The BM series balances are equipped with an RS-232C serial interface and a USB interface to communicate with a computer. While outputting data to a printer using the RS-232C interface, data can be transmitted to a computer at the same time using the USB interface. Using the RS-232C interface and the Windows communication tools software (WinCT) enables bidirectional communication between a Windows computer and the balance. The current version of the WinCT can be downloaded from the A&D website.
- Connecting the USB interface of the balance by a cable to a Windows computer allows transmission of the weighing data to Excel or Word.
- When multiple balances have the BM-08 Ethernet interface installed in place of the USB interface and connected to a LAN, data can be acquired from each of them using the WinCT-Plus software.
- □ When the accessory data logger (AD-1688) is connected, weighing data can be stored without using a computer.

# 1.2. Compliance

# 1.2.1. Compliance with FCC Rules

Please note that this equipment generates, uses and can radiate radio frequency energy. This equipment has been tested and has been found to comply with the limits of a Class A computing device pursuant to Subpart J of Part 15 of FCC rules. These rules are designed to provide reasonable protection against interference when equipment is operated in a commercial environment. If this unit is operated in a residential area, it may cause some interference and under these circumstances the user would be required to take, at his own expense, whatever measures are necessary to eliminate the interference.

(FCC = Federal Communications Commission in the U.S.A.)

# 1.2.2. Compliance with Directives of CE mark

This device features radio interference suppression, safety regulation and restriction of Hazardous Substances in compliance with the following Council Directives

Council directive 2014/30/EU EN61326 EMC directive

Council directive 2014/35/EU EN60950 Safety of Information Technology Equipment

Council directive 2011/65/EU EN50581 Restriction of the use of certain Hazardous Substances

The CE mark is an official mandatory European marking.

Please note that any electronic product must comply with local laws and regulations when sold or used anywhere outside Europe.

5

# **EU Declaration of Conformity**

CE

A&D Instruments Ltd. hereby declare that the following Weighing product conforms to the requirements of the council directives on ...

Electromagnetic Compatibility (EMC) 2014/30/EU,

Low Voltage Equipment (LVD) 2014/35/EU and

Restriction of the use of certain Hazardous Substances (RoHS) 2011/65/EU

provided that they bear the CE mark of conformity.

Model/Series....BM Series

Standards applicable:

#### EN 61326-1:2013

Electrical equipment for measurement, control and laboratory use -EMC requirements Part 1: General requirements

#### EN 60950-1:2006+A11:2009+A1:2010+A12:2011+A2:2013

Safety of Information Technology Equipment

#### EN 50581:2012

Technical documentation for the assessment of electrical and electronic products with respect to the restriction of hazardous substances

6

This Declaration of Conformity is issued under the sole responsibility of the manufacturer A&D instruments, 24-26 Blacklands Way, Abingdon, Oxfordshire OX14 1DY, England

CE Mark first applied 28 January 2011 Signed for A&D Instruments on 02 July 2016

J. Chuman Managing Director


www.aandd-eu.net



# 2. Unpacking the Balance

Unpack the balance carefully. Keep the packing material to be used for transporting the balance in the future. See the illustrations to confirm that everything is included.

#### Accessories for each model

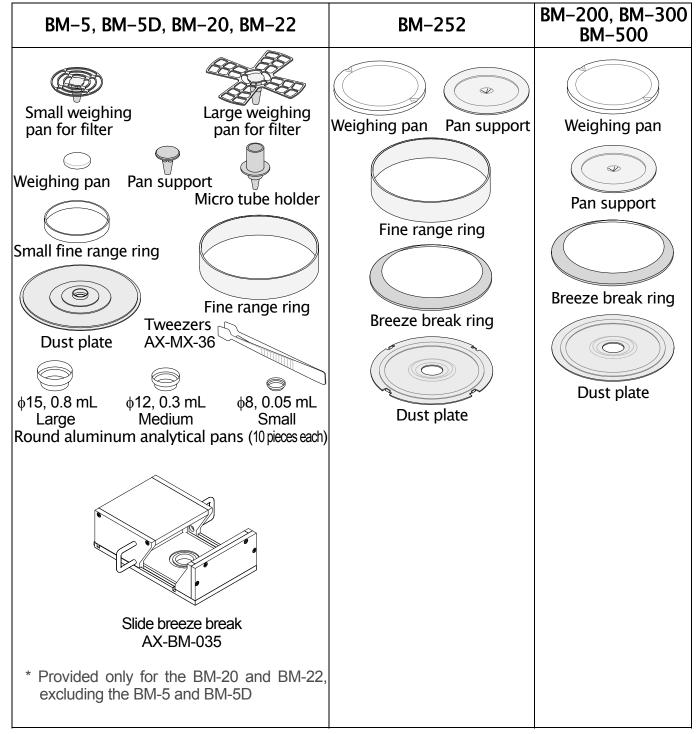

#### Common accessories

Positions to place the AC adapter labels

Attach the AC adapter labels to the AC adapter as shown in the illustration.


識別シール A Cアダブタ、コード等に貼って下さい Identification Sticker Attach Label to the AC Adapter and its Cable			
BM BM			
ВМ	BM		

AC adapter

AC adapter labels

#### **Notes**


- Please confirm that the AC adapter type is correct for your local voltage and receptacle type.
- Please use the dedicated AC adapter specified for the balance.
- Do not use the AC adapter provided with the balance for other models or equipment with which the AC adapter may not be compatible.
- If you use the wrong AC adapter, the balance and other equipment may not operate properly.

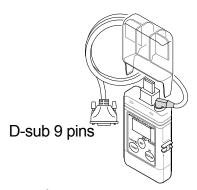

Data logger AD-1688

The weighing data can be stored when connected to the RS-232C interface of the balance.

AD.1689 SUS.304

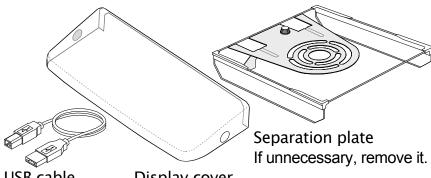


USB cable (1.5 m or 1.8 m)

Display cover (PET resin)

Tweezers AD-1689

Tweezers for handling a weight.


Use it to avoid temperature changes due to having your hand in the weighing chamber, thus to perform accurate calibration.

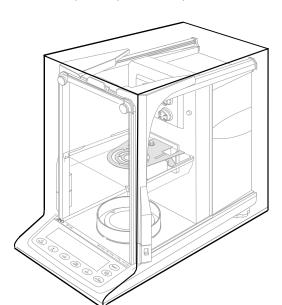

**Body cover** 

Cover during storage to prevent dust and dirt contamination.


The cover can be used for the balance with the AC adapter and other cables connected.

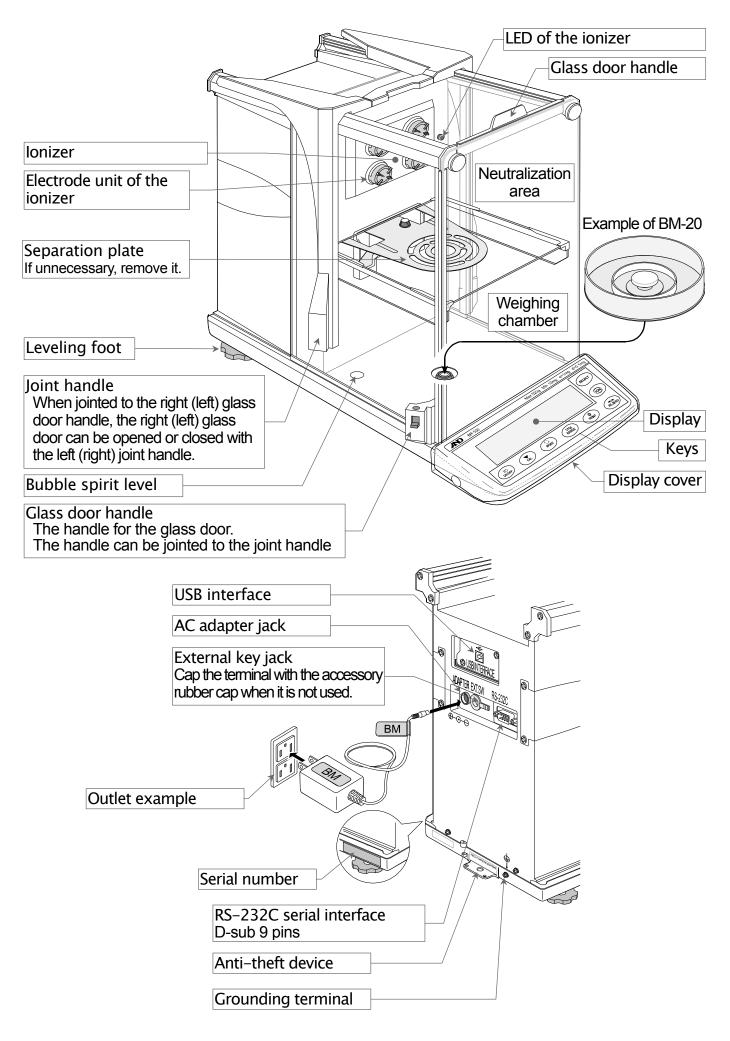

# 3. Installing the Balance and Precautions

# 3.1. Installing the Balance

- 1 Place the balance on a solid weighing table. Refer to "3.2. Precautions Before Use" for details on a place to install the balance.
- 2 Assemble the weighing pan and other parts in accordance with the model and use as shown below.

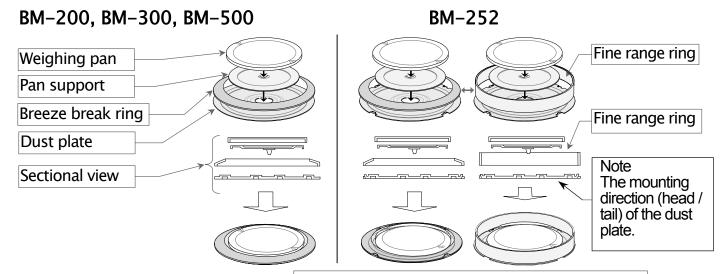


Use the fine range ring in place of the breeze break ring to avoid errors caused by drafts when weighing with a readability of 0.01 mg.


#### BM-5, BM-5D, BM-20, BM-22

Note When using the micro tube holder for the BM-5 and BM-5D:


The micro tube holder is approximately 2 g heavier than the standard pan (weighing pan and pan support). When a micro tube holder is used for the BM-5 and BM-5D, the available weighing range will be up to approximately 3 g including the tare such as a micro tube and test tube.

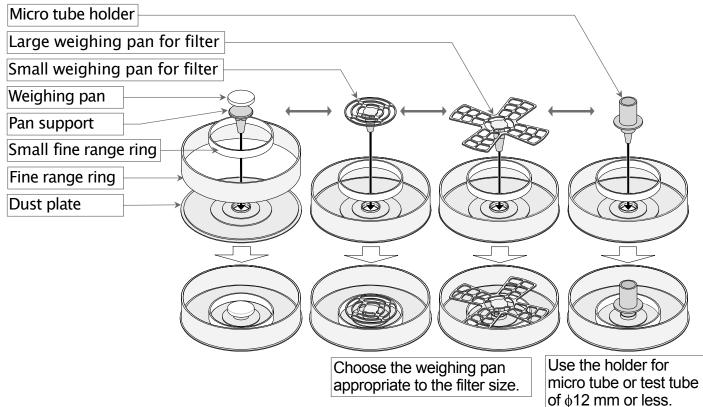

# Assembling the parts in the weighing chamber

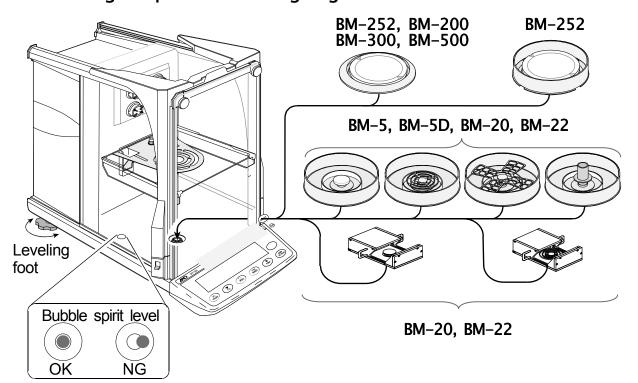


- 3 If the separation plate is not necessary, remove the screw and take the plate off. Refer to "3.1.2. Removing and attaching the separation plate".
- 4 Adjust the level of the balance using the leveling feet. (Refer to "Adjusting the Level" on page 14.)
- 5 Confirm that the AC adapter type is correct for your local voltage and power receptacle type.
- 6 Connect the specified AC adapter to the balance. Warm up the balance for at least one hour with nothing on the weighing pan.

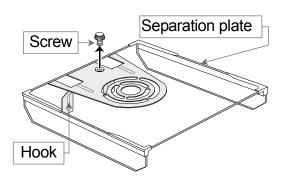

# 3.1.1. Installing the slide breeze break BM-20 / BM-22

The slide breeze break improves weighing precision by reducing the negative effects from the surrounding environment. It is installed inside the weighing chamber of the BM-20 / BM-22.

- 1 Determine positioning of the slide breeze break to align the slide breeze break (base) hole with the weighing pan/pan support mounted section while pressing the sponge, located at the rear of the slide breeze break (base), at an angle to the BM-20 or BM-22 wall surface.
- 2 Attach the weighing pan/pan support and he slide breeze break (door).

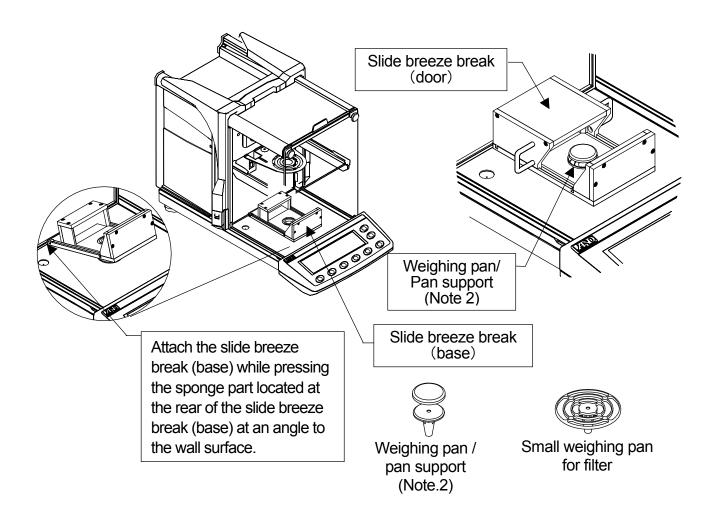

- Note 1 There may be variations in weighing results after attaching the slide breeze break as it has not adjusted to the surrounding temperature. Wait a while before making any measurements.
- Note 2 The weighing pan / pan support or small weighing pan for filter provided with the BM-20 / BM-22 can be used.

# 3.1.2. Removing and attaching the separation plate

Note Take care not to damage the separation plate when handling it.


# Removing the separation plate

- 1 Support the separation plate and remove the screw from the plate.
- 2 Lift the plate and unhook the hooks.
- 3 Rotate the plate on the axis of the front edge. Rotate the plate on the axis of the back edge.
- 4 Remove the plate from the chamber.

#### Attaching the separation plate

Use the arrows in reverse to attach the plate.

- 5 Insert the plate into the chamber.
- 6 Rotate the plate on the axis of the back edge. Rotate the plate on the axis of the front edge. Level the plate.
- 7 Insert the hooks and hook the plate.
- 8 Support the plate and secure the screw.

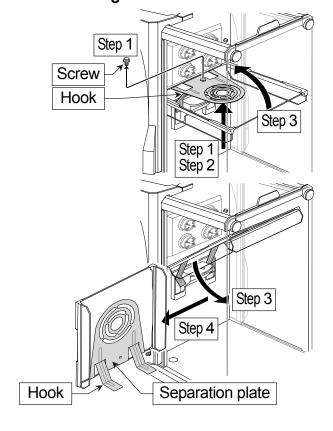


# 3.2. Precautions Before Use

To ensure that you get the most from your balance, please try to follow these conditions as closely as possible, especially for the BM-5, BM-5D, BM-20 and BM-22 that are sensitive.

- The best operating temperature is about 20°C / 68°F at about 50% Relative Humidity.
- The weighing room should be free of dust.
- Clean the weighing chamber and its surroundings to prevent foreign matter from invading, especially when weighing a very small amount of sample.
- □ The weighing table should be solid and free from vibration, drafts (such as frequently opening doors or windows) and as level as possible. We recommend that you use the anti-vibration table (AD–1671) and the remote controller (AD–8922A) for the BM–5, BM–5D, BM–20 and BM–22.
- Do not install the balance where it will be subject to vibration. Corners of rooms on the first floor are best.
- Do not install the balance near a heater, air conditioner, or in a breeze. The influence of breezes can be reduced by using the tabletop breeze break (L) (AD-1672/AD-1672A) or the tabletop breeze break (M) (AD-1676).
- Do not install the balance in direct sunlight. Avoid excessive temperature changes.
- Do not use the balance near other equipment which produces magnetic fields.
- Adjust the level of the balance using the leveling feet.
- Please warm up the balance for at least one hour. Plug in the AC adapter as usual.
- Calibrate the balance if it is installed for the first time or relocated, or when a weighing pan is replaced.
- Ensure a stable power source when using the AC adapter.

⚠ Do not place or use the balance where there is flammable or corrosive gas present.


#### Adjusting the level

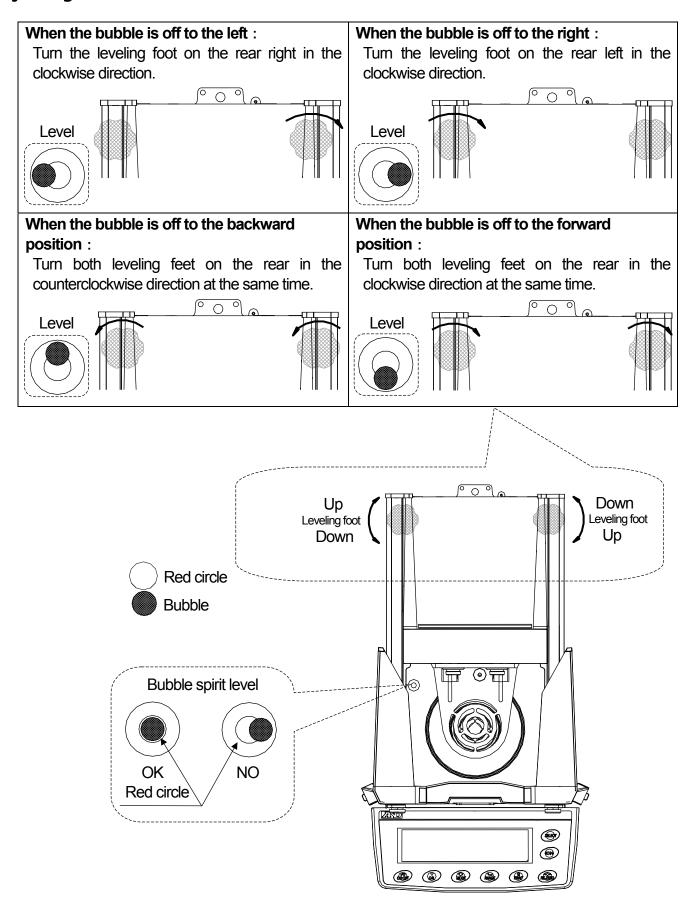

# 3.3. Precautions During Use

Note the following items to get accurate weighing data.

- Discharge static electricity from the weighing sample. When a weighing sample (plastics, insulator, etc.) could have a static charge, the weight value is influenced. Ground the balance, and
  - Eliminate the static electricity using the built-in ionizer.
  - Try to keep the ambient humidity at or above 45% RH in the room.
  - Use a metal shield case.
  - Wipe a charged material (plastic sample, etc.) with a damp cloth.
- This balance uses a strong magnet as part of the balance assembly, so please use caution when weighing magnetic materials. If there is a problem, use the underhook (refer to "13. Underhook") to suspend the material away from the influence of the magnet.
- Eliminate temperature differences between the weighing sample and the environment. When a sample is warmer (cooler) than the ambient temperature, the sample will appear lighter (heavier) than the true mass. This error is due to the rising (falling) draft around the sample. If you touch the sample, the same type of error will occur. Do not touch the sample directly with your hand. Use tweezers or other tools.
- Do not drop things upon the weighing pan, or place a weight beyond the range of the balance on the weighing pan.
- Make each weighing gently and quickly to avoid errors due to changes in the environmental conditions.
- We recommend that the fine range ring and separation plate are used to avoid weighing errors caused by drafts when weighing using the BM-252 (readability: 0.01 mg).
- □ We recommend that the **small fine range ring**, **fine range ring** and **separation plate** are used to avoid weighing errors caused by drafts when weighing using the **BM**−20 or **BM**−22 (readability: 0.01 mg and 0.001 mg). To reduce the influence of breezes, minimize operation of doors of the weighing chamber. In addition, in order to weigh more accurately while reducing the effects of the surrounding environment, attach the slide breeze break to the weighing chamber.
- The weighing value of the BM-5, BM-5D, BM-20 and BM-22 may change after the stabilization indicator is displayed. We recommend that you set a certain amount of wait time (approx. 10 seconds) before reading the value after the stabilization indicator is displayed.
- If contamination may be a problem, perform sampling operations outside the weighing chamber in order to prevent the sample from scattering in the weighing chamber.
- Take into consideration the affect of air buoyancy on a sample when more accuracy is required.
- Do not use a sharp instrument (such as a pencil or ball point pen) to press the keys, use your finger only.

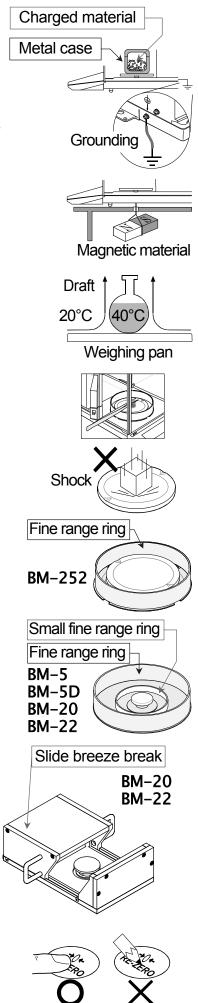

- Press the RE-ZERO key before each weighing to prevent possible errors.
- Keep the balance interior free of foreign matter (dust, liquid or metal fragments).


# 3.4. Precautions After Use

- Avoid mechanical shock to your balance.
- Do not disassemble the balance. Contact your local A&D dealer if your balance needs service or repair.
- Do not use solvents to clean the balance. For best cleaning, wipe with a dry lint free cloth or a lint free cloth that is moistened with warm water and a mild detergent.
- Keep the balance interior free of foreign matter (dust, liquid or metal fragments).

# 3.5. Precautions for Power Supply

- Do not remove the AC adapter while the internal mass is in motion, for example, right after the AC adapter is connected, or during calibration using the internal mass.
   If the AC adapter is removed under the conditions described above, the internal mass will be left unsecured, that may cause mechanical damage when the balance is moved.
   Before removing the AC adapter, press the ON:OFF key and confirm that zero is displayed in the weighing mode.
- After connecting the AC adapter to the balance, warm up the balance for at least one hour before weighing.
- □ When the AC adapter is connected, the balance is in the standby mode if the standby indicator is on. This is a normal state and does not harm the balance. For accurate weighing, we recommend that the AC adapter be kept connected, especially for the BM-5, BM-5D, BM-20 and BM-22.

# 4. Display Symbols and Key Operation

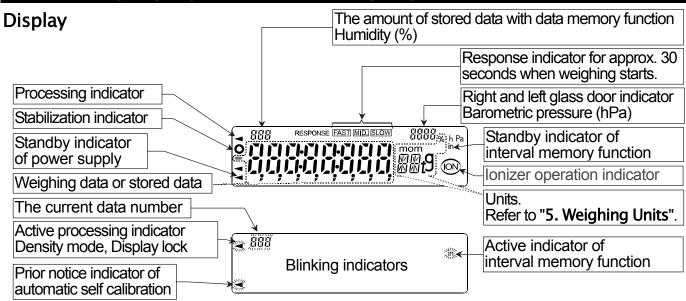

The interval memory function is used to store the weighing data periodically. Refer to "12. Data Memory".

# **Key operation**



Key When pressed and released When pressed and held (approx. 2 seconds) Tums the display on and off. The standby indicator is displayed when the display is tumed off. When the display is turned on, the weighing mode is enabled. If the password function is used, PRSS III is displayed to prompt you to enter a password. Refer to "15.4. Entering" I/O ON:OFF the Password Before Weighing". This key is available anytime. Pressing the key during operation will interrupt the operation and turn the display off. Cancels the operation when performing function settings. CAL Displays other items of the calibration menu. Enters the calibration mode using the internal mass. Switches the preset weighing units stored in the function table. Performs automatic response adjustment. Refer to "5. Weighing Units". Displays the function table menu. In the weighing mode, turns the minimum Refer to "10. Function Table". weighing value on and off. Performs a repeatability test when pressed 1/10d RANGE □ In the counting or percent mode, enters and held for another 2 seconds after the the unit mass storing mode or reference function table menu is displayed. mass storing mode. Refer to "6.7. Repeatability Test". No function at the factory setting. The key to output the weighing data to a By changing the function table: printer or personal computer (or store it in "Title block" and "End block" for GLP / memory) depending on the function table GMP report are output. (Factory setting = output) settings. The data memory menu is displayed. →0← RE-ZERO The key to set the display to zero. (ION) The key to turn the ionizer on and off. The kev to monitor the time and date, and the environmental sensors (temperature, humidity, barometric

SELECT

pressure), for several seconds. Refer to "6.6. Monitoring the Clock and Environmental Sensors".

# 5. Weighing Units

# **5.1.** Units

The units and weighing modes can be selected and stored in the function table as described in "5.2. Storing Active Units".

The sequence of displaying them can be arranged to fit the frequency of use.

They are stored and are maintained in non-volatile memory, even if the AC adapter is removed.

- If the law in your area permits, you may use all of the units. You can disable the units that you don't regularly use. And you are able to turn them back on.
- If a weighing mode (or unit of mass) has been turned off, that mode or unit will be missing in the sequence. Tael has four varieties, one of which can be selected and installed at the factory.
- Press the MODE key to select a unit or mode for weighing.
- For details about the units and modes, see the table below:

Name (unit, mode)	Abbreviation	Display unit	Conversion factor
Gram	g	9	1 g
Milligram	mg	mg	0.001 g
Counting mode	PC	PE	-
Percent mode	%	%	-
Ounce (Avoir)	OZ	OΖ	28.349523125 g
Troy Ounce	ozt	ΩZt	31.1034768 g
Metric Carat	ct	ct	0.2 g
Momme	mom	mom	3.75 g
Pennyweight	dwt	dnt	1.55517384 g
Grain (UK)	GN	<u>G</u> N	0.06479891 g
Tael (HK general, Singapore)			37.7994 g
Tael (HK jewelry)	tl	11	37.429 g
Tael (Taiwan)	U	EL	37.5 g
Tael (China)			31.25 g
Tola (India)	t	t	11.6638038 g
Messghal	m	m	4.6875 g
Density mode	DS	Refer to "14. De	ensity Measurement"

#### Density mode

- To use the density mode, it must be stored in the function table as described on page 22. For details about this mode, refer to "14. Density Measurement".
- To select the density mode, press the MODE key until the processing indictor **\*\*** blinks with the unit "g" displayed.

# Capacity and readability for the BM-252, BM-200, BM-300, BM-500

□ The tables below indicate the weighing capacity and the readability for each balance.

Unit	BM-200	BM-300	BM-500	
Orin	Capacity			Readability
Gram	220	320	520	0.0001
Milligram	220000	320000	520000	0.1
Ounce (Avoir)	7.76	11.2	18.3	0.00001
Troy Ounce	7.07	10.2	16.7	0.00001
Metric Carat	1100	1600	2600	0.001
Momme	58.6	85.3	138	0.0001
Pennyweight	141	205	334	0.0001
Grain (UK)	3395	4938	8024	0.002
Tael (HK general, Singapore)	5.82	8.46	13.7	0.00001
Tael (HK jewelry)	5.87	8.54	13.8	0.00001
Tael (Taiwan)	5.86	8.53	13.8	0.00001
Tael (China)	7.04	10.2	16.6	0.00001
Tola (India)	18.8	27.4	44.5	0.00001
Messghal	46.9	68.2	110	0.0001

Unit	BM-252		
Orint	Capacity	Readability	
Gram	250	0.00001	
Milligram	250000	0.01	
Ounce (Avoir)	8.81	0.000001	
Troy Ounce	8.03	0.000001	
Metric Carat	1250	0.0001	
Momme	66.6	0.00001	
Pennyweight	160	0.00001	
Grain (UK)	3858	0.0002	
Tael (HK general, Singapore)	6.61	0.00001	
Tael (HK jewelry)	6.67	0.00001	
Tael (Taiwan)	6.66	0.000001	
Tael (China)	8.00	0.00001	
Tola (India)	21.4	0.00001	
Messghal	53.3	0.00001	

# Capacity and readability for the BM-5 and BM-5D

□ The **BM-5D** is equipped with the precision range and standard range of the smart range function.

	BM-5		
Unit	Precision range		
	Capacity	Readability	
Gram	5.20	0.000001	
Milligram	5200	0.001	
Ounce (Avoir)	0.183	0.0000001	
Troy Ounce	0.167	0.000001	
Metric Carat	26.0	0.00001	
Momme	1.38	0.000001	
Pennyweight	3.33	0.000001	
Grain (UK)	80.1	0.00002	
Tael (HK general, Singapore)	0.137	0.0000001	
Tael (HK jewelry)	0.138	0.0000001	
Tael (Taiwan)	0.138	0.0000001	
Tael (China)	0.166	0.0000001	
Tola (India)	0.444	0.000001	
Messghal	1.10	0.00001	

	BM-5D				
Unit	Precision range		Standard range		
	Capacity	Readability	Capacity	Readability	
Gram	2.10	0.000001	5.20	0.00001	
Milligram	2100	0.001	5200	0.01	
Ounce (Avoir)	0.0737	0.000001	0.183	0.000001	
Troy Ounce	0.0671	0.000001	0.167	0.000001	
Metric Carat	10.5	0.00001	26.0	0.0001	
Momme	0.560	0.000001	1.38	0.00001	
Pennyweight	1.34	0.000001	3.33	0.00001	
Grain (UK)	32.4	0.00002	80.1	0.0001	
Tael (HK general, Singapore)	0.0551	0.000001	0.137	0.000001	
Tael (HK jewelry)	0.0560	0.000001	0.138	0.000001	
Tael (Taiwan)	0.0560	0.000001	0.138	0.000001	
Tael (China)	0.0671	0.000001	0.166	0.000001	
Tola (India)	0.179	0.000001	0.444	0.000001	
Messghal	0.444	0.000001	1.10	0.00001	

# Capacity and readability for the BM-20 and BM-22

□ The **BM-22** is equipped with the precision range and standard range of the smart range function.

	BM-20		
Unit	Precision range		
	Capacity	Readability	
Gram	22.0	0.000001	
Milligram	22000	0.001	
Ounce (Avoir)	0.776	0.000001	
Troy Ounce	0.707	0.000001	
Metric Carat	110	0.00001	
Momme	5.86	0.000001	
Pennyweight	14.1	0.000001	
Grain (UK)	339	0.00002	
Tael (HK general, Singapore)	0.582	0.0000001	
Tael (HK jewelry)	0.587	0.000001	
Tael (Taiwan)	0.586	0.000001	
Tael (China)	0.704	0.000001	
Tola (India)	1.88	0.000001	
Messghal	4.69	0.000001	

	BM-22				
Unit	Precision range		Standard range		
	Capacity	Readability	Capacity	Readability	
Gram	5.10	0.000001	22.0	0.00001	
Milligram	5100	0.001	22000	0.01	
Ounce (Avoir)	0.179	0.000001	0.776	0.000001	
Troy Ounce	0.163	0.000001	0.707	0.000001	
Metric Carat	25.5	0.00001	110	0.0001	
Momme	1.36	0.000001	5.86	0.00001	
Pennyweight	3.27	0.000001	14.1	0.00001	
Grain (UK)	78.7	0.00002	339	0.0001	
Tael (HK general, Singapore)	0.134	0.000001	0.582	0.000001	
Tael (HK jewelry)	0.136	0.000001	0.587	0.000001	
Tael (Taiwan)	0.136	0.000001	0.586	0.000001	
Tael (China)	0.163	0.000001	0.704	0.000001	
Tola (India)	0.437	0.000001	1.88	0.000001	
Messghal	1.08	0.000001	4.69	0.00001	

# 5.2. Storing Active Units

- The units and modes can be selected and stored in the function table.
   The sequence of displaying them can be arranged to fit the frequency of use.
   The units stored are maintained in non-volatile memory, even if the AC adapter is removed.
- 1 In the weighing mode, press and hold the RANGE key (for approx. 2 seconds) until bR5Fnc is displayed, and then release the key.
- 2 Press the RANGE key several times to display Unit.
- 3 Press the PRINT key to enter the unit selection mode.
- Specify a unit or mode in the order to be displayed using the following keys.

RANGE key ......To display the units sequentially.


RE-ZERO key ······To specify a unit or mode.

The indicator **O** appears when the displayed unit or mode is specified.

Examples

Unit		Display
Gram	g	° lln it 9
Milligram	mg	° Unit mg
Counting mode	PC	° Unit PC
Percent mode	%	° lin it %
Density mode	DS	° Unit 115

- 5 Press the PRINT key to store the units or modes. The balance displays End and then displays the next menu item of the function table.
- 6 Press the CAL key to exit the function table. Then the balance returns to the weighing mode with the selected unit.

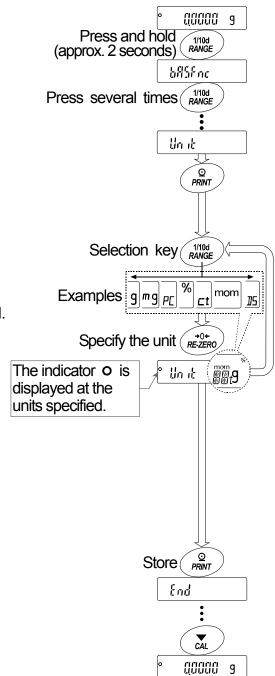


# 5.3. Selecting Units and Weighing Modes

- Preset the sequence of displaying units and weighing modes to fit the frequency of use.
   Refer to "5.2. Storing Active Units".
- Press the MODE key in the weighing mode to display the units and weighing modes in order.
- □ For the BM-5, BM-5D, BM-20 and BM-22, the units preset at the factory, <sup>m</sup>9 (Milligram) and 9 (Gram), are displayed in this order.
- For the BM-252, BM-200, BM-300 and BM-500, the units preset at the factory, <sup>9</sup> (Gram) and <sup>m</sup>9 (Milligram), are displayed in this order.
- Density mode
  - To use the density mode, it must be stored in the function table as described on page 22.
     Refer to "14. Density Measurement" for details about the mode.
  - To select the density mode, press the MODE key until the processing indictor 

    blinks with the unit "g" displayed.

The unit sequence of the factory settings. **Turning the display on.** 

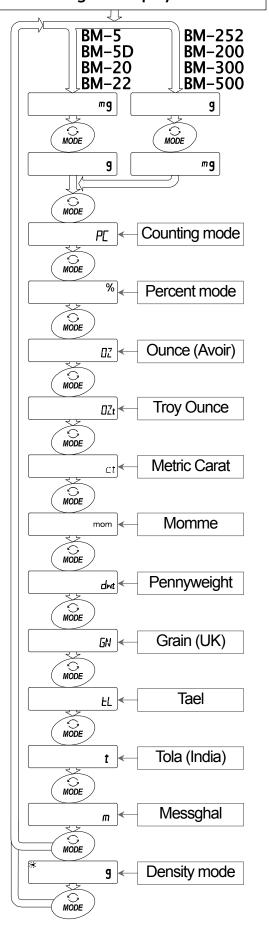


# 5.3.1. Unit Setting Example

The example below sets the units in the order with g (gram) as the first unit followed by PC (counting mode).

- 1 In the weighing mode, press and hold the RANGE key (for approx. 2 seconds) until bR5Fnc is displayed, and then release the key.
- 2 Press the RANGE key several times to display Unit.
- 3 Press the PRINT key to enter the unit selection mode.
- 4 Press the RE-ZERO key to specify the unit of g.

  The stabilization indicator 

  appears when the unit is specified.
- 5 Press the RANGE key several times to display Unit PE.
- 6 Press the RE-ZERO key to specify the unit of PC.
  The stabilization indicator **o** appears when the unit is specified.
- 7 Press the PRINT key to store the units.
  The balance displays End and then displays the next menu item of the function table.
- 8 Press the CAL key to exit the function table. Then the balance returns to the weighing mode with g, the unit selected first.
- 9 Press the MODE key to switch between g and PC.

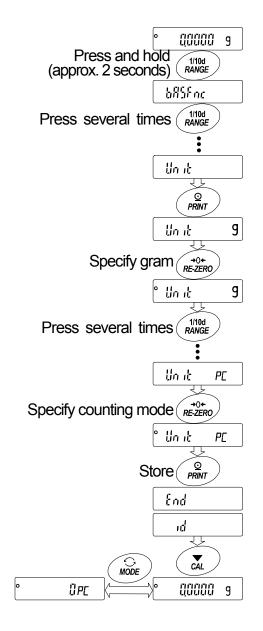

# 6. Weighing

# Precautions for the weighing operation

- □ Press the RE-ZERO key each time, before placing a sample on the weighing pan, to prevent possible errors.
- Place a sample in the center of the weighing pan gently.
- Temperature changes during measurement may cause weighing errors.
- Shorten the operation time as much as possible. (Opening and closing door, placing and removing sample)
   Material with an electrostatic charge or that is magnetic may cause weighing errors.
- Use a pair of tweezers to avoid a temperature change due to having your hand in the weighing chamber.
- Do not drop things on the pan, or place a weight on the pan that is beyond the weighing range.
- Keep the area clean and dry.
- Do not press keys with a sharp instrument (such as a pencil or ball point pen).
- □ For precision weighing, keep the AC adapter connected to the balance.
- Calibrate periodically to maintain weighing accuracy. Refer to "8. Calibration".
- □ For the BM-5, BM-5D, BM-20 and BM-22, perform pre-weighing of the container (tare) so that precision weighing can be performed.
- Consider section "3. Installing the Balance and Precautions" for the weighing operation.

# **6.1. Basic Operation** (Gram Mode)

Refer to "4. Display Symbols and Key Operation" before operation.

Note When turning the balance on with a tare weight placed on the pan, the balance automatically displays the zero display.

I/O ON:OFF 1 Press the ON:OFF key to turn the display on. Press the key to select a unit, 2 Press the MODE key to select a unit, if necessary. MODE if necessary. 00000 Container 3 Place a container (tare) on the weighing pan, if necessary. (tare) Press the RE-ZERO key to cancel the tare weight. Weighing pan Then zero is displayed. →0← RE-ZERO Tare: A vessel placed on the pan, but not Stabilization to be included in the weighing data. indicator 000000 9 Example: Container. Sample Zero display 4 Place a sample on the pan or in the container. The glass door Close the door. is not closed 150972 <sup>600</sup>5 5 Wait for the stabilization indicator o to be Close the door displayed, and then read the value. Stabilization Warning 190972 9 indicator 6 Remove the sample and container from the pan. Remove

#### Zero operation, tare operation and weighing range

#### When weighing is started

The balance will determine the reference zero point when the ON:OFF key is pressed to enter the weighing mode.

Depending on the load condition at that time, the balance will automatically judge whether to perform zero or tare operation. The condition for determining which is used is "power on zero range", and when power on zero range is exceeded, the tare operation is performed.

#### Re-zero operation

By pressing the RE-ZERO key, the display can be set to zero.

The re-zero operation with the RE-ZERO key will automatically judge whether to perform zero or tare operation.

The condition for determining which is used is "zero range", and when zero range is exceeded, the tare operation is performed.

#### Weighing range

The range that the balance can weigh varies depending on the model.

When the total amount (net weight + tare weight) displayed for each model exceeds the maximum display, E is displayed to indicate that the weighing range is exceeded.

When exceeded in the negative direction, -E is displayed.


Model	Power on zero range	Zero range	-E display range	
BM-5	Approx. ±0.1 g	Approx3.0 g to +0.1 g	Loca than approx 2 a	
BM-5D	Арргох. ±0. г у	Approx3.0 g to +0.1 g	Less than approx3 g	
BM-20	Approx ±2.2 g	Approx 2.0 a to ±0.4 a	Less than approx3 g	
BM-22	Approx. ±2.2 g	Approx3.0 g to +0.4 g		
BM-252	Approx. ±25 g	Approx37 g to +5.0 g		
BM-200	Approx. ±22 g	Approx37 g to +4.4 g	Loca than approve 27 a	
BM-300	Approx. ±32 g	Approx37 g to +6.4 g	Less than approx37 g	
BM-500	Approx. ±52 g	Approx37 g to +10.4 g		

# 6.2. Smart Range

- □ The smart range for the BM-5D and BM-22 consists of the standard range and precision range (high resolution).
- Smart range function
  - The range switches automatically, depending on the value displayed.
  - Pressing the RE-ZERO key allows weighing in the precision range, regardless of the tare value.
  - The range can be fixed to the standard range by pressing the RANGE key.

#### **Example**

- 1 Press the RE-ZERO key to use the precision range. The balance displays zero.
- 2 Place a container (tare) on the weighing pan. When the value of the container (tare) exceeds the precision range, the balance automatically switches to the standard range.
- 3 Press the RE-ZERO key to use the precision range. The balance displays zero.
- 4 Place the weighing sample in the container (tare). When the sample is within the precision range, the weighing value can be read in high resolution.

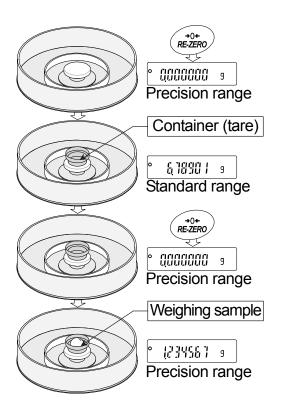

# Precision range and standard range

Model	Unit	Precision range		Standard range	
BM-5D	Milligram	0.000 <b>mg</b> to	2100.009 <b>"9</b>	2100.01 <b>"9</b> to	5200.08 <b>mg</b>
	Gram	0.000000 9 to	2.100009 9	2.10001 9 to	5.20008 9
BM-22	Milligram	0.000 <sup>m</sup> <b>9</b> to	5100.009 <b>"9</b>	5100.01 <b>"9</b> to	22000.08 <b>mg</b>
	Gram	0.000000 9 to	5.100009 9	5.10001 9 to	22.00008 9

Refer to pages 20 and 21 for other units.

# 6.3. Counting Mode (PC)

This is the mode to determine the number of objects in a sample based on the standard sample unit mass. The unit mass means an average mass of the samples. The smaller the variation in the samples is, the more accurate the count will be. The balance is equipped with the Automatic Counting Accuracy Improvement (ACAI) function to improve the counting accuracy.

#### **Notes**

- Use samples with a unit mass of 1 mg or more for counting.
- If the sample unit mass variable is too large, it may cause a counting error.
- □ To improve the counting performance, use the ACAI function frequently or divide the samples into several groups and count each group.

# Selecting the counting mode

1 Press the MODE key to select the unit FC (counting mode).

# Storing a sample unit mass (Weighing input mode)

- 2 Press the RANGE key to enter the sample unit mass storing mode.
- 3 Select the number of samples using the RANGE key. It may be set to 10, 25, 50 or 100.

Note A greater number of samples will yield more accurate counting result.

- 4 Place a container (tare) on the weighing pan, if necessary.

  Press the RE-ZERO key to cancel the weight (tare).


  The number specified in Step 3 appears.
  - Example: 25 0 PC is displayed if 25 is selected in Step 3.
- 5 Place the number of samples specified on the pan. In this example, 25 pieces.
- 6 Wait for the stabilization indicator to be displayed.
  Press the PRINT key to calculate and store the unit mass.
  Then the balance displays 25 PC and is set to count samples with this unit mass. To improve the accuracy of the unit mass, proceed to Step 8.

#### **Notes**

- If the balance judges that the mass of the samples is too light to acquire accurate weighing, it displays an error requiring the addition of more samples to the specified number.
  - Example: 50 PC appears, requiring 25 more samples. Add 25 samples and press the PRINT key. When the unit mass is stored correctly, the balance proceeds to the counting mode.
- If the balance judges that the mass of the samples is too light (under 0.0001g) and cannot be stored as the unit mass, it displays <a href="Lo">Lo</a>.
- The sample unit mass is stored in non-volatile memory, and is maintained even if the AC adapter is removed.

# **Counting operation**

7 Place the samples to be counted on the pan.Read the result and remove the samples from the pan.

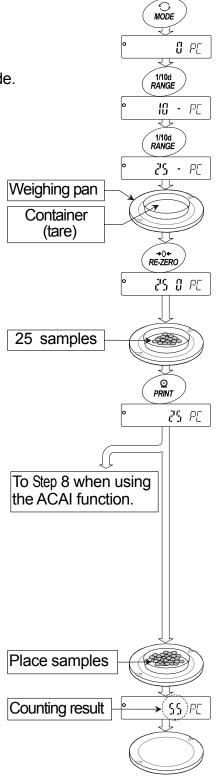

#### Counting mode using the ACAI function

The ACAI is a function that improves the accuracy of the unit mass automatically by increasing the number of samples as the counting process.


**ACAI**: Automatic Counting Accuracy Improvement

Proceed to Step 8 after storing a unit mass on the previous page.

- 8 If a few more samples are added, the processing indicator turns on. To prevent an error, add three or more. The processing indicator does not turn on if overloaded. Try to add the same number of samples as displayed.
- 9 The balance re-calculates the unit mass while the processing indicator is blinking. Do not touch the balance or samples on the pan until the processing indicator turns off.
- 10 Counting accuracy is improved when the processing indicator turns off.

Each time the above operation is performed, a more accurate unit mass will be obtained. There is no definite upper limit to the ACAI range for the number of samples exceeding 100. Try to add the same number of samples as displayed.

11 Remove all the samples used in ACAI and proceed with the counting operation using the improved unit mass.

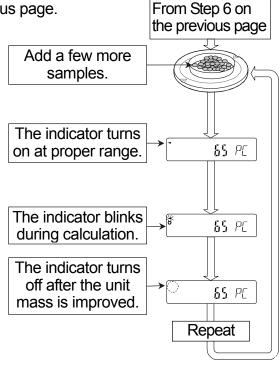

# 6.4. Percent Mode (%)

The percent mode displays the weight value in percentage compared to a 100% reference mass and is used for target weighing or checking the sample variance.

#### Selecting the percent mode

1 Press the MODE key to select the unit % (Percent mode). If the percent mode cannot be selected, refer to "5. Weighing Units".

# Storing the 100% reference mass

- 2 Press the RANGE key to enter the 100% reference mass storing mode.
- 3 Place a container (tare) on the weighing pan, if necessary. Press the RE-ZERO key to cancel the weight (tare). The balance displays 100 0 %.
- 4 Place the sample to be set as the 100% reference mass on the pan or in the container.
- 5 Press the PRINT key to store the reference mass. The balance displays 100,00 %.

#### **Notes**


□ The readability depends on the 100% mass.

BM-252, BM-200 BM-300, BM-500		BM-5, BM-5D BM-20, BM-22		
100% mass	Readability	100% mass	Readability	
0.0100 g to 0.0999 g	1 %	0.00100 g to 0.0099 g	1 %	
0.1000 g to 0.9999 g	0.1 %	0.01000 g to 0.0999 g	0.1 %	
1.0000 g to	0.01 %	0.10000 g to	0.01 %	

- □ If the balance judges that the mass of the sample is too light (under 0.01g) to be used as a reference, it displays \[ \begin{align\*} \lambda \cdot \end{align\*} \].
- The 100% reference mass can be stored in the non-volatile memory and is maintained even if the AC adapter is removed.
- 6 Remove the sample.

# Reading the percentage

7 Place a sample to be compared to the reference mass on the pan. The displayed percentage is based on the 100% reference mass.

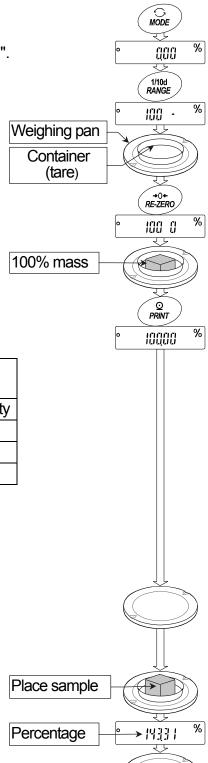

#### **6.5. Ionizer** (Neutralization Device)

The **BM** series is equipped with four ionizer electrode units to neutralize static electricity from a charged weighing sample.

Neutralizing static electricity before weighing will improve stability, reducing errors.

Neutralizing ion: The discharge electrodes of the ionizer continuously generate bipolar ions by corona

discharge. Applying these ions to the weighing sample neutralizes the static charge.

Static electricity: In general, when the ambient humidity is less than 45% RH, nonconductors

such as powders, paper and plastics easily become charged with static electricity. The influence of the static electricity may cause a weighing error of

several milligrams. The ionizer effectively neutralizes the electrical charge.

# Operation

- 1 Place the sample to be weighed on the center (of the metallic circle) of the separation plate to neutralize the static charge.
- 2 Press the ION key to start the neutralization. The ION mark and the LED blink.

  The neutralization will stop after the preset time of "Neutralization time (IDD)" in the function table (the factory setting is 3 seconds).
- □ The readability of the BM-5, BM-5D, BM-20, BM-22 and BM-252 is switched to 0.1 mg while performing the neutralization.
- Pressing the ION key during neutralization stops the neutralization.

# Notes

- Keep the weighing sample away from the electrodes while neutralizing. Placing the sample too close to the electrodes may cause the sample to become charged.
- Remove any obstacle between electrodes and the weighing sample.

# Ionizer Sample Separation plate Press In neutralization Weighing mode QQQQQQQ 9 Weighing mode

Example: Display of BM-20

# 6.5.1. Maintenance of the Electrode Units


- When using the ionizer for a long time, dust and stains may stick to the electrodes. Clean them periodically to maintain performance.
- When needles of the electrodes are worn down and neutralization ability is not restored after cleaning them, replace all four electrodes with new ones. The standard life of an electrode unit is approximately 10000 hours.

# Replacing the electrode units

- 1 Turn the electrode units 45 degrees counterclockwise and remove them from the ionizer.
- 2 Replace all four units with new ones.

Note Do not replace, remove and clean the electrodes while the ionizer is in operation as there is a danger of electrical shock.


3 Turn the electrode units 45 degrees clockwise and insert them into the ionizer.

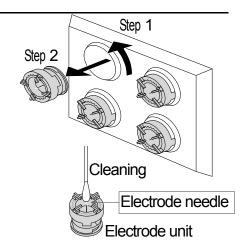

# 6.5.2. Controlling the Ionizer Externally

It is possible to control the BM ionizer function by connecting the remote controller (AD-8922A) or the foot switch (AX-SW137-PRINT, AX-SW137-REZERO).

□ When specifying "□n" to "AD-8922A control (8922)" in the function table (□nFnc 8922 □n), the balance assigns the function of the □ON key to the MODE key on the AD-8922A.

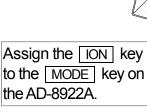

Uhen specifying "External control input (Ε-5)" in "Ionizer function ( μαη Εης)", the following functions can be assigned to AX–SW137–PRINT (option) or AX–SW137–REZERO (option) terminals.

Ionizer function	External control input	AX-SW137-PRINT	AX-SW137-REZERO
■ IonFnc	E-5 O	PRINT key	RE-ZERO key
ionFnc	E-5	PRINT key	ION key
ionFnc	E-5 2	ION key	RE-ZERO key

Factory settings

# 6.6. Monitoring the Clock and Environmental Sensors

In the weighing mode, the built-in clock and environmental sensors (temperature, humidity, barometric pressure) can be monitored and temperature data can be output in the function settings.

Specifications of the environmental sensors


Sensor	Resolution	Measurement range	Applicable range	
Temperature	±1.5 °C	5 °C to 40 °C		
Humidity	±10 %	0 % to 100 %	5 °C to 40 °C	
Barometric pressure	±10 hPa	300 hPa to 1100 hPa		

#### **Notes**

- The environmental sensors are built into the balance casing. Therefore, these values are not in accordance with the environmental condition.
- Refer to "20. Specifications" for the balance operating conditions.

#### Operation

- Press the SELECT key in the weighing mode. The current time, date and sensor values (temperature, humidity, barometric pressure) are displayed for several seconds in order. The balance returns to the weighing mode automatically.
- When pressing the SELECT key while the current data is displayed, the display proceeds to the next data.
- When pressing the CAL key while the current data is displayed, the balance returns to the weighing mode.

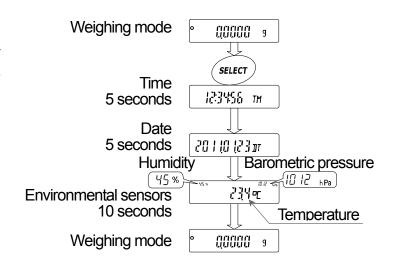

#### **Notes**

- When adjusting the built-in clock, refer to "10.7. Clock and Calendar Function".
- □ When adding the time and date to the output data, refer to "Time/Date output (5-bd)" of "10.2. Details of the Function Table".
- □ When adding temperature data to the output data, refer to "Temperature output  $(5-\ell^p)$ " of "10.2. Details of the Function Table".

# 6.7. Repeatability Test

Repeatability is an index of variation in weighing values when the same mass is repeatedly loaded and unloaded. Standard deviation is generally used for the indicator of repeatability. The repeatability test measures the internal mass 10 times and displays the standard deviation. Use the repeatability test function to inspect the performance of the balance when changing the environment or when moving the balance to a new location.

Example: "Standard deviation = 0.2 mg" means that the results of repeated measurements of the same mass fall within the range of  $\pm 0.2 \text{ mg}$  at a frequency of about 68%.

#### **Notes**


- When the password function is used, the repeatability test function is available only when logged in as an administrator ( ฅժōɪŊ).
- The repeatability test uses the internal mass. Therefore, the obtained standard deviation may disagree with that of the repeatability of "20. Specifications". Regard the standard deviation as a reference.

Model	Internal mass	
BM-5, BM-5D	Approx. 5 g	
BM-20, BM-22	Approx. 20 g	
BM-252, BM-200, BM-300, BM-500	Approx. 190 g	

- 1 Press and hold the RANGE key (for approx. 4 seconds) until rep test is displayed, and then release the key.
- 2 When replaced is displayed, the measurement is automatically started, replaced blinks and the measurement count (0 to 10) is displayed. To cancel the current measurement, press the CAL key.

Note Avoid vibration and breezes for precise measurement during this test.

- 3 At the end of the measurement, the standard deviation of the repeatability is displayed.
- 4 Press the CAL key or PRINT key to return to the weighing mode.

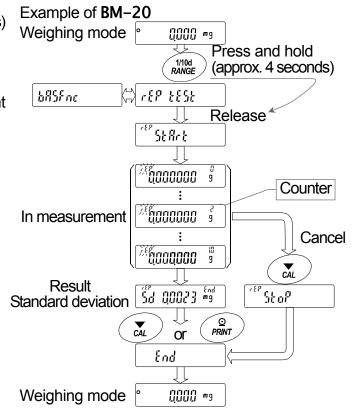


# 7. Response Adjustment

This function stabilizes the weight value by reducing the influence on weighing that is caused by drafts and/or vibration at the place where the balance is installed. The adjustment of the function can be performed by analyzing the environment automatically or hand-operation.

The state of the function has three stages as follows:

Indicator	Parameter	Response	Stability
FAST	Cond O	Fast response,	Sensitive value
MID.	Cond I	1	•
SLOW	Cond 2	Slow response,	Stable value

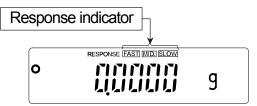

# 7.1. Automatic Response Adjustment

This function automatically updates the response adjustment by analyzing the influence of the environment using the internal mass.


- 1 Press and hold the MODE key (for approx. 2 seconds) until RESPONSE is displayed, and then release the key.
- 2 The balance automatically sets the response characteristic.

# Caution Do not allow vibration or drafts to affect the balance during adjustment.

3 After automatic adjustment, the balance displays *End*, returns to the weighing mode and displays the updated response indicator for about thirty seconds.

#### **Notes**

- If the automatic response adjustment fails, the balance displays LH nL. Check the ambient conditions such as breeze and vibration, also check the weighing pan. Then, perform adjustment again. Press the CAL key to return to the weighing mode.
- If there is anything on the weighing pan, the balance displays [H E]
   Remove the substance from the pan.
   Press the CAL
   key to return to the weighing mode.
- If the automatic response adjustment is not helpful, try"7.2. Manual Response Adjustment".

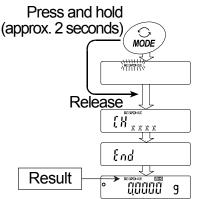

# 7.2. Manual Response Adjustment


- 1 Press and hold the MODE key (for approx. 2 seconds) until RESPONSE is displayed, and then release the key. Press the MODE key again immediately.
- 2 Select a setting for the response adjustment using the MODE key. Either FAST, MID. or SLOW can be selected.
- 3 The balance displays  $\boxed{\mathcal{E}_{nd}}$ , returns to the weighing mode and displays the updated response indicator for about thirty seconds.

#### **Notes**

 When the response adjustment is selected, the "Condition ([and)" parameter of "Environment, Display (bR5Fnc)" in the function table will be changed accordingly as shown below.

Display	[and (Condition)
FAST	0
MID.	1
SLOW	2

□ The response adjustment setting can be changed by the "Condition ([cond)" parameter of "Environment, Display (bff5fnc)" in the function table. For details, refer to "10. Function Table".

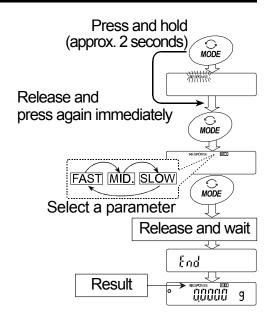

# 8. Calibration

## 8.1. Calibration Group

Since the balance's resolution is high, weighing values may change due to gravity and daily environmental changes. It is necessary to perform calibration (sensitivity adjustment) with the weight in order to keep the weighing values from changing even if gravity or the environment changes.

Calibrate the balance if it is installed for the first time or relocated, or when the weighing values change significantly in daily inspection.

Calibration means to adjust the weighing value of the balance using the reference weight or internal mass. Calibration test means to weigh with the reference weight and compare how much the result deviates from the reference value. (Calibration test does not perform adjustment.)

## Calibration (Sensitivity adjustment)

Automatic self calibration	Automatically adjusts the balance using the internal mass
	depending on the temperature change of the operating
	environment.
Calibration using the internal mass	Using the internal mass, adjusts the balance with a single touch.
Calibration using an external weight	Using an external weight, adjusts the balance.

## Calibration test (Sensitivity check)

Calibration test using the internal mass —— Checks the accuracy of weighing using the internal mass and outputs the result.

\*No adjustment is made.

Calibration test using an external weight— Checks the accuracy of weighing using an external weight (target weight) and outputs the result.

\*No adjustment is made.

#### **Notes**

- Do not allow vibration or drafts to affect the balance during calibration.
- □ Calibration test is available only when "GLP output ( ¬¬¬¬¬)" is set to " /" , "?" or "?".
- □ The calibration and calibration test data can be stored in memory. When using memory, set "Data memory (dRtR)" of the function table. Refer to "12. Data Memory" for details.

## Note on using an external weight

- The accuracy of an external weight can influence the accuracy of weighing.
- Select a weight for calibration and calibration test from the following table.

Model			Usable we	Adjustable range		
BM-5	1 g	2 g	5 g*			
BM-5D	1 g	2 g	5 g*			-3.000 mg to +3.099 mg
BM-20	1 g	2 g	5 g	10 g	20 g*	-3.000 mg to 13.099 mg
BM-22	1 g	2 g	5 g	10 g	20 g*	
BM-252	10 g	20 g	50 g	100 g	200 g*	-15.00 mg to +15.99 mg
BM-200	50 g	100 g	200 g*			
BM-300	50 g	100 g	200 g*	300 g		-30.0 mg to +30.9 mg
BM-500	50 g	100 g	200 g*	300 g	500 g	

**Bold type\***: Factory settings. The weight value can be adjusted within the range above.

#### About the internal mass

The internal mass may change due to corrosion or other damage caused by the operating environment, or due to aging. Check the internal mass periodically and correct the internal mass value if necessary. Refer to "8.7. Correcting the Internal Mass Value".

Display	
◀	This indicator means that the balance is measuring calibration data.
	Do not allow vibration or drafts to affect the balance while the indicator is
	displayed.

# 8.2. Automatic Self Calibration

### Automatic self calibration due to changes in temperature

This function automatically calibrates the balance when the balance detects an ambient temperature change. If GLP output is selected in the function table, the balance outputs the calibration report or stores the data in memory. Automatic self calibration functions even if the display is turned off (standby mode). Refer to "9.1. Permit or Inhibit" for the operation.

#### Caution

- To maintain the calibrated state, keep the weighing pan clear while not in use.
- If something is on the weighing pan, the balance decides that it is in use and does not perform automatic self calibration.
- When performing a long term weighing with something placed on the pan, installing the balance in a system or considering the importance of continuous data, turn off automatic self calibration.


#### **Notes**

	placed on the	on the balance with nothing on the pan, if a sample heavier than 0.5 g is pan, the balance detects the state that a sample is placed on the pan and orm automatic self calibration.
*		Indicates that the balance detects changes in ambient temperature and automatic self calibration will start. If the balance is not used for several minutes with this indicator ( $\triangleleft$ ) blinking, the balance performs automatic self calibration. The blinking duration depends on the operating environment.
<b>▼</b>	LKL	Indicates that the balance is measuring calibration data. Do not allow vibration or drafts to affect the balance while this indicator is displayed. After calibration, the balance returns to indicate the previous display.

The balance can be used while the indicator blinks. But, to maintain the accuracy, stop using the balance and confirm that there is nothing on the pan and allow the balance to perform self calibration.

# 8.3. Calibration Using the Internal Mass

- This function calibrates the balance using the internal mass.
- □ The only operation required is to press the CAL key.
- 1 Connect the AC adapter and warm up the balance for at least one hour with nothing on the weighing pan.
- 2 Press the CAL key to display [FRL 17]
- 3 The balance performs calibration using the internal mass. Do not allow vibration or drafts to affect the balance.
- 4 If GLP output is set,  $\boxed{\text{GLP}}$  is displayed, the calibration report is output to the RS-232C interface and is stored in memory. Refer to "GLP output ( $\neg F_{\alpha}$ )" and "Data memory ( $\exists R \in \mathbb{R}$ )" of the function table in "11.2. GLP Report" and "12. Data Memory".  $\boxed{\mathcal{E}_{\alpha}d}$  is displayed after calibration.
- 5 The balance will automatically return to the weighing mode after calibration.
- 6 Confirm weighing accuracy using calibration test ([[ 17]).

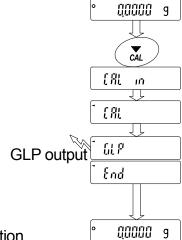

# 8.4. Calibration Test Using the Internal Mass


- This function checks the accuracy of weighing using the internal mass.
- Calibration test does not perform adjustment.
- 1 Connect the AC adapter and warm up the balance for at least one hour with nothing on the weighing pan.
- 2 Press and hold the CAL key until [[[ in ] is displayed, and then release the key.
- 3 The balance measures the zero point.
  Prevent vibration and drafts to affect the balance.
- 4 The measured zero point data is displayed.
- 5 The balance measures the internal mass.
  Prevent vibration and drafts to affect the balance.
- 6 The value of the internal mass is displayed. The normal range of the value is as follows:

Model	Internal mass	Normal range
BM-5, BM-5D	5.00000 g	±0.02 mg
BM-20, BM-22	20.00000 g	±0.02 mg
BM-252, BM-200 BM-300, BM-500	200.0000 g	±0.2 mg

7  $\boxed{\text{GLP}}$  is displayed and the calibration test report is output to the RS-232C interface and is stored in memory. Refer to "GLP output ( InFa)" and "Data memory (dRER)" of the function table.

End is displayed after the calibration test.

8 The balance automatically returns to the weighing mode.

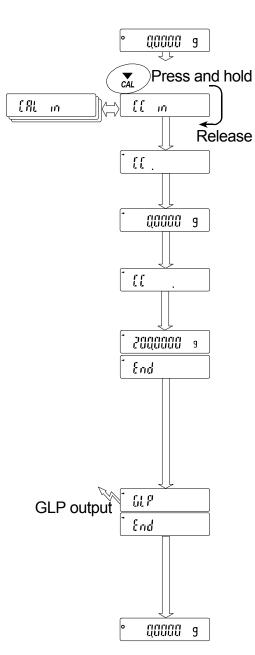

# 8.5. Calibration Using an External Weight

This function calibrates the balance using an external weight.

Note Automatic self calibration (due to changes in temperature) is enabled in the factory setting. Therefore, even after calibration using an external weight, the balance may perform automatic self calibration whenever it detects changes in temperature. When considering the importance of continuous data or managing the balance using the external weight only, refer to "9.1. Permit or Inhibit" to set the function switch to "inhibit"

- 1 Connect the AC adapter and warm up the balance for at least one hour with nothing on the weighing pan.
- 2 Press and hold the CAL key until [FRL out] is displayed, and then release the key.
- 3 When [FRL []] is displayed, proceed as follows:
  - □ If you want to change the calibration weight, press the RANGE key and proceed to Step 4.
  - If you use the calibration weight value stored in the balance, proceed to Step 5.
- 4 Specify the calibration weight value as follows:

RANGE key ......To switch blinking digits.

RE-ZERO (+)key ....To select the calibration weight or adjust the MODE (-)key calibration weight value. Refer to page 37.

PRINT key ......To store the new calibration weight value. Even if the AC adapter is removed, the data is maintained in non-volatile memory.

CAL key ......To cancel the operation and return to

5 Confirm that there is nothing on the pan and press the PRINT key. The balance measures the zero point.


Do not allow vibration or drafts to affect the balance.

The balance displays the calibration weight value.

6 Place the displayed calibration weight on the pan and press the <a href="PRINT">PRINT</a> key. The balance measures the calibration weight. Do not allow vibration or drafts to affect the balance.

7 End is displayed.Remove the weight from the pan.

- 8 If GLP output is set, GLP is displayed and the calibration report is output and stored. Refer to "GLP output ( mFa)" and "Data memory (dRLR)" of the function table in "11.2. GLP Remove the weight Report" and "12. Data Memory".
- 9 The balance will automatically return to the weighing mode.
- 10 Place the calibration weight on the pan and confirm that the value displayed is correct. If it is not within the range, check the ambient conditions such as breeze and vibration, and also check the weighing pan. Then, repeat Steps 2 to 10.

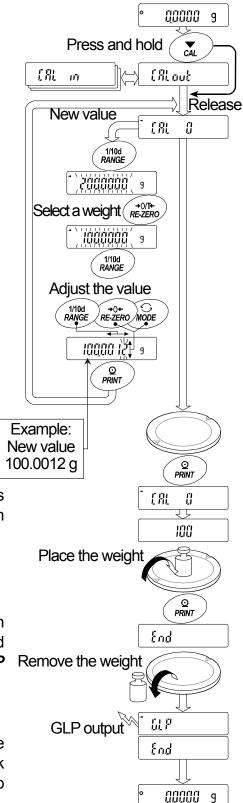

# 8.6. Calibration Test Using an External Weight

- This function checks the accuracy of weighing using an external weight.
- □ This is available only when the "GLP output ( ₁□ F□)" parameter is set to " /", "²" or "³".
- Calibration test does not perform adjustment.
- 1 Connect the AC adapter and warm up the balance for at least one hour with nothing on the weighing pan.
- 2 Press and hold the CAL key until [[ out is displayed, and then release the key.
- 3 When [[[ ]] is displayed, proceed as follows:
  - If you want to change the target weight, press the RANGE key and proceed to Step 4. A list of usable weights is on page 37.
  - □ If the current target weight value is used, proceed to Step 5.
- 4 Specify the target weight value as follows:

RANGE key ...... To switch blinking digits.

RE-ZERO (+)key ... To select the target weight or adjust the MODE (-)key weight value. Refer to page 37.


PRINT key ...... To store the new target weight value.

Even if the AC adapter is removed, the data is maintained in non-volatile memory.

CAL key ..... To cancel the operation and return to


- Confirm that there is nothing on the pan and press the PRINT key. The balance measures the zero point.
   Do not allow vibration or drafts to affect the balance.
   The balance displays the target weight value.
- 6 Place the displayed target weight on the pan and press the PRINT key. The balance measures the target weight. Do not allow vibration or drafts to affect the balance.
- 7 End is displayed.

  Remove the weight from the pan.
- 8 GLP is displayed and the calibration test report is output and stored. Refer to "GLP output (¬¬F¬)" and "Data memory (¬¬F¬)" of the function table in "11.2. GLP Report" and "12. Data Memory".
- 9 The balance will automatically return to the weighing mode.

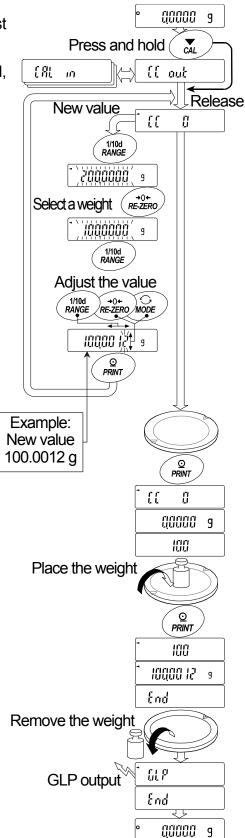

# 8.7. Correcting the Internal Mass Value

The balance can correct the internal mass value within the range shown below. This function corrects the internal mass value to conform to an external weight. The corrected mass value is maintained in non-volatile memory even if the AC adapter is removed. The internal mass value is corrected as follows:

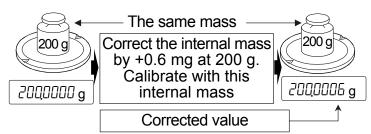

Model	Correction reference value	Correction range
BM-5, BM-5D	5.00000 g	±1.5 mg
BM-20, BM-22	20.00000 g	±1.5 mg
BM-252, BM-200, BM-300, BM-500	200.0000 g	±1.5 mg

**Example**: With the **BM-300**, 200.0000 g is corrected by +0.6 mg (200.0006 g). When performing a +0.6 mg correction with a 100 g external weight, the correction value will be +1.2 mg in accordance with the reference value of 200 g.

- 1 Perform calibration using the internal mass. Weigh an external weight and get the correction value.
- 2 Press the ON:OFF key to turn the display off.
- 3 While pressing and holding the PRINT key and the RANGE key, press the ON:OFF key. P5 is displayed.
- 4 Press the PRINT key to display the function switches. Set the function table switch and internal mass correction switch to " /" as shown to the right using the following keys.

RANGE key.....To select the blinking digit.

RE-ZERO key......To change the value of the blinking digit.

PRINT key.....To store the new value and return to the weighing mode.

CAL key .....To cancel the current operation.

Internal mass correction switch

Function table switch

5 Press and hold the RANGE key (for approx. 2 seconds) to enter the function table and release the key when basen. is displayed.

6 Press the RANGE key several times to display [5 in]

7 Press the PRINT key.

Correct the internal mass value using the following keys.

RE-ZERO (+)key ···· To increase the value.


MODE (-)key ·······To decrease the value.

PRINT key.....To store the new value and display the next

menu item of the function table.

CAL key .....To cancel the current operation and display the next menu item of the function table.

- 8 Press the CAL key to return to the weighing mode.
- 9 Press the CAL key to calibrate the balance using the internal mass.
- 10 Check that the correction has been performed properly with the external weight. If the value is incorrect, repeat the correction.

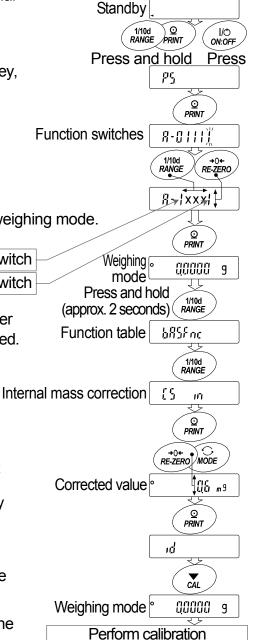

using the internal mass

# **Function Switch and Initialization**

#### Permit or Inhibit 9.1.

- The balance stores parameters that must not be changed unintentionally.
  - Example: Calibration data for accurate weighing, Data for adapting to the operating environment, Control data for the RS-232C interface.
- The balance is equipped with five switches for the purpose of protecting parameters as shown above. Each switch can select either "permit" or "inhibit". "Inhibit" protects parameters against unintentional operations.
  - In addition to the above, when using balances with software version 1.660 or later, the password function can be used to restrict calibration functions.
- 1 Press the ON:OFF key to turn the display off.
- 2 While pressing and holding the PRINT key and the RANGE key, press the ON:OFF key to display
- 3 Press the PRINT key. Then the balance displays the function switches.
- Specify the switches using the following keys.

RANGE key..... To select the blinking digit.

RE-ZERO key...... To change the parameter of the selected switch.

- To inhibit changes. (Cannot be used.)
- To permit changes. (Can be used.)

PRINT key ..... To store the new parameter and return to the weighing mode.

CAL key...... To cancel the operation ( | [L | is displayed ). To return to the weighing mode, press the CAL key once again.

8-81111

(The display shown left indicates the factory settings)

#### Function table

- To inhibit changes to the function table.
  - To permit changes to the function table.

#### Calibration using the internal mass

- Software ver. 1.00 to 1.64 To inhibit calibration using the internal mass. Software ver. 1.660 or later **PRSS** [] To inhibit calibration using the internal mass. To inhibit calibration using the internal mass when
  - See Note 1. logged in as a user (USER) or a guest (UUEST).
  - To permit calibration using the internal mass.

#### Calibration using an external weight

- Software ver. 1.00 to 1.64 To inhibit calibration using an external weight. Software ver. 1.660 or later PR55 [] To inhibit calibration using an external weight. **PRSS** 1.2 To inhibit calibration using an external weight when See Note 1. logged in as a user (USER) or a guest (UUEST).
  - To permit calibration using an external weight.

#### **Automatic self calibration** (for variation of ambient temperature)

- To inhibit automatic self calibration.
- To permit automatic self calibration.

#### Internal mass correction

- To inhibit correction.
- To permit correction.


**Note 1**: The operation is permitted when logged in as an administrator ( $\mathcal{H}d\bar{n}_{IN}$ ).

# 9.2. Initializing the Balance

- This function restores the following parameters to the factory settings.
  - Calibration data
  - Function table (excluding the time and date)
  - Sample unit mass value (counting mode),
     100% reference mass value (percent mode)
  - The data that is stored in the balance using the data memory function
  - External weight value
  - Function switch settings ("9.1. Permit or Inhibit")
  - Liquid density and temperature in the density mode

Note Be sure to calibrate the balance after initialization.

- 1 Press the ON:OFF key to turn the display off.
- 2 While pressing and holding the PRINT key and the RANGE key, press the ON:OFF key to display P5
- 3 Press the RANGE key to display [[Lr]
- 5 Press the RE-ZERO key to display [Lr [ig]
- 6 Press the PRINT key to initialize the balance.
  The balance will automatically return to the weighing mode.

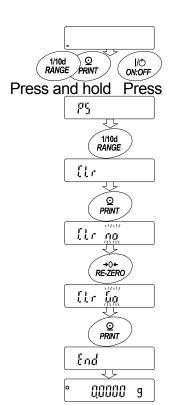

# 10. Function Table

The function table reads or rewrites the parameters that are stored in the balance.

These parameters are maintained in non-volatile memory, even if the AC adapter is removed.

The function table menu consists of two layers. The first layer is the "Class" and the second layer is the "Item".

# 10.1. Setting the Function Table

## Display symbol and keys

<b>o</b>	The symbol "o" indicates that the parameter displayed is in effect.
1/10d RANGE	When pressed and held (for approx. 2 seconds) in the weighing mode, enters the function table mode.  Selects the class or item in the function table mode.
→0← RE-ZERO	Changes the parameter.
Q PRINT	When a class is displayed, enters an item in the class. When an item is displayed, stores the new parameter and displays the next class.
CAL	When an item is displayed, cancels the new parameter and displays the next class. When a class is displayed, exits the function table mode and returns to the weighing mode.


## Setting procedure

- 1 In the weighing mode, press and hold the RANGE key (for approx. 2 seconds) until bR5Fnc is displayed, and then release the key.
- 2 Press the RANGE key to select a class.
- 3 Press the PRINT key to enter the class.
- 4 Press the RANGE key to select an item.
- 5 Press the RE-ZERO key to select a parameter for the selected item.
- 6 Repeat Steps 4 and 5 to select other items of the class.
- 7 To store the parameters of the selected class, press the PRINT key.
  Then the next class is displayed.
  To cancel the current operation, press the CAL key. Then the next class is displayed.
- 8 To specify parameters for another class, proceed to Step 2.

  To finish the setting, press the CAL key to return to the weighing mode.

### Setting example

This example sets the "Data memory (dRER)" parameter to "I" (To store the weighing data) and the "Interval time (IRE)" parameter to "5" (1 minute).

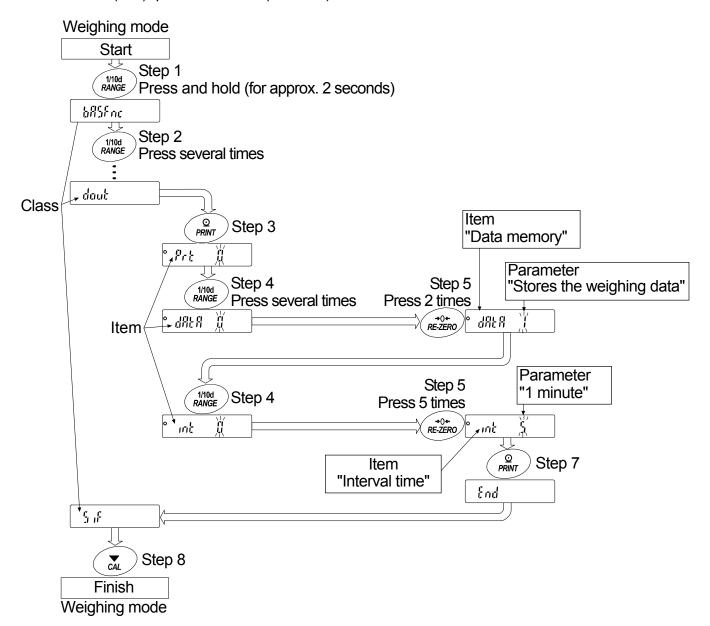

# 10.2. Details of the Function Table

Class	Item and Parameter		Description		
	Eand Condition	• I	Fast response, sensitive value  Slow response, stable value	Can be changed by response adjustment.	
	5と-6 Stability band width	• I	Stable range is ±3 digits	The stabilization indicator illuminates when the display fluctuation is within the range.	
	PL oc Display lock function	■ oFF on	OFF ON	Display lock function to measure the volume dispensed from the pipette.	
₽₩₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽₽	לרכ Zero tracking	• I 2	OFF Normal Strong Very strong	Keeps zero display by tracking zero drift.	
	5P님 Display refresh rate	<b>-</b> []	5 times/second 10 times/second	Period to refresh the display	
	Pnt Decimal point	<b>-</b> []	Point (.) Comma (,)	Decimal point format	
	P-an Auto display-ON	<b>■</b> aFF an	OFF ON	Turns on the weighing mode display when the AC adapter is connected.	
	ьеер	aFF ■ an	OFF ON	Sounds a beep when a key is touched.	
Clock	Refer to "10.7. Clo	ck ar	nd Calendar Function"	The time and date are added to the output data.	
	Prt Data output mode	<b>-</b> []	Key mode	Accepts the PRINT key only when the display is stable.	
		 2	Auto print mode A Reference = zero Auto print mode B Reference = last stable value	Outputs data when the display is stable and conditions of #P-P, #P-b and the reference value are met.	
		3	Stream mode / Interval mode	With dfltfl [], outputs data continuously; with dfltfl l, uses interval memory.	
dอนt Data output	RP-P Auto print polarity	• 0 !	Plus only Minus only Both	Displayed value > Reference Displayed value < Reference Regardless of displayed value	
	RP-b Auto print difference		10 digits 100 digits 1000 digits	Difference between reference value and displayed value #2	
	ਰਸ਼ੁਸ਼ Data memory	• 0 ! 2	Not used Stores the weighing data Stores the calibration data	Related items: Prt, int, d-na, 5-td, inFa	

<sup>• :</sup> Factory settings.

Example: If 1 mg display is selected using the RANGE key for the BM-300, 1 mg is one digit.

Example: In gram display, one digit is 0.01 mg for the BM-252 and 0.1 mg for the BM-300.

<sup>#1</sup> The unit of readability is digit.

<sup>#2</sup> Usable readability of the balance is one digit.

Class	Item and Parame	eter	Description		
	<b>•</b> ()		Every measurement		
			2 seconds		
		2	5 seconds		
		3	10 seconds	Interval time in the interval	
	int	4	30 seconds	memory mode when using	
	Interval time	5	1 minute	PrŁ 3, dRŁR I	
		5	2 minutes		
		7	5 minutes		
		8	10 minutes		
	d-na	<b>■</b> oFF	No output	- C	
	Data number output		Output	Refer to "12. Data Memory"	
	'	<b>-</b> []	No output	Selects whether or not the time or	
	5-Ed		Time only	date is added to the weighting	
	Time/Date output		Date only	data. Refer to "10.7. Clock and	
dout		<u>-</u>	Time and date	Calendar Function" for details.	
Data output	5- 18		No output	Selects whether or not the ID	
	ID number output		<u> </u>	number is output.	
		<b>■</b> oFF	No pause	·	
	Data output pause		Pause (1.6 seconds)	Selects the data output interval.	
	RE-F	■ oFF	Not used	Selects whether or not automatic	
	Auto feed		Used	feed is performed.	
	7 1010 1000	<b>-</b> []	No output	Refer to "11. ID Number and	
	GLP output	<u>-</u>	AD-8121 format	GLP Report" for details.	
		<u></u>		#3 Available for balances with software	
		3	General (external device clock data) #3	version 1.660 or later.	
	Rr - d	<u>-</u> _FF	Not used	Adjusts zero automatically after	
	Zero after output		Used	data is output	
	5- <i>LP</i>		Not output		
	Temperature output	ص م		Adds temperature to data output	
	, , , , , , , , , , , , , , , , , , ,	0	600 bps		
			1200 bps		
	6PS	• 2	2400 bps		
	Baud rate	3	4800 bps		
		4	9600 bps		
		5	19200 bps		
		<b>-</b> []	7 bits, even		
S iF	占ととこ Data bit, parity bit	1	7 bits, odd		
Serial	Data Dit, Parity Dit	2	8 bits, none		
interface	[rLF	<b>-</b> []	CR LF	CR: ASCII code 0Dh	
	Terminator		CR	LF: ASCII code 0Ah	
		<b>-</b> []	A&D standard format		
		1	DP format		
	EYPE Data format	2	KF format	Refer to "10.5. Description of	
		3	MT format	Refer to "10.5. Description of the Item "Data Format" ".	
		<del>-</del> 4	NU format		
		<u>'</u>	CSV format	} 	
			nctony softings	<u>I</u>	

■ : Factory settings.

The balance may not transmit the data completely at the specified refresh rate, depending on the baud rate or data added to the weighing data such as time, date and ID number. Caution

Class	s Item and Parameter		Description		
5 iF	는 - UP Timeout	oFF ■ on	No limit 1 second	Selects the wait time to receive a command.	
Serial	Er[d		No output	a command.	
interface	AK, Error code	-u''	Output	AK: ASCII code 06h	
dS Fnc	,	<b>-</b> ()	Water temperature	A 11.1. 1 1 1 1	
Density	Ld in Liquid density input			Available only when density mode is selected	
function	Liquid del isity il iput	1	Liquid density	mode is selected	
Աո ₁Է Unit			Refer to "5. Weighing Units".		
[5] Internal mas	e correction		Displayed only when the inter switch is set to 1. Refer to "8.		
	3 COITECTION		SWITCH IS SET TO T. INCICL TO U.	Cambration .	
ਾਰ ID number s	etting		Refer to "11. ID Number and	GLP Report".	
		<b>-</b> []	3 seconds		
	Neutralization time	1	10 seconds	Refer to "6.5. lonizer".	
_	TVCdtt diiZdtiOff tiffic	2	Manual operation (max. 10 min.)		
Ionizer	8922	■oFF	The ION key is not used.		
function	AD8922A control	٦٥	Assigns the ON key to the	MODE key on AD-8922A.	
Tariotion	E-5 External control input	<b>-</b> []	Assigns the PRINT key and RE-ZERO key.		
			Assigns the PRINT key and ION key.		
	External control input	2	Assigns the ON key and	RE-ZERO key.	
door	d-dE	oFF	Not used		
Door sensor function	Door sensor	<b>.</b> on	Used	Refer to "6.1. Basic Operation"	
LocFnc	0000	<b>.</b> []	Not used	Defects #15 December	
Password	PR55 #4 Password function	1	Used (Limits weighing operations)	Refer to "15. Password Function".	
function	i assword fullcillit	2	Used (Allows basic weighing)	i diledoli .	
	Adāj).		Password input by the administrator		
Lac na.	Loc no.		Password input by user 1		
Password			i		
		USE <b>R</b> 10	Password input by user 10	#5	

#### : Factory settings.

- #4 Only "Not used ([])" and "Used ( |)" are available for balances with software version 1.50 to 1.64. All the parameters are available for balances with software version 1.660 or later.
- #5 Available for balances with software version 1.660 or later.

  For balances with software version 1.50 to 1.64, only the password input by the administrator is available.

## 10.3. Description of the Class "Environment, Display"

## Condition ([and)

This parameter is for sensitive response to the fluctuation of a mass value. Used for powder target weighing, weighing a very light sample or when quick response weighing is required. After setting, the balance displays FAST.

This parameter is for stable weighing with slow response. Used to prevent a weight value from drifting due to vibration or drafts. After setting, the balance displays SLOW.

Note In automatic response adjustment, this parameter is selected automatically.

### Stability band width (5L-b)

This item controls the width to regard a mass value as a stable value. When the fluctuation per second is less than the parameter, the balance displays the stabilization indicator and outputs or stores the data. The parameter influences the "Auto print mode".

- 5k-b 0 This parameter is used for sensitive response of the stabilization indicator. Used for exact weighing.
- 5½-b 2 This parameter ignores slight fluctuations of a mass value. Used to prevent a mass value from drifting due to vibration or drafts.

### Display lock function (PLoc)

- This function is used to lock the display when measuring the volume dispensed from the pipette. With "Display lock function is used (PLac an)" selected in the function table, when the liquid is dispensed into the weighing bottle from the pipette and a stable weighing value is obtained, the balance calculates an average value and displays the result (locked).
- When using this function for the measurement of a small volume, the influence of liquid evaporation can be reduced.

Caution Even if the weighing bottle (evaporator trap) to reduce the water evaporation is used, at least 0.05 mg (0.05  $\mu\ell$ ) evaporation will occur.

□ When canceling this function, press the RE-ZERO key.



### **Zero tracking** (*Lrc* )

This function tracks zero point drift caused by changes in the environment and stabilizes the zero point. When the weighing data is only a few digits, turn the function off for accurate weighing.

- **Lrc** Π The tracking function is not used. Used for weighing a very light sample.
- Erc | The normal tracking function is used.
- *Erc ≥* The strong tracking function is used.
- *Lrc* ∃ The very strong tracking function is used. Used for stable zero display.

## Display refresh rate (5Pd)

The periodic time to refresh the display. This parameter influences "Baud rate", "Data output pause" and the data output rate of "Stream mode".

## Decimal point (Pnt)

The decimal point format can be selected.

## Auto display-ON (P-an)

When the AC adapter is plugged in, the display is automatically turned on without pressing the ON:OFF key, to display the weighing mode. Used when the balance is built into an automated system. One hour warm up is necessary for accurate weighing.

# 10.4. Description of the Item "Data Output Mode"

The parameter setting of "Data output mode (PrE)" applies to the performance when the "Data memory (dRER)" parameter is set to "I" (to store the weighing data) and when the data is transmitted using the RS-232C interface.

### Key mode

When the PRINT key is pressed with the stabilization indictor turned on, the balance outputs or stores the weighing data and the display blinks one time.

Required setting doub Prt (1) Key mode

### Auto print modes A and B

When the displayed value is stable and the conditions of "Auto print polarity", "Auto print difference" and the reference value are met, the balance outputs or stores the weighing data.

When the PRINT key is pressed with the stabilization indictor turned on, the balance outputs or stores the data and the display blinks one time.

## Auto print mode A

Example For weighing each time a sample is placed and removed, with "8r-d" set

to " /" (to adjust zero after the data is output).

Required setting doub Prb | Auto print mode A (reference = zero)

dout RP-P Auto print polarity
dout RP-b Auto print difference
dout Rr-d on Zero after output

### Auto print mode B

Example For weighing while a sample is added.

Required setting doub Prt 2 Auto print mode B (reference = last stable value)

dout RP-P Auto print polarity
dout RP-b Auto print difference

#### Stream mode

The balance outputs the weighing data continuously regardless of the display condition. When the display refresh rate is set to 5 times / second (5Pd  $\Box$ ), the data output rate is also set to the same 5 times / second. The display does not blink in this mode. The interval memory mode is used when the "Data memory (dRER)" parameter is set to " l" (to store the weighing data).

Example For monitoring data on a computer.

Required setting doub Prt 3 Stream mode

ժոսե ժRER 0 Data memory function is not used

bRSFnc SPd Display refresh rate

5 F 6P5 Baud rate

Caution The balance may not transmit the data completely at the specified refresh rate, depending on the baud rate or data added to the weighing data such as time, date and ID number.

#### Interval mode

The balance outputs the weighing data periodically regardless of the display condition and the display blinks one time. When the "Data memory (dfltfl)" parameter of "Data output (daut)" is set to "!" (to store the weighing data), the weighing data is periodically stored in memory.

Pressing the PRINT key starts output and pressing the PRINT key again during output stops the output.

Example For periodical weighing without a personal computer command and to

output all of the data, to a computer, at one time.

Required setting dout Prt 3 Interval mode

לפטע שאבא Data memory function is used

dout interval time

# 10.5. Description of the Item "Data Format"

#### A&D standard format

S IF LYPE D

This format is used when the peripheral equipment can receive the A&D format.

With an AD-8121B printer, use MODE 1 or 2. With an AD-8127 printer, use the external key printing mode, manual printing mode, auto printing mode or interval printing mode.

- □ This format consists of fifteen or sixteen characters excluding the terminator.
- A header of two characters indicates the balance condition.
- □ The polarity sign is placed before the data with the leading zeros. If the data is zero, the plus sign is applied.
- The unit, consisting of three characters, follows the data.

S	Т	,	+	0	0	0		1	2	7	8		ட	g	CR	LF						
	He	ade	er			D	ata						Ur	nit		Te	rmiı	nato	or			
S	Т	St	abl	e h	eac	der				(	Q	Т	S	Stat	ole I	nea	der	of	СО	unt	ing r	node
U	S	U	nsta	able	e he	ead	ler															
0	L	0	ver	loa	d h	eac	der															
S	Т	,	+	1	0	0		0	1	2	7	8		ш	g	$c_R$	LF					
H	lead	der		8	-nu	me	erica	al c	hai	act	er o	data	ì	Ur	nit		Ter	mir	nat	or		


## DP (Dump print) format

S IF EYPE I

This format is used when the peripheral equipment cannot receive the A&D format.

With an AD-8121B printer, use MODE 3. With an AD-8127 printer, use the dump printing mode.


- This format consists of sixteen characters excluding the terminator.
- A header of two characters indicates the balance condition. No overload header is used.
- □ The polarity sign is placed before the data, with spaces in place of leading zeros, if the data is not zero or overloaded.
- The unit, consisting of three characters, follows the data.

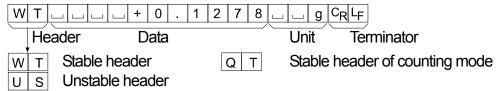

KF format 5 iF EYPE 2

This is the Karl-Fischer moisture meter format and is used when the peripheral equipment can only communicate using this format.

- This format consists of fourteen characters excluding the terminator.
- This format has no header characters.
- The polarity sign is placed before the data, with spaces in place of leading zeros, if the data is not zero or overloaded.
- This format outputs the unit only for a stable value.

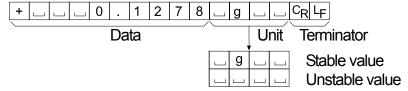

MT format 5 if LYPE 3

This format is used when the peripheral equipment of other manufacturer is connected. Please note that the connection is not guaranteed.

- A header of two characters indicates the balance condition.
- The polarity sign is used only for negative data.
- The weighing data uses spaces in place of the leading zeros.
- The character length of this format changes dependent upon the unit


S				_ 0		1	2	7	8	J	g	C <sub>R</sub> I	-F				
`	He	ader		D	ata				/	Ur	nit	Te	rmin	ato	r		
S		Stab	le he	ader													
S	D	Unst	able	head	der												
S	I	Ove	rload	head	der												
ш		Stab	le he	ader	(W	he	n d	ata	is c	outc	ut	usin	a the	) F	PRINT	٦١	(ev

### **NU** (numerical) format

5 if EYPE 4

This format outputs only numerical data.

- □ This format consists of ten characters excluding the terminator.
- □ The polarity sign is placed before the data with the leading zeros. If the data is zero, the plus sign is used.

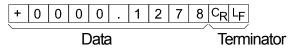

CSV format 5 if EYPE 5

- □ This format separates the data of A&D standard format and the unit by a comma ( , ).
- □ This format outputs the unit even when the data is overloaded.
- □ When a comma ( , ) is selected for the decimal point, the separators are set to semicolon ( ; ).
- □ When the ID number, data number, time and date are added in "Data output (daut)" of the function table, outputs ID number, data number, date, time and weighing data in this order and separates each item by a comma and treats all the items as one group of data.

```
LAB-0123, No,012, 2010/11/01, 12:34:56, ST,+0000.1278,...g<CR><LF>
| D number | Data number | Date | Time | Weighing data | | | | | | | | | | | | | |
| S | T | , | + | 0 | 0 | 0 | . | 1 | 2 | 7 | 8 | , | ... | g | C<sub>R</sub> L<sub>F</sub> |
| O | L | , | + | 9 | 9 | 9 | 9 | 9 | 9 | E | + | 1 | 9 | , | ... | g | C<sub>R</sub> L<sub>F</sub> |
```

Note When the balance is connected to a personal computer via the USB interface, the ID number, data number, time and date cannot be added to the weighing data.

ID number dout 5- id on

The number to identify a specific balance.


□ This format consists of eight characters excluding the terminator.

L A B - 0 1 2 3 C<sub>R</sub>L<sub>F</sub>

Data number doub done on

This format outputs the data number just before the data is transmitted using the RS-232C interface.

- □ This format consists of six characters excluding the terminator.
- □ When CSV format (5 ,F Ł YPE 5) is selected, the period ( . ) is replaced with a comma ( , ).

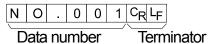

Date

dout 5-td 2 or 3

□ The date output order can be changed in "Time/Date output (5-½d)" and "Clock (½ ਕਰਹ)". This format outputs the year in a four-digit format.

```
2 0 1 0 / 1 1 / 0 1 C<sub>R</sub>L<sub>F</sub>
```

Time

dout 5-td | or 3

□ This format outputs the time in a 24-hour format.

```
1 2 : 3 4 : 5 6 C<sub>R</sub>L<sub>F</sub>
```

**Temperature** 

dout 5-tP on

□ This format outputs the temperature. The unit is "C".

Note Humidity and barometric pressure are not output.

```
+ 0 2 3 . 4 _ C C<sub>R</sub>L<sub>F</sub>
```

# 10.6. Data Format Examples

## Stable

° 01218 9

S	Т	,	+	0	0	0		1	2	7	8	ш		g	$C_{R}$	LF	
S	Т	,	+	٦	0	0	•	0	1	2	7	8	J	J	g	$c_R$	LF
W	Т		J			+	0		1	2	7	8			g	$c_R$	LF
+	Ţ	]	]	0		1	2	7	8		g			$c_R$	LЕ		
S	Ţ	]	J	]		0		1	2	7	8		g	$c_R$	LF		
+	0	0	0	0		1	2	7	8	$C_R$	LБ						

## Unstable

- 183690 g

U	S	,	_	0	1	8		3	6	9	0			g	$C_{R}$	LF	
U	S	,	1	1	0	1		8	3	6	9	0			g	$c_R$	LF
U	S		]		-	1	8		3	6	9	0			g	$c_R$	ᅱ
-	]	]	1	8		3	6	9	0	]				$c_R$	۱F		
S	О	]	]	ı	1	8		3	6	9	0		g	$c_R$	LF		
-	0	0	1	8		3	6	9	0	$C_{R}$	LF						

# Overload

Positive error

**£** 9

	0	L	,	+	9	9	9	9	9	9	9	Е	+	1	9	$C_{R}$	LF	
ι	L			]	]				Е						J		$C_{R}$	LF
ı	L	]					Н				ш	ш	ш		$c_R$	LF		
	ഗ	ı	+	$c_{R}$	ᆦ													
	+	9	9	9	9	9	9	9	9	9	$c_R$	۱F						

## Overload

Negative error

**- ⊱** g

0	L	,	-	9	9	9	9	9	9	9	Е	+	1	9	$C_{R}$	LБ	
							-	Е	J	]	]	J	]			$c_R$	봔
ш						L	]			]		]	]	$c_R$	LБ		
S	ı	-	$c_R$	LF													
-	တ	တ	9	9	9	9	9	9	9	$C_{R}$	LF						

 $\begin{tabular}{lllll} $\square$ & Space, & ASCII & 20h \\ $^{C}_{R}$ & Carriage Return, & ASCII & 0Dh \\ $^{L}_{F}$ & Line Feed, & ASCII & 0Ah \\ \end{tabular}$ 

# Units

		A&D	D.P.	KF	MT
g	9		g		□ g
mg	™g	m   g	m   g	$\square$ m g $\square$	$\square$ m g
Counting mode	PE	□ P C	□ P C	□ p c s	□ P C S
Percent mode	%	<u> </u>	<b>□</b> □ <b>%</b>	<u></u>  %  _   _	<u></u> %
Ounce (Avoir)	OΖ	0 Z	_ 0 Z	⊔ O Z ⊔	⊔ O Z
Troy Ounce	0Zt	o z t	o z t	⊔ o z t	⊔ o z t
Metric Carat	ct	∟ c t	∟ c t	□ c t □	□ c t
Momme	mom	m o m	m o m	$\square m o m$	□ m o
Pennyweight	dnt	d w t	d w t	d w t	ᆸ d w t
Grain	БN	□GN	⊔GN	g r	uGN
Tael (HK general, Singapore)	EL	山 t I	ഥ t l	ᆸ t l s	山 t l
Tael (HK, jewelry)	EL	山 t l	山 t l	山 t l h	∟ t I
Tael (Taiwan)	EL	山 t l	ш t I	⊔ t l t	山 t l
Tael (China)	EL	山 t l	ഥ t l	⊔ t I c	山 t l
Tola (India)	t	t	t	니 t O I	山 t
Messghal	m	m e s	m e s	uМSu	<u></u> m
Density	115	□ D S	□ D S		□ D S

□ Space, ASCII 20h

# 10.7. Clock and Calendar Function

The balance is equipped with a clock and calendar function. When "GLP output ( יח २ ם)" and "Time/Date output (5 - ৮ d)" are used, the time and date can be added to the output data. The operation is as follows:

00000 9 Press and hold

basfac

#### Operation

- (approx. 2 seconds) 1 In the weighing mode, press and hold the RANGE key (for approx. 2 seconds) until basence is displayed, and then release the key. Press several times
- 2 Press the RANGE key several times to display [[L RdJ]].
- 3 Press the PRINT key to enter the time and date setting mode.

## Checking the current time

- 4 The current time is displayed with all the digits blinking.
  - □ To change the time, press the RE-ZERO key and proceed to Step 5.
  - □ When the time is correct, press the RANGE key and proceed to Step 6.
  - □ To exit the mode, press the CAL key and proceed to Step 8.

## Storing a new time

5 Set the time in a 24-hour format using the following keys.

RE-ZERO (+)key... To increase the value by one.

MODE (-)key ...... To decrease the value by one.

RANGE key...... To select the digit to change the value.

PRINT key..... To store the new time and proceed to Step 6.

CAL key..... To cancel the operation and proceed to Step 6.

## Checking the current date

- 6 The current date is displayed with all the digits blinking.
  - $\Box$  To change the display order of year ( $\Box$ ), month ( $\bar{\cap}$ ) and day (d), press the MODE key. The date is output in the order as specified.
  - □ To change the date, press the RE-ZERO key and proceed to Step 7.
  - □ When the date is correct, press the CAL key and proceed to Step 8.
  - □ To change the time again, press the RANGE key and proceed back to Step 4.

## Storing a new date

Set the date using the following keys.

RE-ZERO (+)key ·· To increase the value by one.

MODE (-)key ...... To decrease the value by one.

RANGE key ...... To select the digits to change the value.

PRINT | key ...... To store the new date and proceed to Step 8.


CAL key ..... To cancel the operation and proceed to Step 8.

## Quitting the operation

8 The balance displays the next menu item of the function table. Press the CAL key to exit the mode and return to the weighing mode.

Note Do not enter invalid values such as non-existing time and date in this function. When the clock backup battery has been depleted, the balance displays | rec PF |. Under this condition, press any key and set the time and date. The dead battery only affects the clock and calendar function. Even so, the function works normally as long as the AC adapter is connected to the balance.

58

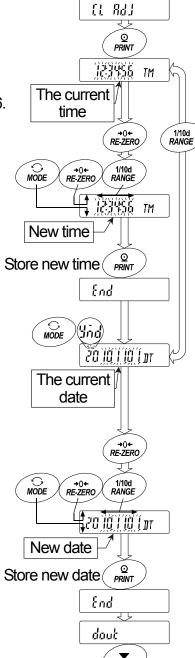

# 11. ID Number and GLP Report

- The ID number is used to identify the balance when Good Laboratory Practice (GLP) or Good Manufacturing Practice (GMP) is used.
- The output format for GLP/GMP compliant report is selected in "GLP output (nFa)" of the function table and can be output to a personal computer or printer using the RS-232C serial interface.
- The GLP/GMP compliant report includes the balance manufacturer, model, serial number, ID number, date, time and space for signature. The calibration report and the calibration test report include the weight used and the result.
- □ The balance can output the following reports for GLP / GMP using the RS-232C serial interface.
  - "Calibration report" of the automatic self calibration and the calibration using the internal mass.
  - "Calibration report" of the calibration using an external weight.
  - "Calibration test report" of the calibration test using the internal mass.
  - "Calibration test report" of the calibration test using an external weight.
  - "Title block" and "End block" for the weighing data.
- Calibration report and calibration test report can be stored in memory and several reports can be output at the same time. Refer to "12. Data Memory" for details.
- The ID number is maintained in non-volatile memory even if the AC adapter is removed.
- Refer to "10.7. Clock and Calendar Function" for checking and storing the time and date.
- When an AD-8127 printer is connected to the balance and outputs the GLP report, the time and date can be added to the report using the clock function of the printer. (Available for balances with software version 1.660 or later) This is very useful for managing the time and date using the AD-8127 password lock function to prevent them from being falsified.

To output the GLP/GMP compliant report, set the AD-8127 printer to the dump printing mode.

# 11.1. Setting the ID Number

	Title Setting the 15 Manibel
1	Press and hold the RANGE key (for approx. 2 seconds) until BRSFnc is displayed, and ther release the key.
2	Press the RANGE key several times to display .
3	Press the PRINT key. Set the ID number using the following keys.  RE-ZERO key·····To set the character of the digit selected.  Refer to the display character table shown below.
	RANGE key······To select the digit to change the character.  PRINT key······To store the new ID number and display basen.

4 When bR5Fnc is displayed, press the CAL key to return to the weighing mode.

CAL key.....To cancel the new ID number and display 6855 nc

Display character table

0	1	2	3	4	5	6	7	8	9	-	ш	Α	В	С	D	Ε	F	G	Н	I	J	K	L	М	N	0	Ρ	Q	R	S	Т	U	V	W	Χ	Υ	Z
[]	1	2	3	4	5	5	7	8	3	-		R	Ь	Ľ	₫	E	F	ប្រ	Н	ı	ц	2	<u>L</u>	ñ	<i>[]</i>	٥	P	7	r	5	Ŀ	Ц	Ľ,	u -	11	4	<u>,</u>
													Sr	ac	e																						

## 11.2. GLP Report

To output the GLP/GMP compliant report to an AD-8121B printer, AD-8127 printer or to a personal computer, set the "GLP output ( \( \in F\_a \) \)" parameter of the function table to " \( \in \)" (AD-8121 format), "\( \in \)" (General format using the balance built-in clock data) or "\( \in \)" (General format using the external device clock data).

### Notes on outputting the GLP/GMP compliant report to a printer

- □ Refer to "16.2. Connecting to Peripheral Equipment" for connection to a printer.
- □ With an AD-8121B printer, use MODE 3. If MODE1 is used, select temporarily the dump print mode by pressing the STAT. key of the AD-8121B printer.


  With an AD-8127 printer, use the dump printing mode. If the external key printing mode is used, press and hold the ENT key of the AD-8127 printer (for approx. 2 seconds) to switch between the external key printing mode and the dump printing mode. If the data memory function is used (except dRER □), the GLP output cannot be output automatically.
- □ If the time and date are not correct when the balance built-in clock data is output ( ,□F□ 1, 2), set the correct time and date in "Clock (EL RdJ)" of the function table.
- □ " יח + □ ∃" is available for balances with software version 1.660 or later.

## Calibration report using the internal mass

Note If the data memory function is used (except dfltfl (1)), the "Title block" and "End block" cannot be output.

### Key operation

- 1 Press the CAL key to display [FRL 17]. The balance performs calibration using the internal mass.
- 2 If GLP output is used, LGLP is displayed and the calibration report is output.
- 3 The balance returns to the weighing mode.

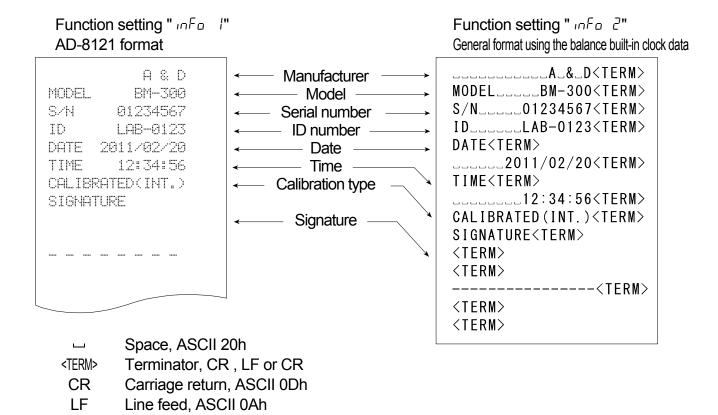


#### Outputting the clock data of the external device " ¬F□ ∃"

When the "GLP output (mFa)" parameter of the function table is set to " $\exists$ ", the clock data of external devices such as a printer or a personal computer can be used instead of the balance built-in clock data for the GLP/GMP compliant report. This is useful for standardizing the clock data using the clock function of an external device.

#### **Notes**

- To output the clock data of an external device, the external device such as an AD-8127 printer or Windows communication tools software WinCT RsCom must have a function to output the time and date when <ESC>D or <ESC>T is received.
- □ Even with InFa 3 set, the data memory function saves the balance built-in clock data for the calibration report.

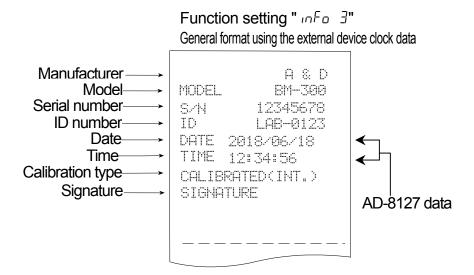

### Calibration test report using the internal mass

Note 

Calibration test does not perform adjustment.

□ If the data memory function is used (except dfltfl 0), the "Title block" and "End block" cannot be output.

#### **Key operation**

- 1 Press and hold the CAL key to display [[[ In ]]] and release the key.
- 2 [[ is displayed and the balance is checked automatically.
- 3 The zero point is measured and the weight value is displayed for a few seconds.
- 4 The internal mass is weighed and the weight value is displayed for a few seconds.
- 5 If GLP output is used, [GLP] is displayed and the calibration test report is output.
- 6 The balance returns to the weighing mode.

#### Command


CR

LF

This calibration test can be performed with the TST command.

Carriage return, ASCII 0Dh

Line feed, ASCII 0Ah

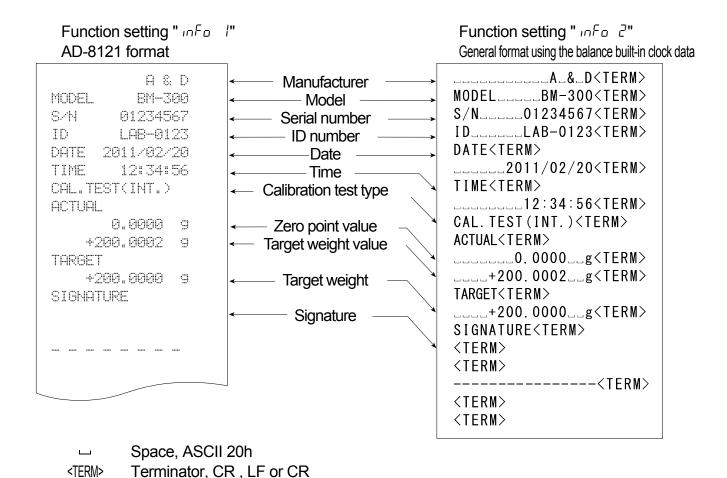

### Calibration report using an external weight

Note If the data memory function is used (except dfltfl (1)), the "Title block" and "End block" cannot be output.

#### Key operation

- 1 Press and hold the CAL key to display [FRL out] and release the key.
- 2 [FIL [] is displayed.
- 3 To update the calibration weight value, press the RANGE key and proceed to Step 4. To use the preset calibration weight value, proceed to Step 5.
- 4 Specify the calibration weight value using the following keys.


RANGE key.....To select the blinking digit.

RE-ZERO (+)key...To increase the value of the blinking digit.

MODE (-)key ······To decrease the value of the blinking digit.

PRINT key .....To store the new value.

- 5 Press the PRINT key. The zero point is measured and the weight value is displayed for a few seconds.
- 6 Place the displayed weight on the pan. Press the PRINT key to measure the weight and the weight value is displayed for a few seconds.
- 7  $[E \cap d]$  is displayed. Remove the weight from the pan.
- 8 If GLP output is used, [GLP] is displayed and the calibration report is output.
- 9 The balance returns to the weighing mode.

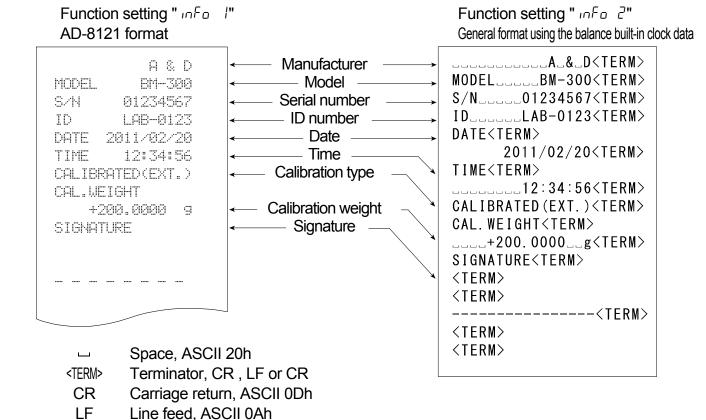

### Calibration test report using an external weight

- Note 

  Calibration test does not perform adjustment.
  - □ If the data memory function is used (except dfltfl ①), the "Title block" and "End block" cannot be output.

#### **Key operation**

- 1 Press and hold the CAL key to display [[ and release the key.
- 2 [FL [] is displayed.
- 3 To update the target weight value, press the RANGE key and proceed to Step 4. To use the preset target weight value, proceed to Step 5.
- 4 Specify the target weight value using the following keys.


RANGE key .....To select the blinking digit.

RE-ZERO (+)key ···· To increase the value of the blinking digit.

MODE (-)key ...... To decrease the value of the blinking digit.

PRINT key.....To store the new value.

- 5 Press the PRINT key. The zero point is measured and the weight value is displayed for a few seconds.
- 6 Place the displayed target weight on the pan. Press the PRINT key to measure the weight. The weight value is displayed for a few seconds.
- 7  $\boxed{\mathcal{E} \cap d}$  is displayed. Remove the weight from the pan.
- 8 If GLP output is used, [GLP] is displayed and the calibration test report is output.
- 9 The balance returns to the weighing mode.

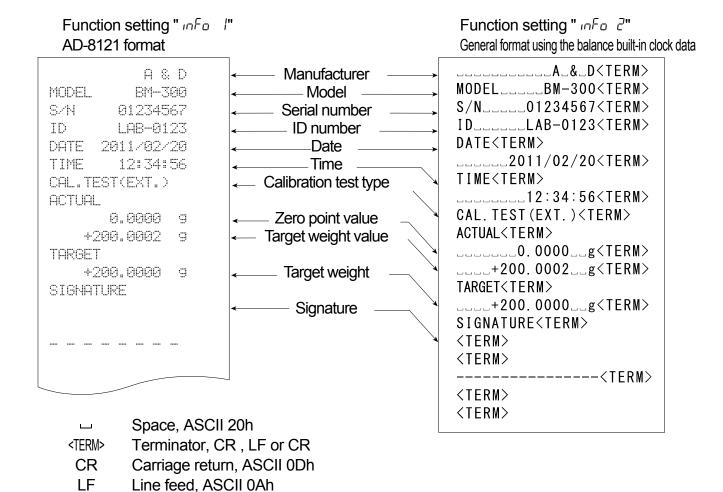


#### Title block and end block

When the weight values are recorded as GLP data, a "Title block" is inserted at the beginning and an "End block" is inserted at the end of a group of weight values in the GLP report.

Note If the data memory function is used (except dALA 0), the "Title block" and "End block" cannot be output.

#### **Key operation**

- 1 With the weighing data displayed, press and hold the PRINT key (for approx. 2 seconds) until 5£8r£ is displayed, and then release the key. The "Title block" is output.
- 2 The weighing data is output according to the parameter setting of the data output mode ( $P_r t$ ) of the function table.
- 3 Press and hold the PRINT key (for approx. 2 seconds) until release the key. The "End block" is output.

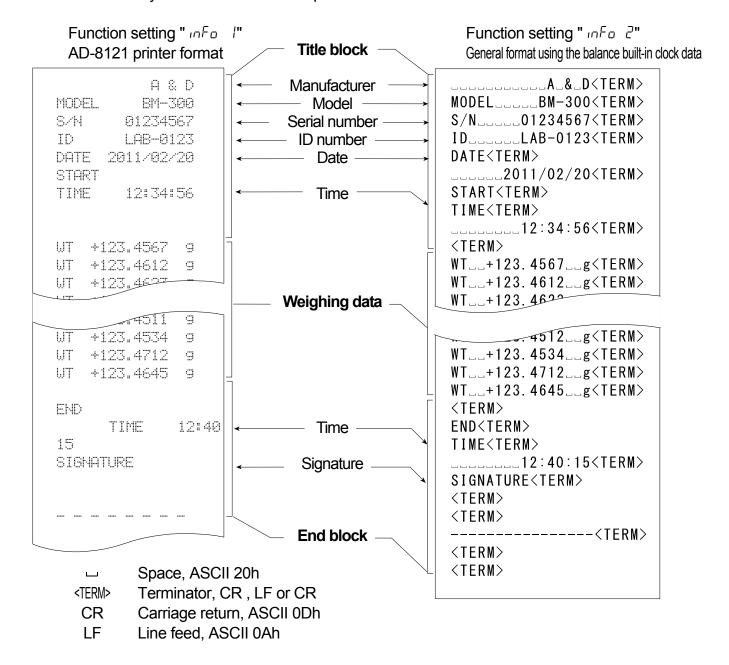

# 12. Data Memory

Data memory is a function to store weighing data and calibration data in memory. The data stored in memory are available for outputting at one time to a printer or personal computer.

Weighing data	Excluding the time and date	Up to 200 sets							
Including the time and date  Up to 100 sets									
•	ng the internal mass and an external weight comatic self calibration	Last 50 sets							
	Using the internal mass and an external weight	2401 00 0010							

# 12.1. Notes on Using Data Memory

- □ To use the memory function, set the "Data memory (dฅ೬ฅ)" parameter and the "Time/Date output (5-៤d)" parameter of the function table. Refer to "10. Function Table" for details on setting the data memory.
- □ For weighing data, the data contents to be stored and the storage capacity depend on the "Time/Date output (5-৮d)" parameter setting.
- If a different type of data exists in memory, "[Lr" blinks in the upper left of the display to prompt you to delete the data.
  For example: You want to store weighing data but calibration data or unit mass data remains in memory. "[Lr" blinks.
- If the stored data is corrupted, "Εςς" blinks in the upper left of the display.
   Under such a condition, before storing data, delete the data in memory as follows:

## Releasing "[Lr" or "Err"

- 1 Press and hold the PRINT key (for approx. 2 seconds) until Ltr no with "no" blinking is displayed, and then release the key.
- 2 Press the RE-ZERO key to display [[Lr [[g]]] with "[g]" blinking.
- 3 Press the PRINT key to delete all the data in memory.
- 4 The balance displays  $\lceil \mathcal{E} \cap d \rceil$  and returns to the weighing mode.

# 12.2. Data Memory for Weighing Data

#### **Features**

- □ The balance can store 200 sets of weighing data in memory (if the time and date are added, the balance can store 100 sets). Even if the AC adapter is removed, the data is maintained in the memory.
- It is not necessary to connect a printer or personal computer to the balance continually, because the balance stores the weighing data in memory.
- The data memory function does not occupy a printer or personal computer while weighing.
- □ The data in memory can be displayed on the balance for confirmation.
- The data in memory can be output at one time to a printer or personal computer.
   Comment: Select items for addition to the output data, such as the ID number, data number, time and date, in the function settings.

Upper left of the display

I.L.

### Storing the weighing data

Note If "[[r"] blinks in the upper left of the display, delete the data in memory.

- 1 Set the "Data memory (dRER)" parameter to ";" (To store the weighing data).
- 2 Specify whether or not to add the time and date using the "Time/Date output (5-6d)" parameter.
- 3 The storing mode depends on the "Data output mode (Prt)" parameter setting. Four modes are available to store data.

Key mode ......When the PRINT key is pressed and the displayed value is stable, the balance stores the weighing data.

Auto print mode A ......When the displayed value is stable and the conditions of "Auto print polarity", "Auto print difference" and zero point (reference value) are met,

the balance stores the weighing data.

Auto print mode B ......When the displayed value is stable and the conditions of "Auto print polarity", "Auto print difference" and last stable data (reference value)

are met, the balance stores the weighing data.

Interval mode......Weighing data is stored at an interval specified in "Interval time ( יחל )".

Press the PRINT key to start and stop this mode.

## **Enabling the data memory function**

1 Press and hold the RANGE key (for approx. 2 seconds) until b85Fnc is displayed, and then release the key.

2 Press the RANGE key several times to display doub.


3 Press the PRINT key.

4 Press the RANGE key several times to display dRLR.


5 Press the RE-ZERO key to display dRER !

6 Press the PRINT key to store the setting.

7 Press the CAL key to return to the weighing mode.

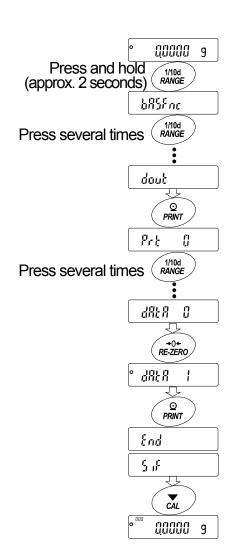

### Display symbols for data memory

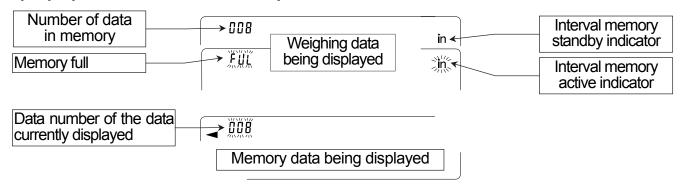

#### **Notes**

- When weighing data is being stored in memory, the data can be output to a personal computer using the RS-232C interface at the same time.
- □ "FUL" indicates that memory is full or the memory capacity has been reached. More data cannot be stored unless the memory data is deleted.
- Automatic self calibration cannot be used while the interval memory mode is active.
- The following commands cannot be used during data storage.

## Setting the function table

Parameter settings for each output mode are as follows:

Item	Data output	Auto print polarity,	Data memory	Interval time
Mode	mode	difference	function	interval time
Key mode	Prt O	Not used	JAFA I	
Auto print mode A	Prt I	AP-A O to 2	dRER I	Not used
Auto print mode B	Prt 2	ЯР-Ь О <b>to</b> 2	dRER I	
Interval mode	Prt 3	Not used	dRER I	ınĿ 0 <b>to</b> 8

Additional parameter settings, as follows:

Data number	No	d-no off
Data Humber	Yes	d-no on
ID number	No	5- id off
ID Humber	Yes	5- id on

Time and date	No	5-F9 0	200 sets
	Time only	5-Ed 1	
	Date only	5-Ed 2	100 sets
	Both	5-Ed 3	

### Recalling the memory data

Confirm that the "Data memory (dRLR)" parameter is set to " /".

- 1 Press and hold the PRINT key (for approx. 2 seconds) until FERLL is displayed, and then release the key. The type of data appears in the upper left of the display as shown to the right.
- 2 Press the PRINT key to enter the memory recall mode. Recall the data in memory using the following keys. RE-ZERO key ······To proceed to the next data set. MODE key ·······To go back to the previous data set.

PRINT key······To transmit the current data using the RS-232C interface.

With RANGE held down, press the CAL key

To delete the current data.

CAL key.....To exit the memory recall mode.

3 Press the CAL key to return to the weighing mode.

Indicators

-d
Weighing data
without time and date


Weighing data with time and date

## Transmitting all memory data at one time

Confirm that the "Serial interface (5 ,F)" parameters are set properly. Refer to "10. Function Table" and "16.2. Connecting to Peripheral Equipment".


- 1 Press and hold the PRINT key (for approx. 2 seconds) until FEERLL is displayed, and then release the key.
- 2 Press the RANGE key to display out
- 3 Press the PRINT key to display aut no with "no" blinking.
- 4 Press the RE-ZERO key to display συξ μως with "Γω" blinking.
- 5 Press the PRINT key to transmit all data using the RS-232C interface.
- 6 The balance displays [LLERr] when all data is transmitted.

Press the CAL key to return to the weighing mode.

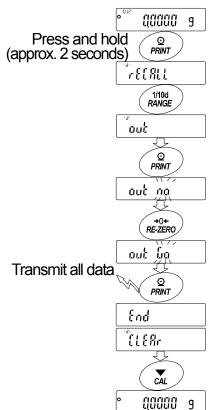

### Deleting all memory data at one time

- 1 Press and hold the PRINT key (for approx. 2 seconds) until recent is displayed, and then release the key.
- 2 Press the RANGE key several times to display
- 3 Press the PRINT key to display [Lr no with "no" blinking.
- 4 Press the RE-ZERO key to display [[Lr [μ]]] with "[μ] blinking.
- 5 Press the PRINT key to delete all data.
- 6 The balance displays  $\boxed{\textit{End}}$  and returns to the weighing mode.

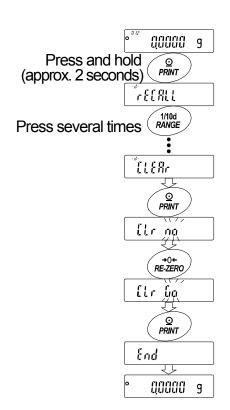

# 12.3. Data Memory for Calibration and Calibration Test Data

#### **Features**

- Calibration data (when and how it is performed) and calibration test data can be stored in memory.
- All the data in memory is available to be output at one time to a printer or personal computer.
- □ Up to 50 data sets of the latest calibration or calibration test can be stored. When the memory capacity has been reached, "FUL" is displayed in the upper left of the display.

Indicator FUL

## Storing the calibration and calibration test data

Note If "[Lr" appears blinking in the upper left of the display, delete the data in memory.

Store the calibration and calibration test data as follows:


- 1 Set the "Data memory (♂RER)" parameter to "?" (to store the calibration data).
- 2 Set the "GLP output ( InFa)" parameter to " !", "2" or "3".
- 3 With the settings above, each time calibration or calibration test is performed, the data is stored automatically.

# Transmitting the memory data Notes

- □ Confirm that the "Serial interface (5 ,F)" parameters are set properly.


  Refer to "10. Function Table" and "16.2. Connecting to Peripheral Equipment".
- □ Confirm that the "Data memory (dALA)" parameter is set to "2".
- 1 Press and hold the PRINT key (for approx. 2 seconds) until aut is displayed, and then release the key.
- 2 Press the PRINT key to display out no with "no" blinking.
- 3 Press the RE-ZERO key to display συλ μω with "Γο" blinking.
- 4 Press the PRINT key to transmit all memory data using the RS-232C interface.
- 5 The balance displays [LERr] when all memory data is output.

  Press the CAL key to return to the weighing mode.

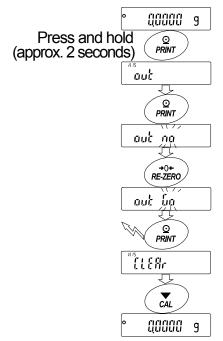

## Deleting data stored in memory

- 1 Press and hold the PRINT key (for approx. 2 seconds) until aut is displayed, and then release the key.
- 2 Press the RANGE key to display [LERr]
- 3 Press the PRINT key to display [Lr ng with "no" blinking.
- 4 Press the RE-ZERO key to display [Lr Lo] with "Lo" blinking.
- 5 Press the PRINT key to delete all data.
- 6 The balance displays *End* and returns to the weighing mode when all the data has been deleted.

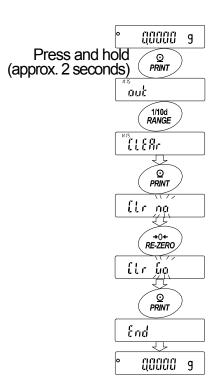

# 13. Underhook


The underhook can be used for magnetic materials or density measurement. The built-in underhook is revealed by removing the cap on the bottom of the balance. Use the underhook as shown below.

## Caution

- Do not apply excessive force to the underhook.
- When not in use, attach the cap to prevent dust from getting into the balance.
- Do not push the underhook upward
- When turning the balance over, parts and weighing sample in the weighing chamber will fall off. (Parts: the weighing pan, pan support, breeze break ring, fine range ring and dust plate)

Remove them first.

- When the cap is removed, a weighing error may occur, because of drafts entering into the internal portion of the balance. Arrange the room condition.
- 1 Remove the cap on the bottom of the balance.
- 2 Hang your weighing pan on the underhook. Place the balance on the solid table.

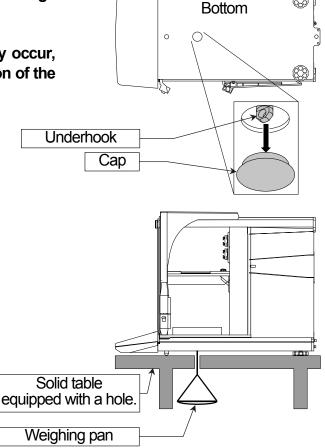

# 14. Density Measurement (Specific gravity)

The balance is equipped with a density mode. It calculates the density of a solid using the weight of a sample in air and the weight in liquid. We recommend the use of the AD-1653 Density Determination Kit.

#### **Notes**

- The density mode was not selected for use when the balance was shipped from the factory.
   To use this mode, change the function table settings to activate the density mode.
   Refer to "5.2. Storing Active Units".
- When the density mode is selected, the response adjustment function cannot be used.
- □ The readability is 0.0001 g while using the density mode.

## Formula to obtain the density

The density can be obtained by the following formula.

 $\rho = \frac{A}{A - B} \times \rho_0$  Where  $\rho : Density of a sample A : Weight value of a sample in air B : Weight value of a sample in liquid$ 

## Prior to measurement: Changing the function table

Prior to measurement, change the function table as follows:

1 Selecting the density mode.

The density mode is not available in the factory setting. To use the mode, select it (unit of Unit 15).) in the function table. Refer to "5.2. Storing Active Units". Then, the density mode can be selected as one of the weighing units by pressing the MODE key.

2 Selecting a way to set the density of a liquid.
Select the liquid density input method from the function table below.

Note The following function table is displayed only when the unit of  $[U \cap I \vdash IS]$  is active. The table of the density function  $(dS \vdash F \cap C)$  is displayed after  $S \vdash F$ .

Class	Item and Parameter		Description
dS Fnc	Ld in		Water temperature
Density function	Liquid density input	1	Liquid density

Factory setting.

## Entering the density of a liquid

Two ways to set the density of a liquid are available in the function table, "Liquid density input (Ld In): by entering the water temperature or by entering the density directly.

1 Press the MODE key as necessary to select the density mode. Press and hold the MODE key (for approx. 2 seconds) to select to "water temperature" or "liquid density". In the density mode: The unit is "g (gram)". The processing indicator (◄) blinks. Response adjustment is off. If you cancel the density mode, press the MODE key.



00000 9

ŀ

Press and hold

(approx. 2 seconds)

## Entering the water temperature (Ld in 11)

2 Press and hold the MODE key (for approx. 2 seconds). The water temperature currently set (unit: °C, factory setting : 25°C) is displayed. Use the following keys to change the value.

RE-ZERO (+)key .... To increase the temperature by one degree. (0°C is displayed after 99°C)

MODE (-)key ......... To decrease the temperature by one degree. (99°C is displayed after 0°C)

PRINT key......To store the new water temperature, display End and return to the density mode. Proceed to Step 1.

CAL key.....To cancel the change and return to the density mode. Proceed to Step 1.

The relation between the water temperature and density

°C	+0	+1	+2	+3	+4	+5	+6	+7	+8	+9
0	0.99984	0.99990	0.99994	0.99996	0.99997	0.99996	0.99994	0.99990	0.99985	0.99978
10	0.99970	0.99961	0.99949	0.99938	0.99924	0.99910	0.99894	0.99877	0.99860	0.99841
20	0.99820	0.99799	0.99777	0.99754	0.99730	0.99704	0.99678	0.99651	0.99623	0.99594
30	0.99565	0.99534	0.99503	0.99470	0.99437	0.99403	0.99368	0.99333	0.99297	0.99259
40	0.99222	0.99183	0.99144	0.99104	0.99063	0.99021	0.98979	0.98936	0.98893	0.98849
										a/om <sup>3</sup>

g/cm

00000 9

## Entering the density directly (Ld in 1)

Press and hold the MODE key (for approx. 2 seconds). The density currently set (unit: g / cm<sup>3</sup>, factory setting: 1.0000g / cm<sup>3</sup>) is displayed.

Use the following keys to change the value.

The range to set the density is  $0.0000g / cm^3$  to  $1.9999g / cm^3$ .

RE-ZERO (+)key..... To increase the value of the digit selected.

MODE (-)key ......To decrease the value of the digit selected.

RANGE key.....To select the digit to change the value.

PRINT key......To store the change, display End and return to the density mode. (Step 1).

CAL key ......To cancel the change and return to the density mode. (Step 1)

Press and hold MODE (approx. 2 seconds) a jjaaaa

## Measuring the density (specific gravity)

Note If the liquid temperature or the type of liquid is changed during measurement, input the value of the liquid density again as necessary. The density display is four decimal places. The readability cannot be changed with the RANGE key.

The density (or specific gravity) is displayed after "Weighing in air" and "Weighing in liquid". The procedure of each measurement is as follows:


- 1 Enter the density mode. (g lights. blinks.) Place nothing on both pans and press the RE-ZERO key to display zero.
- 2 Place the sample on the pan in air (g lights. ■ blinks.) and wait for the weight value to be stable. To store or output the weight value, press the PRINT key. Press the RANGE key to confirm the weight value in air and proceed to the next step.

Note If negative value or  $\boxed{\mathcal{E}}$  (out of range) is displayed, the  $\boxed{\mathsf{RANGE}}$  key is inactive.

Note If  $\mathcal{E}$  (out of range) is displayed, the RANGE key is inactive.

- 4 To store or output the density value, press the PRINT key. (Unit: №. g turns off. ✓ lights.)

  Press the RANGE key to measure another sample and proceed to Step 2.
- 5 If the liquid temperature or the type of liquid is changed during measurement, input the value of the liquid density again as necessary.
- 6 Press the MODE key to proceed to other modes.

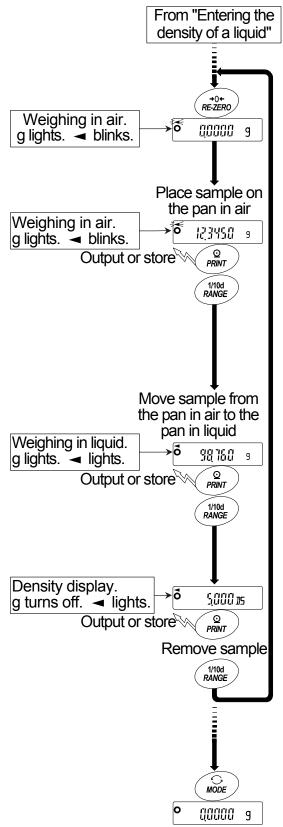

# 15. Password Function

The password function can be used to restrict balance use or functions. It is useful for preventing the time and date setting from being falsified and preventing users from changing other function settings.

Password: Four digits (4 x 4 x 4 x 4 = 256 outcomes) using the MODE, RANGE, PRINT and RE-ZERO keys.

At the factory setting, the password function has been disabled.

Settings related to the password function are performed in the function table.

Note The password function varies depending on the balance software version.

# 15.1. Balance Software Version 1.50 to 1.64

Two settings are available in "Password function (LocFnc)" of the function table.

PRSS 0	No password function.
PRSS	Entering a password is required before starting weighing.

## PR55 [] (No password function)

No password is required. Anybody can perform weighing, use all the balance functions and make changes to them.

## **PR55** (Entering a password is required before starting weighing.)

The administrator (Ranin) sets a unique password to restrict balance users. Only one password by the administrator can be set.

The correct password must be entered for the balance to enter the weighing mode.

# 15.2. Balance Software Version 1.660 or later

Three settings are available in "Password function (LocFnc)" of the function table.

PRSS [		No password function.
PRSS 1		Entering a password is required before starting weighing.
PRSS 2	•	Logging in by the administrator password is required before making changes.

## **PRSS** (No password function)

No password is required. Anybody can perform weighing, use all the balance functions and make changes to them.

## **PR55** (Entering a password is required before starting weighing.)

The administrator ( $Rd\tilde{n}IN$ ) sets unique passwords to restrict balance users. Pressing the ON:OFF key before weighing will prompt you to enter a password.

The correct password must be entered for the balance to enter the weighing mode.

Two login levels are available: administrator (#dn]N) and user (USER III to III).

Administrator (#dnIN)	All the functions and settings can be used.		
Autilistator (nanin)	A password can be set for each of 10 users.		
User (USER ()) to (())	Initialization and changes to the settings are restricted, including the clock setting.		
No password The balance cannot be used.			

## **PR55** 2 (Logging in by the administrator password is required before making changes.)

Anybody can perform weighing, but initialization and changes to the settings are restricted, including the clock setting. Even if the ON:OFF key is pressed to start weighing, a password will not be required.

Two login levels are available: administrator (AdnIN) and guest ([[][[5]]).

Administrator (RdnIN)	All the functions and settings can be used.				
Guest ([]][5])	Initialization and changes to the settings are restricted,				
*No password	including the clock setting.				

When weighing is started by pressing the ON:OFF key while holding down the CAL key in the standby mode, the balance will prompt you to enter the administrator password.

## Items restricted depending on the login level

	Weighing				
Login level	Entering password before weighing	Calibration	Changes to the settings *1		
Administrator (RdnIN)	Required	Available	Available		
User (USER ()) to (())	Required	Available or not available	Not available		
Guest ( <b>[]][</b> [5])	Not required	* <b>2</b>			

<sup>\*1:</sup> Response adjustment, repeatability test, function switch and initialization, function settings (clock setting)

<sup>\*2:</sup> Available under ordinary circumstances, but not available to users and guests when the administrator sets the function switch to "inhibit".


# 15.3. Using the Password Function


The password function can be switched among "Not used ( $\mathcal{I}$ )", "Used ( $\mathcal{I}$ )" and "Used ( $\mathcal{I}$ )" in "Password function ( $L a \in \mathcal{F} \cap \mathcal{I}$ )" of the function table.

- 1 In the weighing mode, press and hold the RANGE key (for approx. 2 seconds) until base is displayed, and then release the key.
- 2 Press the RANGE key several times to display LocFnc
- 4 Press the RE-ZERO key to switch the numbers. Display " *l*" or "∠".
- 5 Press the PRINT key to display 5ur E: 9E5 Na. with "No" blinking (selected).
- 6 Press the RE-ZERO key to switch between "YES" and "No".
- 7 Display <u>โรนายี เนีย์วิห</u> with "<u>ป</u>ียีรี" blinking (selected).
- 8 Press the PRINT key to store the new setting. The password function can now be used.
- 9 Lac na is displayed. To store or change the password, proceed to Step 3 in "15.6. Storing or Changing the Password". Not to store or change the password, press the CAL key to return to the weighing mode.


Note When the balance software version is 1.50 to 1.64, only "Not used (I)" or "Used (I)" is available.

78

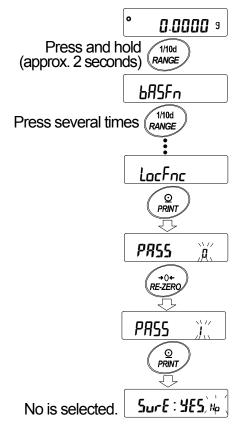

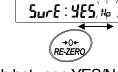

Switch between YES/No

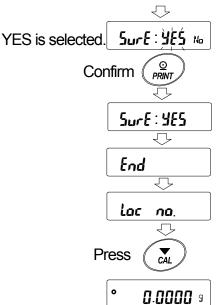

# 15.4. Entering the Password Before Weighing

## 15.4.1. PR55 (Entering a password is required before starting weighing)

1 In the standby mode, press the ON:OFF key.

2 After PRSS NII, ---- PRSS is displayed to prompt you to enter a password.

3 Enter a four-digit password using the following keys. The display will turn off automatically after ten minutes of non-operation.

MODE key......To enter  $\overline{n}$ RANGE key.....To enter  $\overline{n}$ PRINT key.....To enter  $\overline{n}$ RE-ZERO key.....To enter  $\overline{n}$ 


CAL key .....To go back one character

4 When the password is correct, the balance displays the login level, turns all the display segments on, and then enters the weighing mode.

When the administrator password is entered, you are logged in as an administrator.

The default password set at the factory is (press the RE-ZERO) key four times), for administrator level.

When the password is incorrect, the balance displays FRIL, sounds the buzzer three times and turns the display off.

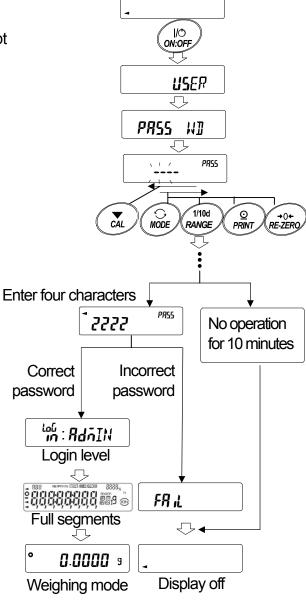

Note When the balance software version is 1.50 to 1.64, the login level is not displayed when logged in.

When the balance software version is 1.660 or later, enter the administrator password to log in as an administrator.

# 15.4.2. PR55 2 (Logging in by the administrator password is required before making changes)

# Logging in as an administrator (AdnIN)

- 1 In the standby mode, while holding down the CAL key, press the ON:OFF key.
- 2 Enter a four-digit password using the following keys. The display will turn off automatically after ten minutes of non-operation.

MODE key......To enter  $\bar{n}$ RANGE key......To enter  $\bar{n}$ PRINT key......To enter  $\bar{p}$ RE-ZERO key......To enter  $\bar{c}$ CAL key......To go back one character


3 When the password is correct, the balance displays the login level, turns all the display segments on, and then enters the weighing mode.


The default password set at the factory is [2222] (press the RE-ZERO] key four times), for administrator level.

When the password is incorrect, the balance displays FRIL, sounds the buzzer three times and turns the display off.

# Logging in as a guest ([][][5])


- 1 In the standby mode, press the ON:OFF key.
- 2 After LUE, the balance enters the weighing mode.

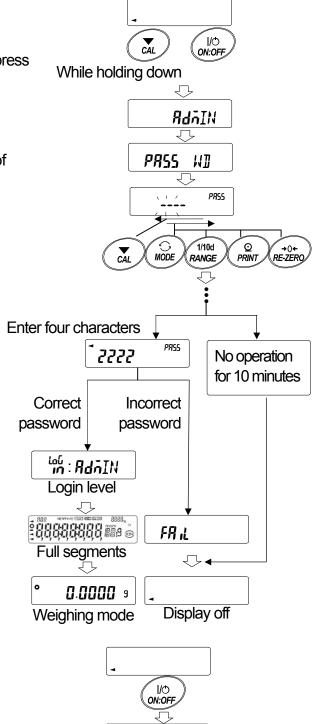

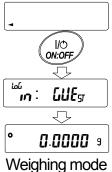

# 15.5. Logging Out

1 To log out, press the ON:OFF key to turn the display off. With the function PR55 / set, the password will need to be entered again to enter the weighing mode from the standby mode.

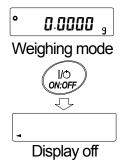


# 15.6. Storing or Changing the Password

The password can be changed in "Password (Lac na)" of the function table.

- 1 In the weighing mode, press and hold the RANGE key (for approx. 2 seconds) until basence is displayed, and then release the key.
- 2 Press the RANGE key several times to display Lac na.
- 3 Press the PRINT key to display the login level (Adn IN).
- 4 Press the RANGE key to display the login level to be changed (RdnIN/USER UI to UI).

  The stabilization indicator appears at a login level where a password has been stored. (The password can be changed.)
- 5 To change the password, press the PRINT key. For details, refer to "15.7. Changing the Password".

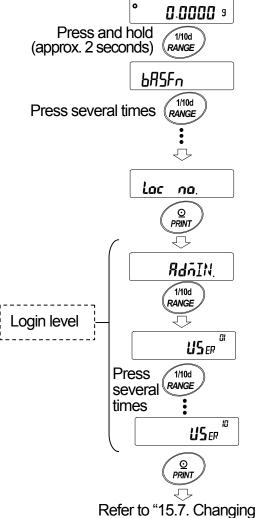

Refer to "15.7. Changing the Password".

#### **Notes**

- To log out, press the ON:OFF key to turn the display off.
- □ With the function PRSS 2 set, the administrator password is required to log in as an administrator. No password needs to be stored for USER UH to U.
- □ When the balance software version is 1.50 to 1.64, only one password can be stored. While the balance displays 

  L□c □□□, press the PRINT key to display the current password.

# 15.7. Changing the Password


- 1 Refer to "15.6. Storing or Changing the Password" to display the login level to be changed.
- 2 Press the PRINT key to display the current password. The default password set at the factory is [2222] (press the RE-ZERO key four times), for administrator level.
- 3 Enter a new four-digit password using the following keys. The display will turn off automatically after ten minutes of non-operation.

MODE key......To enter no ente

- 4 The new password is displayed.
- 5 <u>SurE : YES No.</u> is displayed with "No" blinking (selected).

  Press the <u>CAL</u> key to return to the fourth character entry position.
- 6 Press the RE-ZERO key to display Sur E: YES No with "YES" blinking (selected).
- 7 While "YES" is blinking, press the PRINT key to store the new password.
- When setting is complete, End and then the next login level is displayed.
   To set a password for the next level, follow the procedure starting from Step 2.

To finish the operation, press the CAL key twice to return to the weighing mode.

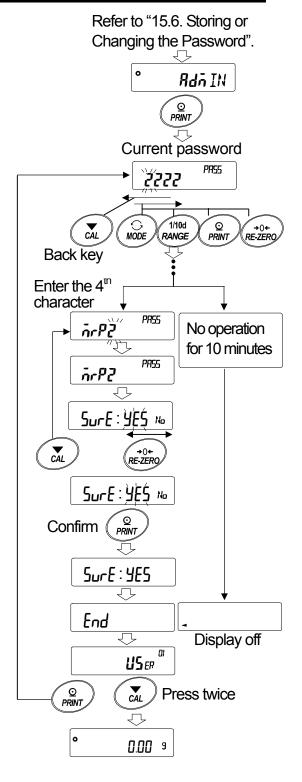

## **Notes**


- If you have forgotten the password, the balance cannot be used. Make a record of the password and keep it where you can refer to it if needed.
- □ The same password as the administrator password cannot be used for USER II to II.

# 15.8. Deleting the Password (USER CI)

- 1 Refer to "15.7. Changing the Password" to select the user whose password to be deleted and display the password.
- 2 Press and hold the CAL key (for approx. 2 seconds) to display \(\begin{aligned} \tilde{\mathcal{E}} \tilde{\mathcal{E}} \mathcal{F} \\ \tilde{\mathcal{E}} \end{aligned}\) (blinking).
- 3 Press the PRINT key to display [[LERr] ]
- 4 Press the RE-ZERO key to switch between " $G_{\alpha}$ " and " $N_{\alpha}$ ".
- 5 While [[LERr] is displayed, press the PRINT] key to delete the password.

  The balance displays [End] and then displays the login level.

Refer to "15.7. Changing the Password".

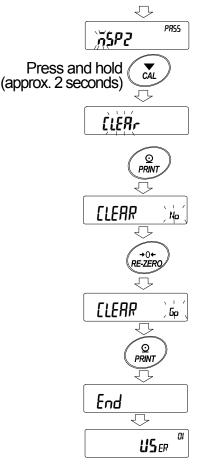

Note The administrator password cannot be deleted, but it can be changed.

Refer to "15.6. Storing or Changing the Password" and "15.7. Changing the Password" to change it.

# 15.9. Forgotten Password

If you have forgotten the password, the balance cannot be used. Contact your local A&D dealer to undo the password.

# 16. Standard Input and Output Interface

# 16.1. RS-232C Interface

The balance is a DCE device. Connect the balance to a personal computer (DTE) using a straight through cable.

Transmission system : EIA RS-232C

Transmission form : Asynchronous, bi-directional, half duplex

Transmission rate : 10 times/second or 5 times/second (same as data refresh rate)


Data format : Baud rate : 600, 1200, 2400, 4800, 9600, 19200 bps

Data bits : 7 or 8 bits

Parity: Even, Odd (Data bits 7 bits)


None (Data bits 8 bits)

Stop bit : 1 bit
Code : ASCII RS-232C

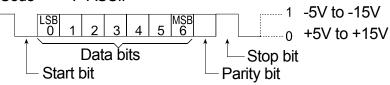


D-Sub 9 pin assignments

Pin No.	Signal name	Direction	Description
1	_	_	No connection
2	RXD	Output	Transmit data
3	TXD	Input	Receive data
4	_	_	No connection
5	SG	_	Signal ground
6	DSR	Output	Data set ready
7	RTS	Input	Request to send
8	CTS	Output	Clear to send
9	_	_	No connection

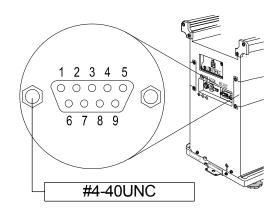

All signal names mean name for the DTE side.

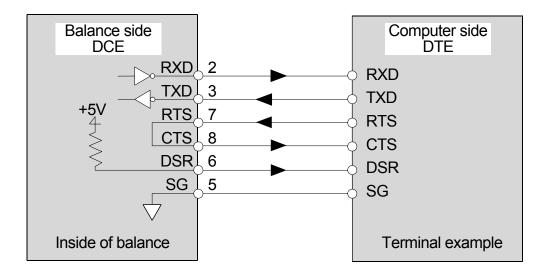

# 16.2. Connecting to Peripheral Equipment

# 16.2.1. Connecting to the AD-8121B Printer

When connecting the AD-8121B printer to the balance for printing weighing data, configure the printer and the balance as follows

Class	Item and Parameter	Factory settings	AD-8121B MODE 1	AD-8121B MODE 2	AD-8121B MODE 3	
	Prt Data output mode	0	0, 1, 2	3	0, 1, 2	
	RP-P Auto print polarity	0	Ш4	NI (C	11.4	
	RP-ь Auto print difference	1	#1	No setting	#1	
dout	5- id ID number output	oFF	oFF	oFF	oFF, on	
Data output	d-no Data number output	oFF	oFF	oFF	oFF, on	
	5-とd Time/Date output	0	0	0	0, 1, 2, 3	
	PUSE Data output pause	oFF	oFF	oFF	oFF, on #2	
	RE-F Auto feed	oFF	oFF	oFF	oFF, on	
	երը Baud rate	2	2	2	2	
5 ,F Serial interface	ьեРг Data bit, parity bit	0	0	0	0	
	ErlF Terminator	0	0	0	0	
	ੁ ੁੁੁੁੂੁੁੂੂ Data format	0	0	0	1	

<sup>#1</sup> Set the parameters when auto print mode A or B ( $Prt \mid$  or 2) is selected.

#### **Notes**

Refer to "11.2. GLP Report" for print samples.

Settings of AD-8121B

MODE	AD-8121B DIP switch	Description
MODE 1	MODE \$ \$ \$	Print at receiving data. Standard mode, statistical calculation mode
MODE 2	MODE	Print by DATA key operation or built-in timer. Standard mode, interval mode, chart mode
MODE 3	MODE TO STATE OF THE PROPERTY	Print at receiving data. Dump print mode

DIP switch No.3: Handling unstable data

ON To print
OFF Not to print

DIP switch No.4: Data input specifications (Interface selection)

ON Current loop OFF RS-232C





<sup>#2</sup> Set an when multiple lines are printed. Example: When adding ID number, set an.

# 16.2.2. Connecting to the AD-8127 Printer

When connecting the AD-8127 printer to the balance for printing weighing data, configure the printer and the balance as follows according to the usage examples.

Function settings of the printer

	AD-8127
Usage example	Multi printer
	Printing mode settings
When printing the balance weighing data	
with the balance's PRINT key or its	EXT.KEY
"Auto print" mode.	
When printing the balance weighing data	MANUAL
with the printer's "Printing" key or its	AUTO
"Interval printing" mode.	TIMER
When printing charts with the printer.	CHART
When printing the balance's statistical	
calculation results.	DUMP
When printing the balance's GLP report.	

<sup>□</sup> Refer to the instruction manual of the AD-8127 printer for how to change the function settings of the AD-8127 printer.

Function settings of the balance

Usage example	Balance ア <sub>r</sub> Ł Data output mode	Balance Ł YPE Data format
When printing the balance weighing data with the balance's PRINT key or its "Auto print" mode.	0,1,2	0
When printing the balance weighing data with the printer's "Printing" key or its "Interval printing" mode. When printing charts with the printer.	3	0
When printing the balance's statistical calculation results. When printing the balance's GLP report.	0, 1, 2	1

<sup>□</sup> Refer to "10. Function Table" for how to change the function settings of the balance.

# 16.2.3. Connecting to a Computer Using the RS-232C Interface

## **Features**

The balance is of the DCE type (Data Communication Equipment), which can be connected to a personal computer using the RS-232C interface. Before connection, read the personal computer manual thoroughly. Use a standard DCE cable for connection (cable type: straight-through).

## Using Windows Communication Tools Software (WinCT)

When Windows is used as an operating system in a personal computer, the WinCT software can be used to transmit the weighing data to the personal computer. The current version of the WinCT can be downloaded from the A&D website. Confirm the windows version supported by this software when downloading. Refer to the WinCT instruction manual in the A&D website. The WinCT software has three communication methods: "RsCom", "RsKey" and "RsWeight".

## RsCom

- RsCom can transmit commands to control the balance.
- RsCom can make bi-directional communication between the balance and a personal computer using the RS-232C interface.
- RsCom can display or store the data using a text file format. RsCom can also print the data using a printer connected to the personal computer.
- □ When several ports of a personal computer have balances connected, the personal computer can communicate with each balance simultaneously.
- RsCom can share a personal computer with other application software.
- RsCom can receive the balance GLP report.

## RsKey

- RsKey can transmit the weighing data output from the balance directly to other application software such as Microsoft Excel.
- RsKey can be used with most application software.
- RsKey can receive the balance GLP report.

## RsWeight

- Can retrieve the weight data from the balance and display the data as a graph in real time.
- Can calculate and display the maximum, minimum, average, standard deviation and coefficient of variation.

# Using the WinCT software, the balance can do the following:

Analyzing the weighing data and the statistics input by "RsKey"

The weighing data can be input directly into an Excel worksheet. Then, Excel can analyze the data to obtain total, average, standard deviation, maximum and minimum values, and display them in a graph.

Controlling the balance using commands from a personal computer

By using "RsCom", the personal computer sends commands such as "re-zero" or "send weighing data" to the balance and controls the balance.

Printing the balance GLP report using your printer

The balance GLP report can be printed using a printer connected to the personal computer.

Receiving weighing data at a certain interval

The weighing data can be received at a certain interval and data characteristic with elapsed time can be obtained.

Using the balance memory function

The data can be stored in the balance's memory. Of the data stored, the weighing data and calibration data can be transmitted to a personal computer at one time.

Using a personal computer as an external indicator

With the "RsKey" test mode function, a personal computer can be used as an external weight indicator for the balance. (To do this, set the balance data output mode to stream mode.)

## 16.2.4. Connecting to a Computer Using the USB Interface

## **Features**

- The weighing data can be transmitted to a personal computer when connecting the balance and personal computer via the USB interface. The applicable operating system is Windows 7 or later, as of December 2018.
- The USB interface uses standard device driver software and does not require installation of special driver software. By simply connecting, weighing data can be transmitted.
- □ The data can be transmitted to most application software (ex.: EXCEL, WORD, Notepad).

#### **Notes**

- The balance can transmit the weighing data only without the header and unit.
- The USB interface of the balance cannot receive the command that controls the balance. When using the command, connect to the RS-232C interface of the balance. When connecting the USB interface of a personal computer to the RS-232C interface of the balance, use the USB converter (AX-USB-9P).
- Do not use the screensaver or suspend mode of a personal computer.
- Even if stream mode is used, output will be performed approximately 2 times/second due to key input operation.
- □ To add the ID number, data number, time and date to the weighing data or to output the GLP report, set the "Data output pause (PUSE)" parameter of "Data output (daut)" to "an" (Pause 1.6 seconds). The time and the date are separated by a dot when output.
- □ When the "Data format (⊱ℲℙΕ)" parameter is set to "5" (CSV format), the ID number, data number, time and date cannot be added to the weighing data.

## Using the USB interface

- Start the Windows of the personal computer.Connect the AC adapter to the balance and press the ON:OFF key to turn the display on.
- 2 Connect the personal computer to the balance using the USB cable.

Note Do not use stream mode of data output mode (doub: Prt 3).

- 3 At the first connection only, the USB driver software is automatically installed.
- 4 Open the application (ex.: EXCEL) that is to receive the weighing data. Move the cursor onto the position to input the weighing data.

Case 1: Transmitting the weighing data with key operation

- 5 Press the PRINT key to input the weighing data at the position of the cursor.
- 6 To finish the operation, remove the USB cable.

## **Examples**

□ Case 2: Transmitting the stable weighing data  Data output mode of the balance: Auto print mode A and B,  When a sample is placed and the stabilization indicator o  displayed, the weighing data is transmitted. This mode is useful	t:Prt O
When a sample is placed and the stabilization indicator o is	
· · ·	Ł lor∂
displayed, the weighing data is transmitted. This mode is useful	
when repeating the weighing.	

□ Case 3: Transmitting the weighing data periodically

Data output mode of the balance: Interval mode,

The weighing data is transmitted periodically.

The interval time: 2, 5, 10, 30 seconds, 1, 2, 5, 10 minute(s).

# 16.3. Commands

# 16.3.1. Command List

Note Before transmitting a command, add the terminator that is specified in the "Terminator  $(f \cap LF)$ " parameter of "Serial interface  $(f \cap F)$ " in the function table.

Commands to query weighing data			
С	Cancels the S or SIR command.		
Q	Requests the weighing data immediately.		
S	Requests the weighing data when stabilized.		
SI	Requests the weighing data immediately.		
SIR	Requests the weighing data continuously.		

Commands to control the balance				
CAL	Same as the CAL key.	Calibration using the internal mass.		
OFF	Turns the display off.			
ON	Turns the display on.			
P	Same as the ON:OFF key.	Turning the display on or off.		
PRT	Same as the PRINT key.			
R	Same as the RE-ZERO key.	Zero display. #1		
RNG	Same as the RANGE key.	Changing the readability.		
TR	Tares the balance.	Canceling the container's weight. #1		
TST	Performs calibration test.	Calibration test using the internal mass.		
U	Same as the MODE key.	Changing the unit.		

Commands to request data		
?ID	Requests the identification number.	
?PT	Outputs the tare value.	#1
?SN	Requests the serial number of the balance.	
?TN	Requests the model name of the balance.	

Command to set data		
PT:***.*** பப g	Sets the tare value.	#1

- #1 
  Command R assumes the point as zero and sets the display to zero.
  - Command TR cancels the tare weight when the displayed value is greater than zero.
     The canceled tare value can be read using the command ?PT.
  - - An example of this command sets a negative target value and fills with the sample until the display becomes zero. For details, refer to page 94.
    - Step 1 Place a container.
    - Step 2 Set the display to zero using the R command.
    - Step 3 Set a target value using the PT: command.

## 16.3.2. Acknowledge Code and Error Codes

When the "AK, Error code  $(E \cap E d)$ " parameter of "Serial interface function  $(5 \cap F)$ " is set to "an", the balance outputs <AK> code or error code to each command as follows:

<a>AK> (06h)</a> Acknowledge in ASCII code.

 $\ \square$  When the balance receives a command to request data and cannot process it, the balance transmits an error code ( $\ \square \ \square \ \square$ ).

When the balance receives a command to request data and can process it, the balance outputs the data.

When the balance receives a command to control the balance and cannot process it, the balance transmits an error code (EC, Exx).

When the balance receives a command to control the balance and can process it, the balance transmits the acknowledge code.

<a>AK> (06h)</a> Acknowledge in ASCII code.

Among commands to control the balance, the following transmit the acknowledge code both when the balance receives the command and when the balance has accomplished the command. If the command cannot be processed properly, the balance transmits an error code (EC, Exx). This error can be resolved using the CAL command. xx is error code number.

CAL command (Calibration command using internal mass)

ON command (Display ON command)

P command (Display ON/OFF command)

R command (RE-ZERO command)

TR command (Tare command)

TST command (Calibration test command)

When a communication error has occurred due to external noise, or a parity error has occurred due to transmission error, the balance transmits an error code (EC, Exx). In this case, send the command again. xx is error code number.


## 16.3.3. Settings Related to RS-232C

Concerning the RS-232C, the balance has two functions: "Data output (daut)" and "Serial interface  $(5 \ F)$ ". Set each function as necessary.


# 16.3.4. Command Examples

This example uses "Er[d an" of "5 F" so that the <AK> (06h) code is output.


# ON command (turning the display on)

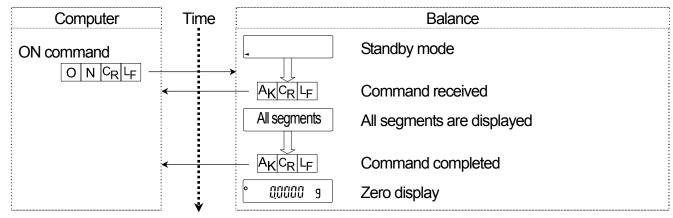

## R command (Re-zeroing the display)

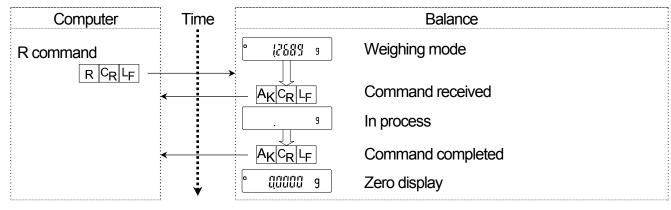


## CAL command (Calibration with internal mass)

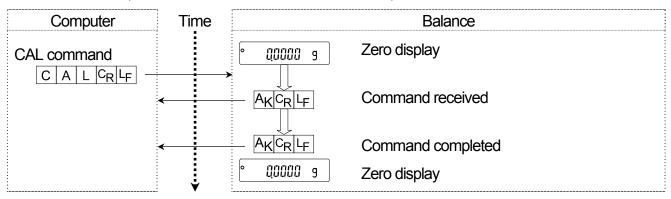


## Error code

This example is of an error using the R command. "ErEd an" is used. The balance transmits an error code when the received command cannot be achieved.

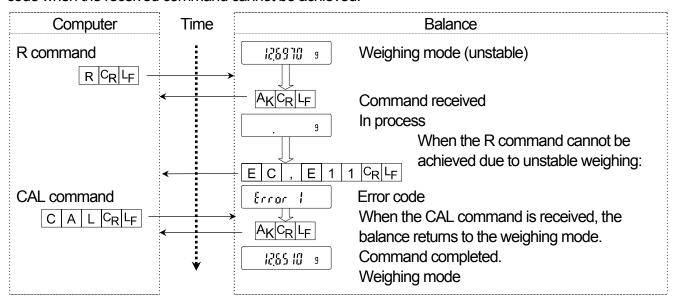


## Weighing with a tare

This example uses " $\mathcal{E}_{\mathcal{F}}[d \circ \mathcal{F}]$ " of "5  $\mathcal{F}$ " so that the <AK> (06h) code is output.

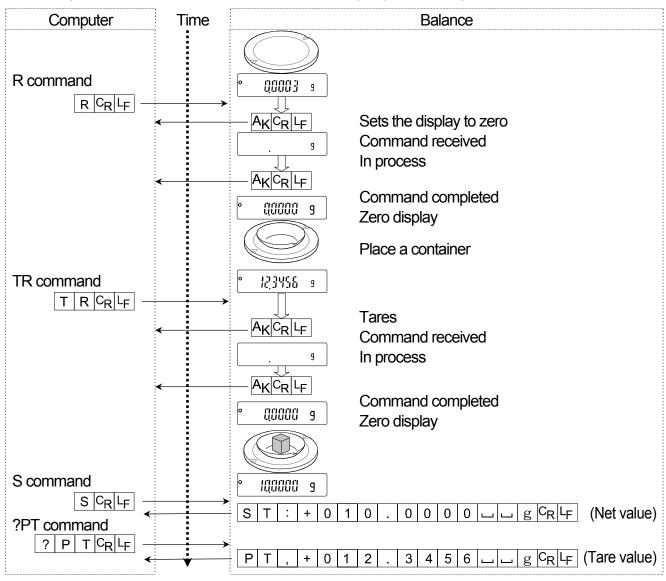

# Setting a negative target value and filling with a sample until the display becomes zero

This example uses "Er[d an" of "5 F" so that the <AK> (06h) code is output.

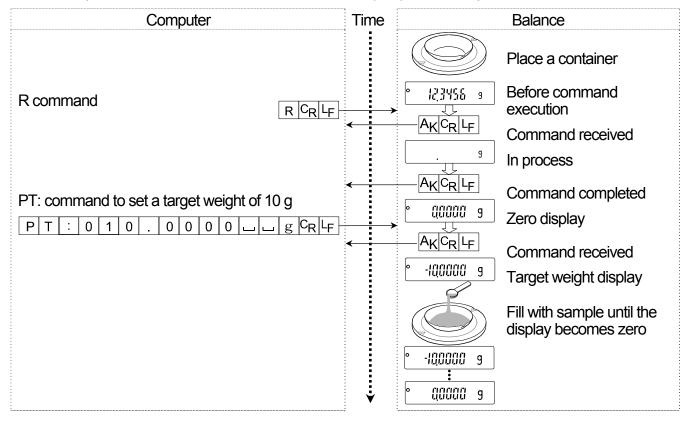

# 17. Key Lock Function

Sending specified commands to the balance will lock the keys on the balance to disable the key functions. This is useful for controlling the balance only by an external device such as a personal computer.

- Even if the keys are locked, key operations using commands are available. For details on commands to operate the keys, refer to "16.3. Commands".
- The key lock state can be checked using the command to request the state of the balance keys.
- □ The key lock state is maintained until the balance receives the key unlocking command or the balance is turned off by removing the AC adapter.
- The key lock function (KL command and LK command) is available for balances with software version
   1.660 or later.

# 17.1. Locking All the Keys

Using the KL command, all the keys on the balance can be disabled.

Command string	Description		
KL:***	KL:000 Unlocks all the keys.		
(Enter 000 or 001 in place of ***.)	KL:001 Locks all the keys		
?KL	Requests the state of the all key lock function.		
	KL,000 All keys unlocked.		
	KL,001 All keys locked.		

# 17.2. Locking the Specified Keys

Entering numerical values \*\*\*\*\* to the LK command will lock specified keys on the balance to disable their key functions. Set the numerical values \*\*\*\*\* by adding the decimal numbers converted from bits assigned to each key as shown below.

Bits	Decimal	Keys
0	1	ON:OFF
1	2	CAL
2	4	MODE
3	8	RANGE
4	16	PRINT
5	32	RE-ZERO
6	64	SELECT
7	128	ION

# Example 1: Locking all the keys excluding the PRINT key

- Add the decimal numbers corresponding to the keys to be locked.
   1 (ON:OFF) + 2 (CAL) + 4 (MODE)+ 8 (RANGE) + 32 (RE-ZERO) + 64 (SELECT) +128 (ION) = 239
- 2 Send the LK command to the balance with the sum that is obtained in Step 1. LK: 00239

## Example 2: Unlocking all the keys


1 Send the LK command to the balance with 0, which indicates no keys to be locked. LK: 00000

Command string	Description			
LK:****	Locks the specified keys.			
(Enter the numerical values	e.g.: LK: 00239 Locks all the keys excluding the PRINT key.			
00000 <b>to</b> 00255 <b>in place</b>				
of ****.)				
?LK	Requests the state of the specified key lock function.			
	e.g.1: LK, 00239 Specified keys locked.			
	e.g. 2: LK, 00000 All keys unlocked.			

# 18. Checking the Balance Software Version

The balance specifications vary depending on the balance software version. Follow the steps below to check the balance software version.

- 1 Unplug the AC adapter from the balance and plug it back in.
- 2 P-\*.\*\*\* is displayed with the software version in place of \*\*\*\*

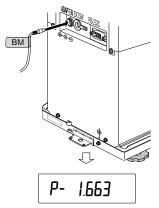

# 19. Maintenance

# 19.1. Cleaning and Handling the Balance

- Clean the balance with a lint free cloth that is moistened with warm water and a mild detergent.
- Do not use organic solvents to clean the balance.
- □ Do not disassemble the balance. Contact the local A&D dealer if the balance needs service or repair.
- Use the original packing material for transportation.
- Consider "3. Installing the Balance and Precautions" when handling the balance.

# 19.2. Error Codes

Display	Error code	Description		
Error 0	EC,E00	Internal error If this error is displayed continuously, contact your local A&D dealer to repair the balance.		
Error 1	EC,E11	Stability error The balance cannot stabilize due to an environmental problem. Check around the pan. Prevent vibration, drafts, temperature changes, static electricity and magnetic fields, from influencing the balance.  Press the CAL key to return to the weighing mode.		
Error 6	EC,E16	Internal mass error Applying the internal mass does not yield a change in the mass value as specified. Confirm that there is nothing on the pan and perform the weighing operation from the beginning again.		
Error 7	EC,E17	Internal mass error The internal mass application mechanism does not function properly. Perform the weighing operation from the beginning again.		
[ FRL F	EC,E20	Calibration weight error The calibration weight is too heavy. Confirm the calibration weight value. Press the CAL key to return to the weighing mode.		
-686-8	EC,E21	Calibration weight error The calibration weight is too light. Confirm the calibration weight value. Press the CAL key to return to the weighing mode.		
<b>£</b> 9		Overload error A sample beyond the balance weighing capacity has been placed on the pan. Remove the sample from the pan.		
- <b>£</b> g		Weighing pan error The mass value is too light. Confirm that the weighing pan and pan support are properly installed. Press the ON:OFF key two times to return to the weighing mode. If the error still persists, calibrate the balance.		
Lo		Sample mass error The balance cannot store the sample for the counting mode or for the percent mode because it is too light. Use a larger sample.		

Display Error code	Description
25 - PC 50 - PC 100 - PC	Unit mass error The sample unit mass for the counting mode is too light. Storing and using it for counting will cause a counting error. Add samples to reach the specified number and press the PRINT key. Pressing the PRINT key without adding samples will shift the balance to the counting mode. But, for accurate counting, be sure to add samples.
<u> K</u> <u>K</u>	Automatic response adjustment zero error  The automatic response adjustment cannot be performed because there is something on the pan. Clear the pan. Press the CAL key to return to the weighing mode.
(Check NG)	Automatic response adjustment unstable error The automatic response adjustment cannot be performed because the mass value is unstable. Check the ambient conditions such as breeze, vibration and magnetic fields, also check the weighing pan. Press the CAL key to return to the weighing mode.
rte Pf	Clock battery error The clock backup battery has been depleted. Press any key and set the time and date. The clock and calendar function works normally as long as the AC adapter is connected to the balance. If this error appears frequently, contact the local A&D dealer.
Lb	Low battery error Voltage of the power supply drops. Confirm whether the correct AC adapter is used.
医III (Blinking)	Memory full The amount of weighing data in memory has reached the maximum capacity. Delete data in memory to store new data. For details, refer to "12. Data Memory".
ि ६॥ (Illuminated)	Memory full The amount of calibration or calibration test data in memory has reached the maximum capacity (50 sets). The data in memory will be deleted automatically to store new data. For details, refer to "12. Data Memory".
	Memory type error Type of memory set in the function table and type of data stored are different. For details, refer to "12. Data Memory".
<u>Err</u>	Memory data error A fatal error occurs in the stored data. To use the memory function, delete all the old data and resolve the error. For details, refer to "12.1. Notes on Using Data Memory".
EC,E00	Communications error A protocol error occurred in communications. Confirm the format, baud rate and parity.
EC,E01	Undefined command error An undefined command was received. Confirm the command.

Display Error code	Description
EC,E02	Not ready A received command cannot be processed.  Examples:  □ The balance received a "ℚ" command, but not in the weighing mode.  □ The balance received a "ℚ" command while processing a RE-ZERO command.  Adjust the delay time to transmit a command.
EC,E03	Timeout error  When the timeout parameter is set to "Ł-UP an", the balance did not receive the next character of a command within the time limit of one second. Confirm the communication.
EC,E04	Excess characters error The balance received excessive characters in a command. Confirm the command.
EC,E06	Format error A command includes incorrect data. Example:  □ The data is numerically incorrect. Confirm the command.
EC,E07	Parameter setting error The received data exceeds the range that the balance can accept. Confirm the parameter range of the command.
Other error code	If an error described above cannot be resolved or other errors are displayed, contact the local A&D dealer.

# 19.3. Other Display



When this indicator blinks, automatic self calibration is required. The indicator blinks when the balance detects changes in ambient temperature. If the balance is not used for several minutes with this indicator blinking, the balance performs automatic self calibration. The blinking duration depends on the operating environment.

Note The balance can be used while this indicator is blinking. We recommend that you perform automatic self calibration for precision weighing.



This is the display to prompt you to enter a password. When the password function is used, the correct password must be entered for the balance to enter the weighing mode.

For details, refer to "15. Password Function".

# 19.4. Checking the Balance Performance and Environment

The balance is a precision instrument. When the operating environment or the operating method is inadequate, correct weighing cannot be performed. Place a sample on the pan and remove it, and repeat this several times. If the balance seems to have a problem with repeatability or to perform improperly, check as described below. If improper performance persists after checking, contact the local A&D dealer for repair.

## Checking that the operating environment or weighing method is proper Operating environment

- Is the weighing table solid enough?
- Is the balance level?
- Is the operating environment free from vibration and drafts?
- Is there a strong electrical or magnetic noise source such as a motor near the balance?

## Weighing method

- Does the weighing pan edge touch anything? Is the weighing pan assembly installed correctly?
- □ Is the RE-ZERO key pressed before placing a sample on the weighing pan?
- Is the sample placed in the center of the weighing pan?
- □ Is the fine range ring installed for weighing with a readability of 0.01 mg for the BM-252?
- □ Is the fine range ring and small fine range ring installed for weighing with a readability of 0.001 mg for the BM-5, BM-5D, BM-20 and BM-22?
  - Is the slide breeze break installed for weighing with a readability of 0.001 mg for the **BM-20** and **BM-22**?
- Has the balance been calibrated using the internal mass?
- Has the balance been warmed up for one hour before weighing?

## Sample and container

- Has the sample absorbed or lost moisture due to the ambient conditions such as temperature and humidity?
- Has the temperature of the container been allowed to equalize to the ambient temperature?
- Is the sample charged with static electricity? Is the built-in ionizer (page 31) used to discharge the sample charged with static electricity.
- Is the sample a magnetic material such as iron? Cautions are required when weighing magnetic materials.

# Checking that the balance performs properly

- Check the balance repeatability using an external weight. Be sure to place the weight in the center of the weighing pan.
- Check the balance repeatability, linearity and calibrated value using an external weight with a known value.
- Check whether the calibration is performed correctly. If an error occurs and the calibration cannot be performed quickly, the balance may not measure correctly.

# 19.5. Asking for Repair

If the balance needs service or repair, contact your local A&D dealer.

The balance is a precision instrument. Use much care when handling the balance and observe the following when transporting the balance.

- Use the original packing material for transportation.
- Remove the weighing pan, pan support, rings and dust plate from the balance.

# 20. Specifications

(Typical at FAST) 10 seconds Approx. 10 seconds App			BM-5	BM-5D	BM-20	BM-22	
Auximum display   5.200084 g   5.20008 g   22.000084 g   1 5.100009 g   1 5.1000009 g   1 5.100009 g   1 5.100009 g   1 5.100009 g   1 5.100009 g   1 5.10	Weighing canacity		52a	_	22 a		
Readability  Repeatability  0.001 mg  10.0025 mg  10.002 mg  10.001 mg  10.0	vveigning capacity		5.2 g		22 y		
Page	Mavin	num display	5 200084 a	5.20008 g	22 000084 a	22.00008 g	
Repeatability (Standard deviation)   Repeatability   \$\frac{\pmathbb{\text{20.001} mg}{\pmathbb{\text{20.004} mg}} \frac{\pmathbb{\text{20.004} mg}{\pmathbb{\text{20.004} mg}} \frac{\pmathbb{\text{20.004} mg}{\pmathbb{\text{20.010} mg}} \frac{\pmathbb{\text{20.010} mg}}{\pmathbb{\text{20.010} mg}} \frac{\pmathbb{\text{20.000} mg}}{\pmathbb{\text{20.010} mg}} \frac{\pmathbb{\text{20.010} mg}}{\text{20.0	IVIGAIII	nam display	3.20000+ g	*1 2.100009 g	22.00000+ g	* <b>1</b> 5.100009 g	
Repeatability (Standard deviation)   **2	Reads	ahility	0.001 mg	0.01 mg	0.001 mg	0.01 mg	
Continued   Cont	rtcaut	ability		0.001 mg	9	0.001 mg	
Linearity				0.01 mg		0.01 mg	
Stabilization time (Typical at FAST)	(S	tandard deviation)	0.001211Ig	*2 0.004 mg	0.0025 mg	* <b>2</b> 0.004 mg	
Stabilization time (Typical at FAST) Approx. Approx. 8 seconds (Typical at FAST) 10 seconds Approx. 10 secon	l inea	ritv	+0.010 ma	±0.02 mg	+0.010 ma	±0.02 mg	
(Typical at FAST) 10 seconds Approx. 10 seconds App	LiiiCai	rity	±0.010111g	±0.010 mg	±0.010111g	±0.010 mg	
Sensitivity drift  Display refresh rate  Environmental sensors  Minimum unit mass  Number of samples  Readability  Display refreence mass  Readability  Display refresh rate  External calibration  Signaphy and some supply & Confirm that the adapter type  Power supply & Confirm that the adapter type  Signaphy refresh rate  10 no or 0.1 mg  10 mg	Stabili	ization time	Approx.	Approx. 8 seconds		Approx. 8 seconds	
Operating environment Display refresh rate  Environmental sensors  Resolution: Temperature ±1.5 °C, Humidity ±10 %, Barometric pressure ±10 hPa Applicable range: 5 °C to 40 °C  Minimum unit mass Number of samples  Headability  O.01 %, O.1 %, 1 % (Depends on the reference mass stored.)  Interface  External calibration weight  Size of loading area for weighing mm  Power supply & AC adapter tyree  Operating environmental 5 °C to 40 °C, 85 % RH or less (No condensation)  Stimes/second or 10 times/second  Resolution: Temperature ±1.5 °C, Humidity ±10 %, Barometric pressure ±10 hPa Applicable range: 5 °C to 40 °C  O.1 mg  O.1 mg  O.1 mg  O.1 mg  O.1 mg  O.25, 50 or 100 pieces  USB and RS-232C  USB and RS-232C  External calibration  Oxide the second of the second of the AC adapter of the local voltage and power of the local voltage		(Typical at FAST)	10 seconds	Approx. 10 seconds	10 seconds	Approx. 10 seconds	
Display refresh rate  Environmental sensors  Resolution: Temperature ±1.5 °C, Humidity ±10 %, Barometric pressure ±10 hPa Applicable range: 5 °C to 40 °C  Minimum unit mass  Number of samples  Minimum 100% reference mass  Readability  0.01 %, 0.1 %, 1 % (Depends on the reference mass stored.)  Interface  USB and RS-232C  External calibration weight  Size of loading area for weighing mm  Size of loading area for weighing mm  Power supply & AC adapter tyree  Power supply & Confirm that the adapter type is correct for the local voltage and power	Sensi	tivity drift	±2 ppm/°C	(10 °C to 30 °C, A	utomatic self calibr	ation: OFF)	
Environmental sensors  Resolution: Temperature ±1.5 °C, Humidity ±10 %, Barometric pressure ±10 hPa Applicable range: 5 °C to 40 °C  O.1 mg  Minimum unit mass  Number of samples  Number of samples  Number of samples  Nound of samples  Number	Opera	ating environment	5 °C t	o 40 °C, 85 % RH c	or less (No condens	sation)	
Applicable range: 5 °C to 40 °C    Document	Displa	ay refresh rate		5 times/second or	10 times/second		
Minimum 100% reference mass   10 mg	. ,		Resolution: Temperature ±1.5 °C, Humidity ±10 %, Barometric pressure ±10 hPa Applicable range: 5 °C to 40 °C				
Minimum 100% reference mass   10 mg	の Minimum unit mass		-				
Interface  USB and RS-232C   5 g 5 g 20 g 20 g  2 g 10 g 10 g  1 g 5 g 5 g  2 g 2 g 10 g  2 g 2 g 10 g  5 g 5 g  5 g 2 g 2 g  1 g 5 g 5 g  2 g 2 g 2 g  1 g 1 g 5 g  2 g 2 g  1 g 1 g  Size of loading area for weighing mm  Weighing pan: \$\phi 25\$  Filter weighing: \$\phi 50\$, \$\phi 95\$  Micro tube holder: \$\phi 12\$  External dimensions  Power supply & Power consumption: Approx. 30 VA (supplied to the AC adapter)  Confirm that the adapter type is correct for the local voltage and power	Cour		10, 25, 50 or 100 pieces				
Interface  USB and RS-232C   5 g 5 g 20 g 20 g  2 g 10 g 10 g  1 g 5 g 5 g  2 g 2 g 10 g  2 g 2 g 10 g  5 g 5 g  5 g 2 g 2 g  1 g 5 g 5 g  2 g 2 g 2 g  1 g 1 g 5 g  2 g 2 g  1 g 1 g  Size of loading area for weighing mm  Weighing pan: \$\phi 25\$  Filter weighing: \$\phi 50\$, \$\phi 95\$  Micro tube holder: \$\phi 12\$  External dimensions  Power supply & Power consumption: Approx. 30 VA (supplied to the AC adapter)  Confirm that the adapter type is correct for the local voltage and power	rcent			10	mg		
External calibration weight    5 g   5 g   20 g   20 g   10 g   10 g   10 g   10 g   5 g   2 g	Pe	Readability	0.01 %, 0.1	%, 1 % (Depends	on the reference n	nass stored.)	
External calibration weight    2g	Interfa	ace	· · · · · · · · · · · · · · · · · · ·				
External calibration weight    2g			5 g	5 g	20 g	20 g	
weight    Size of loading area for weighing mm   Weighing pan: φ25	Evtorr	nal calibration	2 g	2 g	_	_	
Size of loading area for weighing mm  Weighing pan: \$\phi 25\$ Filter weighing: \$\phi 50\$, \$\phi 95\$ Micro tube holder: \$\phi 12\$  External dimensions  Power supply &  AC adapter type  Power danger type  Power consumption: Approx. 30 VA (supplied to the AC adapter )  Confirm that the adapter type is correct for the local voltage and power			ı g	ı g		5 g	
Size of loading area for weighing mm  Weighing pan: \$\phi 25\$ Filter weighing: \$\phi 50\$, \$\phi 95\$ Micro tube holder: \$\phi 12\$  External dimensions  259(W) x 466(D) x 326(H) mm  Power supply &  AC adapter type  Confirm that the adapter type is correct for the local voltage and power						2 g	
Filter weighing: \$\phi 50\$, \$\phi 95\$  Micro tube holder: \$\phi 12\$  External dimensions  Power supply &  AC adapter type  Filter weighing: \$\phi 50\$, \$\phi 95\$  Micro tube holder: \$\phi 12\$  259(W) x 466(D) x 326(H) mm  Power consumption: Approx. 30 VA (supplied to the AC adapter )  Confirm that the adapter type is correct for the local voltage and power				\A/a i albi in arra arra		ı g	
Micro tube holder: φ12  External dimensions  259(W) x 466(D) x 326(H) mm  Power supply &  AC adapter type  Confirm that the adapter type is correct for the local voltage and power			,				
External dimensions  259(W) x 466(D) x 326(H) mm  Power supply &  Confirm that the adapter type is correct for the local voltage and power	weighing mm						
Power supply & Power consumption: Approx. 30 VA (supplied to the AC adapter ) Confirm that the adapter type is correct for the local voltage and power	External dimensions		'				
Confirm that the adapter type is correct for the local voltage and power							
receptacie type.							
Mass Approx. 10 kg	Mass	Approx. 10 kg					

- □ The operating environment with no excessive changes in ambient temperature and humidity, no vibration and drafts, no influence by magnetic fields and static electricity.
- The internal mass may change due to corrosion or other damage caused by the operating environment, or due to aging. Check the internal mass periodically and correct the internal mass value if necessary.
- \*1: Even when a tare weight beyond the precision range is placed, the weighing sample can be weighed in the precision range if using the smart range.
- \*2: Repeatability using a 1 gram weight within the weighing capacity.


		BM-252	BM-200	BM-300	BM-500			
Weigh	ning capacity	250 g	220 g	320 g	520 g			
Maximum display		250.00084 g	220.0084 g	320.0084 g	520.0084 g			
Readability		0.01 mg	0.0 mg					
	atability tandard deviation)	*3 0.03 mg	0.1 mg 0.2 mg					
Linear	rity	±0.10 mg	±0.2 mg	±0.3 mg	±0.5 mg			
Stabili	zation time (Typical at FAST)	Approx. 8 seconds		Approx. 3.5 seconds				
Sensit	tivity drift	±2 ppm/°C	(10 °C to 30 °C, A	utomatic self calibra	ation: OFF)			
Opera	ating environment	5 °C t	o 40 °C, 85 % RH o	or less (No condens	ation)			
Displa	y refresh rate	5 times/second or 10 times/second						
Enviro	onmental sensors	Resolution: Temperature ±1.5 °C, Humidity ±10 %, Barometric pressure ±10 hPa Applicable range: 5 °C to 40 °C						
Counting mode	Minimum unit mass	0.1 mg						
Coul	Number of samples	10, 25, 50 or 100 pieces						
Percent mode	Minimum 100% reference mass		10	mg				
Pe	Readability	0.01 %, 0.1	%, 1 % (Depends	on the reference m	ass stored.)			
Interfa	ace		USB and	RS-232C				
External calibration weight		200 g 100 g 50 g 20 g 10 g	200 g 100 g 50 g	300 g 200 g 100 g 50 g	500 g 300 g 200 g 100 g 50 g			
Size o	of loading area for ing mm	Weighing pan: φ90 Micro tube holder: φ12 (Option)						
Exterr	nal dimensions	259(W) x 466(D) x 326(H) mm						
Power supply & AC adapter type		Power consumption: Approx. 30 VA (supplied to the AC adapter ) Confirm that the adapter type is correct for the local voltage and power receptacle type.						
Mass		Approx. 10 kg						

The operating environment with no excessive changes in ambient temperature and humidity, no vibration and drafts, no influence by magnetic fields and static electricity.

<sup>□</sup> The internal mass may change due to corrosion or other damage caused by the operating environment, or due to aging. Check the internal mass periodically and correct the internal mass value if necessary.

<sup>\*3:</sup> Repeatability using a 100 gram weight within the weighing capacity.

# 20.1. External Dimensions

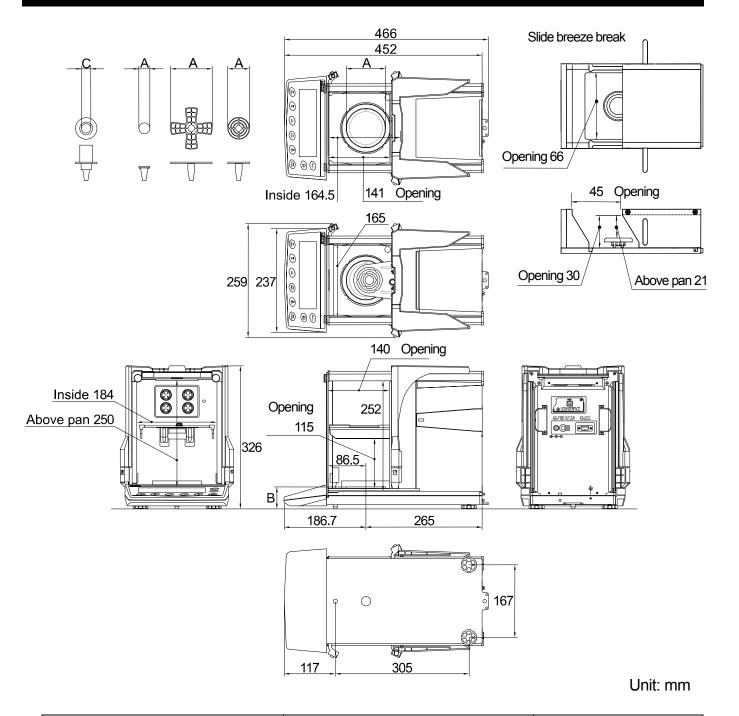

	BM-5, BM-5D BM-20, BM-22	BM-252, BM-200 BM-300, BM-500
A: Diameter of the weighing pan	Weighing pan: φ25 Weighing pan for filter: φ50, φ95	Weighing pan: φ90
B: Height of the weighing pan	54	50
C: The inside diameter of the test tube holder	φ12	φ12 (Option)

# 20.2. Options and Peripheral Equipment

## Ethernet Interface BM-08

- The interface to connect the balance to a LAN.
- Accessory: "WinCT-Plus" data communication software.
  - The software can acquire data from multiple balances connected to a LAN.
  - The software can control these balances with commands.
  - The software can acquire data transmitted from balances. Example: When pressing the PRINT key of the balance, data is output and is acquired by the computer.
  - The stored data can be used with Microsoft EXCEL (if installed).

# | A&D ComponyLimited | Image: Second | Image:

# Pipette Accuracy Tester BM-014 Note This option cannot be used for the BM-5 and BM-5D.

- This option can be used to inspect the volume dispensed from the pipette.
- The accessory evaporation trap is used to prevent the evaporation of the test liquid inside it.
- The WinCT-Pipette software converts the test liquid mass into a volumetric value.



5 mL, 30 mL Sample cup

## Electrode Unit of the Ionizer AX-BM-NEEDLESET (4 pieces)

- Use these electrode units for replacement parts of the ionizer built into the balance.
- Replace all four units with new ones at the same time.
- Refer to "6.5. lonizer" for the replacement.





## Display Cover AX-BM-031 (5 pieces)


This option is the same as the accessory display cover.

# Micro Tube Holder AX-BM-032 (for the BM-5/5D/20/22, standard accessory) AX-BM-034 (for the BM-252/200/300/500)

This option is used when weighing with the general micro tube or test tube under Φ 12 mm.

# Note When using the micro tube holder for the BM-5 and BM-5D:


The micro tube holder is approximately 2 g heavier than the standard pan (weighing pan and pan support). When a micro tube holder is used for the BM-5 and BM-5D, the available weighing range will be up to approximately 3 g including the tare such as a micro tube and test tube.

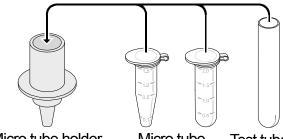

Micro tube holder Micro tube Test tube


## Body Cover AX-BM-033

- Cover during storage to prevent dust and dirt contamination.
- The cover can be used for the balance with the AC adapter and other cables connected.
- When removing the cover, take care so as not to get its edge caught on the balance.
- Specification
  - Material: NylonColor : Light gray

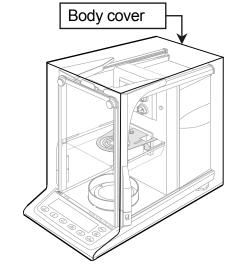

# Slide Breeze Break AX-BM-035 (for the BM-5/5D/20/22, standard accessory)

 This option is attached to the weighing chamber to reduce the effects of the surrounding environment so as to improve weighing accuracy.

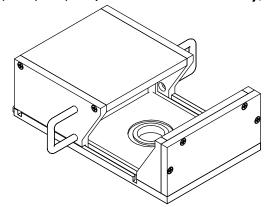

## **Aluminum Pans**

This option is the same as the accessory aluminum pans used for weighing a very small amount of sample.

Name and desc	Order code	View			
Round aluminum analytical pan (L)	φ15	0.8 mL	100 pieces	AX-ROUND-PAN-L	
Round aluminum analytical pan (M)	φ12	0.3 mL	100 pieces	AX-ROUND-PAN-M	
Round aluminum analytical pan (S)	ф8	0.05 mL	100 pieces	AX-ROUND-PAN-S	


## AD-1653 Density Determination Kit

Note This option cannot be used for the BM-5, BM-5D, BM-20 and BM-22.

Weight in air
Weight in air - Weight in liquid x liquid density = sample density

Example  $\frac{10.0000 \text{ g}}{10.0000 \text{ g} - 9.5334 \text{ g}} \times 0.9970 \text{ g/cm}^3 = 21.4 \text{ g/cm}^3$ 

Temperature	Liquid density
0 ℃	0.99984 g/cm <sup>3</sup>
10 °C	0.99970 g/cm <sup>3</sup>
20 °C	0.99820 g/cm <sup>3</sup>
30 °C	0.99565 g/cm <sup>3</sup>

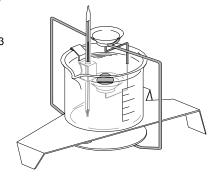

## **Option**

## AD-8127: Multi printer

- Small dot impact printer that connects to the balance via the RS-232C interface.
- Statistical calculation mode, calendar/clock function, interval printing mode, chart printing mode, dump printing mode

## AD-1691: Balance environment analyzer

- During the routine inspection, repeatability and the minimum weighing value can be easily checked.
- Also, this calculates measurement uncertainty and supports environmental evaluation of the balance.
- This analyzer can be easily carried to the balance site, so more than one balance can be controlled using this analyzer.

## AD-1687: Weighing environment logger

- A data logger equipped with 4 sensors for temperature, humidity, barometric pressure and vibration that can measure and store environmental data. When connected to the RS-232C interface of the balance, the AD-1687 can store environmental data along with weighing data. Therefore, it is possible to store data in an environment where a computer cannot be used.
- The stored data can be read to a personal computer using USB. As the AD-1687 is recognized as USB memory, special software is not required to read the data.

## AD-1688: Data logger

■ When connected to the RS-232C interface of the balance, the AD-1688 can store the data in an environment where a personal computer cannot be used.

## AD-8526: Ethernet converter

■ This option can be used to connect the RS-232C interface of the balance to the Ethernet (LAN) port of a computer. This allows management of the balance weighing data with a computer connected to a network.

## AD-8527: Quick USB adapter

■ This option transmits the weighing data directly to software applications such as Excel and Word.

#### AD-8920A: Remote display

Connected to the balance using the RS-232C interface.

#### AD-8922A: Remote controller

■ This option can be connected to the balance using the RS-232C interface and can control the balance remotely. Various options such as comparator output or analog output are available.

## AX-USB-9: USB converter

- Adds a COM port to a personal computer.
- Enables bi-directional communication between the personal computer and the balance when a USB driver is installed.
- Can use serial communication software such as WinCT on a personal computer without COM ports.
- An RS-232C cable is provided to connect the USB converter to the balance.

## AD-1671: Anti-vibration table

- Approximately 27 kg, Artificial stone (Terrazzo).
- Use this table when unstable weighing is caused by vibration coming from the floor.
- Use the AD-8922A remote controller to avoid a weighing error by a slight table tilt that key operation may make.

## AD-1672/AD-1672A Tabletop breeze break (Large)

- Large sized table breeze break
- Protects the balance from wind from sources such as air conditioning or people passing by, thereby reducing balance weighing errors.
- The transparent panel assembly consists of antistatic plastic material that protects the balance from static electricity.
- Samples can be inserted or removed through the opening on the front.
- AD-1672: 680 (W) x 600 (D) x 720 (H) mm, AD-1672A: 680 (W) x 584 (D) x 720 (H) mm

## AD-1676: Tabletop breeze break (Medium)

- A tabletop breeze break of a size best suited for the BM series balances.
- Protects the balance from wind from sources such as air conditioning or people passing by, thereby reducing balance weighing errors.
- The transparent panel assembly consists of antistatic plastic material that protects the balance from static electricity.
- Sliding panels on the right and left side enable samples to be inserted or removed from each side.
- 368(W) x 514(D) x 350(H) mm (Including the handle)

## **AD-1682: Rechargeable Battery**

■ This option allows use of the balance in a place where AC power is not available.

#### AD-1683: DC static eliminator

- This static eliminator eliminates static electricity to prevent weighing errors caused by static electricity during measurement.
- This direct current and windless type of eliminator is the most suitable for precisely weighing powder, etc.
- Small size, light weight

#### AD-1684A: Electrostatic field meter

- This option measures the amount of the static charge on the sample, tare or peripheral equipment and displays the result.
- If those are found to be charged, discharge them using the AD-1683 DC static eliminator.

## AD-1689: Tweezers for calibration weight

■ A pair of tweezers ideally suited for holding calibration weights of 1 g to 500 g.

## AD-8529PC-W: Bluetooth converter (for connecting a personal computer)\*

- This option connects a personal computer to the balance wirelessly (via Bluetooth®) up to 10 m.
- Driver installation is required.

## AD-8529PR-W: Bluetooth converter (for connecting a printer)\*

■ This option connects a printer to the balance wirelessly (via Bluetooth<sup>®</sup>) up to 10 m.

#### \*Note on the AD-8529PC-W and AD-8529PR-W

The two products may require certification for compliance with local laws pertaining to radio broadcast and equipment for wireless communication. Please refer to the A&D website for countries where the products are already certified for use.

## AX-SW137-PRINT: Foot switch print function with a plug

■ The foot switch has the same function as the PRINT key.

## AX-SW137-REZERO: Foot switch rezero function with a plug

■ The foot switch has the same function as the RE-ZERO key.

## **AX-HOLDER-SET: Sample cup holder**

Note This option cannot be used for the BM-5, BM-5D, BM-20 and BM-22.

 A set of holders to secure a sample cup such as a test tube, centrifuge tube or micro tube.



Centrifuge tube secured slantingly



Sample cup holder used as test tube stand



Sample cup holder used as holder for micro tube

# 21. Terms/Index

# 21.1. Terms

Calibration Adjustment of the balance so that it can weigh accurately.

Calibration weight A weight used for calibration.

Data number Numbers assigned sequentially when weighing data or unit weight is stored.

Digit The minimum weighing value available. Used for the balance, one digit is

the smallest mass that can be displayed.

Environment Ambient conditions such as vibration, drafts, temperature, static electricity

and magnetic fields which affect the weighing operation.

External weight A weight that you have.

GLP Good Laboratory Practice.

GMP Good Manufacturing Practice.

Internal mass Built-in calibration weight.

ISO International Organization for Standardization.

Micro tube 2 mL and 1.5 mL micro tubes in the market can

be weighed with the micro tube holder.

1.5 mL 2 mL Micro tube Micro tube holder

Repeatability Variation in measured values obtained when the same mass is placed and

removed repetitively. Usually expressed as a standard deviation.

Example: "Standard deviation = 1 digit" means that a weighing error may occur within

±1 digit in the frequency of approx. 68% when weighing the same mass

repeatedly.

Re-zero To set the display to zero.

Sensitivity drift An affect that a change in temperature causes to the weighing data.

Expressed as a temperature coefficient.

Example: Temperature coefficient = 2 ppm/°C, If a load is 10 g and the temperature

changes by 10°C, the value displayed changes by the following value.

 $0.000\,002\,\text{ppm/°C} \times 10\,\text{°C} \times 10\,\text{g} = 0.0002\,\text{g}$ 

Stabilization time Time required after a sample being placed, until the stabilization indicator

illuminates and the weighing data is displayed.

Tare To cancel the mass value of a container which is not to be included in the

weighing data. Normally, refers to an operation of placing a container and

setting the display to zero.

Target weight An external weight used for calibration test.

Zero point A weighing reference point. Usually refers to the value displayed when

nothing is on the weighing pan.

# 21.2. Index

0	Left, Upper, L	ighting	Stabilization	n indicator	17, 25,	46	Ar-d	Z	ero after	output		49	, 52
<b>⋖</b>	Left, Upper, L	ighting	Process inc				RE-F			•			
***	Left, Upper, E	Blinking	Density mo	ode		17	Auto disp	olay-C	ON			48,	, 51
◀	Left, Lower, L	ighting	Standby inc				Auto print	it				48,	, 52
#	Left, Lower, E	Blinking,	Prior notice of	calibration	17, 38,	99	AX-BM-0	)31	Display	cover			104
door	Right, Upper,	Lighting	Shut side g	lass door		25	AX-BM-0	)32					
in	Right, Upper,	Lighting	Interval mo	de, Stand	by	17	AX-BM-0	)33	Body c	over		8, 1	105
ìņ	Right, Upper,	Blinking	Interval mo	de, Action	· !	17	AX-BM-0	)34	Test tub	e holder		<i>'</i>	104
RESPO	NSE FA	ST MIC	SLOW Res	sponse ind	dicator	17	AX-BM-0	)35	Slide br	eeze breal	k 7	', 12, '	105
(ION)	ION r	mark			17,	31	AX-BM-N	NEEC	LESET	Electroc	de units		104
_	Spac	e mark	<56	, 57, 60, 6	2, 63, 64,	65	AX-HOLE	DER-	SET	Sample	cup holder	·	107
I/O ON:OFF	ON:C	OFF ke	y			17	AX-MX-3	36	Tweez	ers			7
CAL	CAL	key			17,	46	AX-ROUI	ND-F	PAN-L	Aluminum	pan,∳15, 0.8 m	ر '	105
	ION I	key			17,	31	AX-ROUI	ND-F	PAN-M	Aluminum	pan, ∮2, 0.3 ml	ر '	105
MODE	MOD	E key			17, 18,	23	AX-ROUI	ND-F	PAN-S	Aluminum	pan, φ8, 0.05 n	nL′	105
PRINT	PRIN	IT key			17,	46	AX-SW137-	-PRIN	Γ Foot	switch print funct	tion with plug	32, 1	107
1/10d RANGE	RAN	GE ke	у		17,	46	AX-SW137-	-REZE	RO Foot	switch rezero fun	nction with a plug	32, 1	107
◆0◆ RE-ZERO	RE-Z	ERO k	кеу		17,	46	AX-USB-	-9	US	B converte	er	<i>'</i>	106
SELECT	SELE	CT ke	y			17							
8922	AD-8	922A	control		32,	50	Barometr	ric pre	essure			17,	, 33
							685Fnc	Е	nvironn	nent & Disp	olay	46,	, 48
AC a	dapter			8,	9, 11, 13,	16	Baud rate					49,	, 85
ACAI						29	ЬЕЕР	be	ер				. 48
AD-10	653	Densi	ty determina	ation kit	1	05	BM-014	Р	ipette a	ccuracy tes	ster	<i>'</i>	104
AD-10	671	Anti-vi	ibration tabl	e	1	06	BM-08	Е	thernet	interface		4, 1	104
AD-16	72/ AD-	1672A	Tabletop bree	eze break (	L) 13, 1	06	bPS	В	aud rate			49,	, 85
AD-10	676	Tabletop	breeze break	(M)	13, 1	07	Breeze b	reak	ring			7,	, 10
AD-10			argeable ba				<i>ЬЕР</i> г	D	ata bit, p	arity bit		49,	, 86
AD-10	683	DC sta	atic eliminat	or	1	07	Bubble sp	pirit le	evel			.9, 11,	, 14
AD-10	684A	Electro	ostatic field	meter	1	07							
AD-10	687	Weigh	ning environ	ment logg	er 1	06	⁻[RL	Α	utomati	c self calib	ration		. 38
AD-10	688	Data l	ogger		4, 8, 1	06	CRL E	С	alibratio	on error. To	o heavy		. 97
AD-10	689	Tweez	zers for calib	oration we	ight 8, 1	07	-ERL E	С	alibratio	on error. To	o light		. 97
AD-10	691	Weigh	ning environ	ment anal	yzer 1	06	[RL in	C	alibraion u	ising the intern	nal mass	39,	, 60
AD-8	121B	Printe	r		53, 65,	86							
AD-8	127	Multi p	orinter	53, 59, 60	, 61, 87, 1	06							
AD-8	526	LAN c	onverter		1	06					nass weight		
AD-8	527	Quick	<b>USB</b> adapt	er	1	06	Correc	ctina t	the inter	nal mass v	/alue	41	, <del>77</del> . 50
AD-8	529PC	-W B	luetooth cor	nverter	1	07					39, 41, 6		
AD-8	529PR	-W B	luetooth cor	nverter	1	07	Calibratio				,	-,,	
AD-89	920A	Remo	te display		1	06				ng an exter	nal weight.		. 42
AD-89	922A	Remo	te contoroll	er	1	06				-	nal mass		
Air bu	ioyancy	y				15		on tes	•		40, 42, 6		
AK co	ode			9	1, 92, 93,	94	[RLout			•	ternal weight		
Anti-tl	heft de	vice				9	[[ out	-					
Anti-v	vibratior	n table	AD-1671		13, 1	06	[[ in						
ЯР-Ь		Auto p	orint differen	ice	48, 52,	86	EH E					•	•
RP-P		-	orint polarity				[X n[	CHE	ECK NG	}		35,	, 98

[L RdJ Clock	48, 55, 58	Er[d AK, Error code50, 91, 92, 93, 94
ELERr Delete memory data	70, 71	Err Memory data error98
[Lr Initializing the balance	45, 66, 70	Error code50, 91, 93, 97
[Lr Memory type error	98	E-5 External control Input
[Lr [io Initialization, Perform	45	Ethernet
ELr no Initialization, Cancel		Evaporator trap51, 104
Command		External key jack9
Eand Condition		
Container		FAST35, 36
Counting mode		FCC5
CR Carriage return56, 6		Filter
ErLF   Terminator		Fine range ring
	• •	<u> </u>
[5] In Internal mass correction	43, 50	FUL Memory full
	00	
-d- Weighing data		Glass door handle9
d [0000 Initial density		GLP output
d R L R Data memory		GLP 39, 41, 49, 59, 60, 61, 63, 65, 108
Data bits, parity bits		GMP108
Data format	49, 53, 56, 86	Grounding15
Data memory	48, 66	Grounding terminal9
Data memory function	66	
Data number	. 49, 55, 66, 108	Humidity33
Data output	49	·
Data output mode		ıd ID number setting50, 59
Data output pause		ID number
Date		Influence of magnetic fields
d-dt Door Sensor		Influence of static electricity
Decimal point		то Fa GLP output49, 60, 61, 62, 63, 64, 65, 70
Density		Inhibition
Density mode 15		Initializing45
Digit		intelliging
•		Interval mode in
Dimensions	•	, ,
Discharge electrode		Interval time
Display character table		Neutralizing time
Display cover		lon31
Display lock function		ומח <sup>F</sup> חב Ionizer function
Display refresh rate		
d-no Data number		Joint handle9
door Door sensor function		
ժոսե Data output		Key
d5 Fnc Density function		CAL Key
ರ-೬ Weighing data with time and	I date 69	ION Key17, 31 MODE Key17, 18, 23, 25
Dust plate	7, 10	ON:OFF Key17
		PRINT Key17, 46
E Weighing error. Too heavy	97	RANGE Key 17, 46
-E Weighing error. Too light		RE-ZERO Key17, 25, 46
Electrode unit of the ionizer		SELECT Key
Electrostatic field meter AD-1684A		key lock function
Environment		KL command95
Environmental sensor		
	11111 11, 00	Ld in Liquid density input
		110 RM series

Leveling foot9, 11, 14	RsKey Data output mode for WinCT 88
LF Line feed56, 60, 62, 63, 64, 65	רב PF Clock battery error 58, 98
Linearity101, 102	
LK command95, 96	Sensitivity drift101, 102
Lo Sample mass error	Separation plate
Lac na Password input50, 82	Serial interface
LocEnc Password function50, 76	5- <i>id</i> ID number output49, 54, 68, 86
	5 ,F Serial interface49, 52, 53, 86
Magnetic material4, 15, 72	SLOW35, 36
Mass (of the product) 101, 102	Smart range4, 20, 21, 27
Maximum display101, 102	5 <i>Pd</i> Display refresh rate48, 51, 52
Micro tube holder7, 101, 102	Specific gravity73, 75
MID35, 36	Stability
Readability 101, 102	Stability band width48, 51
	Stabilization indicator •
Neutralization31	Stabilization time101, 102, 108
	Standard deviation34
Operating environment101, 102	Standby indicator Left, Lower, Lighting
Report output	5£8r£ Title block of output65
ոսե նո Report output, Start69, 71	Static electricity
ոս Report output, cancel 69, 71	
	5Ł-b Stability band width48, 51
Pan support	5- <i>Ed</i> Time/Date output49, 58, 66, 67, 86
Parity	Stop bit85
PR55 Password	5-ŁP Temperature output
password function 50, 76	Stream mode
Percent mode	75
Permission	£ 25 Initial temperature
Pipette	Tare
PLac Display lock function	Tare weight
Pnt Decimal point	Temperature4, 15, 17, 55, 75, 101, 102
P-an Auto display-ON	Temperature output
Power consumption	<term> Terminator</term>
Power supply	Terminator
Prior notice of calibration	Test tube holder
Processing indicator Left, Upper, Lighting <b>◄</b> 17 Pr L Data output mode48, 52, 66, 86, 87, 89	Time
	Timeout $\xi$ - $U$ P
	とって Zero tracking
PUSE Data output pause49, 86	
rEERLL Display format for memory data 69, 70	Tweezers
,	LJI L Data torrial
rEcEnd End block of output65 Relative Humidity13, 15, 101, 102	Underhook72
Remote Controller AD-8922A	Unit17, 18, 19, 22, 23, 27, 57
rEP LESE Repeatability test	9 Gram17, 18, 19, 22, 23, 27, 37
Repeatability34, 101, 102, 108	<b>"9</b> Milligram18, 19, 23, 57
RESPONSE	PE Counting mode18, 19, 23, 28, 57
Round aluminum pan	% Percent mode18, 19, 23, 30, 57 Ounce (Avoir)18, 19, 23, 30, 57
RS-232C4, 9, 85, 88	©Zt Troy Ounce
RsCom Command mode for WinCT 88	<i>ct</i> Metric carat18, 19, 23, 57
	mom Momme18, 19, 23, 57
RM series 111	

dnt	Pennyweight	
БN	Grain (UK)	
EL	Tael (preset in the factory)	
t	Tola (India)	
m	Messghal	
115	Density mode	22, 57, 73, 75
Un ıE	Unit selection mode	
USB		4, 9, 89
Water eva	aporation	51
Water ter	mperature	74
Weighing	capacity	101, 102
Weighing	data logger AD-1688	4, 8
Weighing	mode	18, 22, 23
Weighing	pan	7
Weighing	pan for filter	7, 10
Weighing	response	35
Weight		
Adiusta	able range	37

42, 108
7, 10
Tools Software 4, 88
inspect pipette 104
connect to LAN 104
58



## A&D Company, Limited

3-23-14 Higashi-Ikebukuro, Toshima-ku, Tokyo 170-0013, JAPAN Telephone: [81] (3) 5391-6132 Fax: [81] (3) 5391-1566

#### A&D ENGINEERING, INC.

1756 Automation Parkway, San Jose, California 95131, U.S.A. Telephone: [1] (408) 263-5333 Fax: [1] (408)263-0119

## A&D INSTRUMENTS LIMITED

Unit 24/26 Blacklands Way, Abingdon Business Park, Abingdon, Oxfordshire OX14 1DY United Kingdom Telephone: [44] (1235) 550420 Fax: [44] (1235) 550485

## A&D AUSTRALASIA PTY LTD

32 Dew Street, Thebarton, South Australia 5031, AUSTRALIA Telephone: [61] (8) 8301-8100 Fax: [61] (8) 8352-7409

## **A&D KOREA Limited** 한국에이.엔.디(주)

서울특별시 영등포구 국제금융로6길33 (여의도동) 맨하탄빌딩 817 우편 번호 07331 (817, Manhattan Bldg., 33. Gukjegeumyung-ro 6-gil, Yeongdeungpo-gu, Seoul, 07331 Korea ) 전화: [82] (2) 780-4101 팩스: [82] (2) 782-4264

#### OOO A&D RUS

## ООО "ЭЙ энд ДИ РУС"

121357, Российская Федерация, г.Москва, ул. Верейская, дом 17 (Business-Center "Vereyskaya Plaza-2" 121357, Russian Federation, Moscow, Vereyskaya Street 17 ) тел.: [7] (495) 937-33-44 факс: [7] (495) 937-55-66

#### A&D INSTRUMENTS INDIA PRIVATE LIMITED

ऐ&डी इन्स्ट्रयमेन्ट्स इण्डिया प्रा0 लिमिटेड

509, उद्योग विहार , फेस -5, गुड़गांव - 122016, हरियाणा , भारत

( 509, Udyog Vihar, Phase-V, Gurgaon - 122 016, Haryana, India ) फोन : 91-124-4715555 फैक्स : 91-124-4715599