

# INSTRUCTION MANUAL

MC-1000 MC-6100



1WMPD4002204A

## This Manual and Marks

All safety messages are identified by the following, "WARNING" or "CAUTION", of ANSI Z535.4 (American National Standard Institute: Product Safety Signs and Labels). The meanings are as follows:

<b>A</b> WARNING	A potentially hazardous situation which, if not avoided, could result in death or serious injury.
	A potentially hazardous situation which, if not avoided, may result in minor or moderate injury.



This is a hazard alert mark.

- This manual is subject to change without notice at any time to improve the product.
- □ The contents of the product specifications and this manual are subject to change without any obligation on the part of the manufacturer.
- Under the copyright laws, the software (program) described in it are copyrighted, with all rights reserved.
   The software may be installed into one computer and may not be installed into other computers without the prior written consent of A&D Company. Copying includes translation into another language, reproduction, conversion, photocopy and offer or loan to another person.

 Microsoft, Windows, Word and Excel are registered trademarks of the Microsoft Corporation.

© 2012 A&D Company, Limited All rights reserved.

No part of this publication may be reproduced, transmitted, transcribed, or translated into any language in any form by any means without the written permission of A&D Company, Limited.

## CONTENTS

## **Basic Operation**

1.	. INTRODUCTION	3
	1-1 About This Manual	
	1-2 Features	
	1-3 Compliance	3
2.	. UNPACKING THE BALANCE	4
	2-1 MC-1000	
	2-2 MC-6100	
	2-3 Installing the Balance	
3.	. PRECAUTIONS	10
	3-1 Before Use	10
	3-2 During Use	11
	3-3 After Use	
	3-4 Power Supply	
4.	. DISPLAY SYMBOLS AND KEY OPERATION	13
5.	. WEIGHING UNITS	14
6.	. WEIGHING	16
	6-1 Selecting a Weighing Unit (Mode)	
	6-2 Basic Weighing	

## Adapting to the Environment

7. RESPONSE ADJUSTMENT / SELF CHECK FUNCTION	18
7-1 Automatic Response Adjustment / Self Check Function 7-2 Manual Response Adjustment	
8. CALIBRATION	20
8-1 Calibration Group	
8-2 Automatic Self Calibration (Calibration due to changes in temperature)	
8-3 Calibration Using the Internal mass (One-touch calibration)	
8-4 Calibration Using an External Weight	
8-5 Calibration Test Using an External Weight	
8-6 Correcting the Internal Mass Value	

### **Functions**

9. FUNCTION SWITCH AND INITIALIZATION	28
9-1 Permit or Inhibit	
9-2 Initializing the Balance	

10. FUNCTION TABLE	30
10-1 Structure and Sequence of the Function Table	30
10-2 Display and Keys	30
10-3 Details of the Function Table	
10-4 Description of the Class "Environment, Display"	35
10-5 Description of the Item "Data output mode"	
10-6 Description of the Item "Data format"	
10-7 Description of the Data Format Added to the Weighing Data	
10-8 Data Format Examples	40
11. ID NUMBER AND GLP REPORT	41
11-1 Setting the ID Number	41
11-2 GLP Report	42

## **RS-232C Serial Interface**

12. RS-232C SERIAL INTERFACE / EXTERNAL INPUT	45
13. CONNECTION TO PERIPHERAL EQUIPMENT	47
13-1 Connection to the AD-8121B Printer	
13-2 Connection to a Computer	
13-3 Using Windows Communication Tools (WinCT)	
14. COMMANDS	50
14-1 Command List	
14-2 Acknowledge Code and Error Codes	
14-3 Control Using CTS and RTS	
14-4 Settings Related to RS-232C	
15. EXTENDED FUNCTION	53
15-1 Description of "Averaging range" and "Averaging time"	

### Maintenance

16. MAINTENANCE	56
17. TROUBLESHOOTING	56
17-1 Checking the Balance Performance and Environment 17-2 Error Codes 17-3 Asking For Repair	
18. SPECIFICATIONS	
19. OPTIONS	62
20. EXTERNAL DIMENSIONS	63
21. TERMS/INDEX	64

## 1. INTRODUCTION

This manual describes how the MC Series Mass Comparator Balances, MC-1000 and MC-6100, work, and how to get the most out of them in terms of performance.

Read this manual thoroughly before using the balance and keep it at hand for future reference.

For other functions and operations that this manual does not describe, refer to the GX series instruction manual.

## **1-1 About This Manual**

This manual consists of the following five parts:

Basic operation	Describes precautions on handling the balance, balance construction and basic balance operation.
Adapting to the environment	Describes response adjustment, calibration and calibration test.
Functions	Describes various functions of the balance.
RS-232C serial interface	Describes the interface which transmits data and controls the balance.
Maintenance	Describes maintenance, error codes, troubleshooting, specifications and options.

### **1-2 Features**

- Display resolution, one digit greater than a standard balance. This allows management of OIML class F1 weights.
- Capable of weighing small amounts of powdery or liquid material, even with a massive tare.
- When used as a mass comparator, the balance can achieve even more precise weighing, by using the optional auto-centering pan (sold separately), which reduces eccentric loading errors.

## **1-3 Compliance**

#### **Compliance with FCC Rules**


Please note that this device generates, uses and can radiate radio frequency energy. This device has been tested and has been found to comply with the limits of a Class A computing device pursuant to Subpart J of Part 15 of FCC rules. These rules are designed to provide reasonable protection against interference when this device is operated in a commercial environment. If this unit is operated in a residential area, it may cause some interference and under these circumstances the user would be required to take, at his own expense, whatever measures are necessary to eliminate the interference.

(FCC = Federal Communications Commission in the U.S.A.)

## 2. UNPACKING THE BALANCE

- The balance is a precision instrument. Unpack the balance carefully. Keep the packing material to be used for transporting the balance in the future.
- The packing contents depend on the balance model. See the illustrations to confirm that everything is contained..


## 2-1 MC-1000

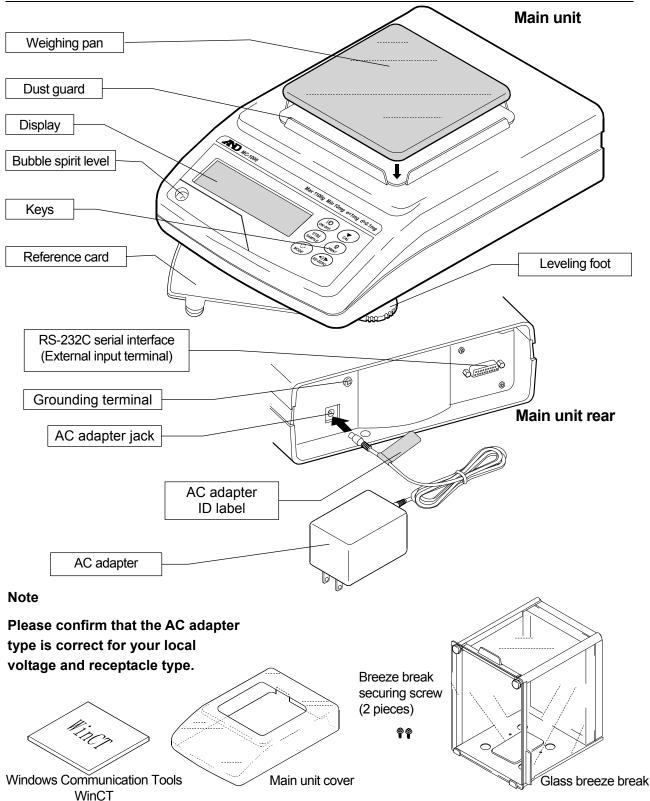


#### Glass breeze break installation procedure

#### Caution: During installation, do not apply excessive force to the balance.


 Remove the weighing pan, pan support receptors, and dust guard from your balance. To remove the dust guard, remove the two screws that secure it to the balance.
 Weighing pan
 Pan support receptor (4 pieces)

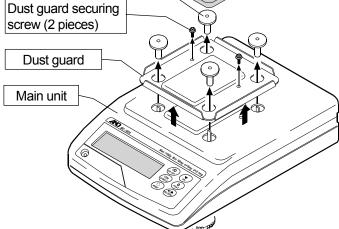

 Secure the breeze break to the main unit using two screws provided with the breeze break.
 Replace the pan support receptors and weighing pan.

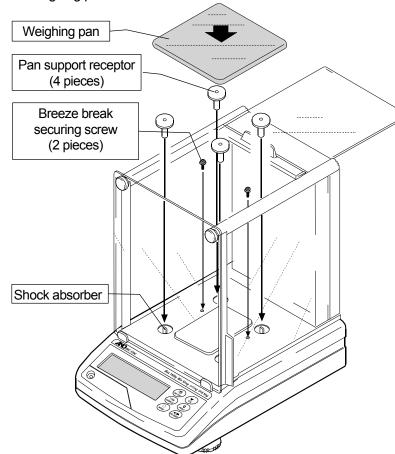

2. Place the breeze break on the main unit

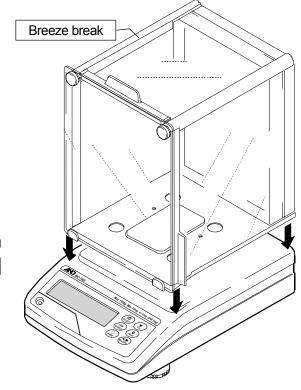

• When installing the pan support receptors:

If the pan support receptors are hard to push in, hold the head of the receptor as shown below and push it in while rotating the head to the right and left.




- Keep the removed dust guard and dust guard securing screws (2 pieces).
- To reinstall the dust guard after removing the breeze break, use the two screws to secure the dust guard.
- The dust guard securing screws and breeze break securing screws are different in length.


#### Auto-centering pan (AX-MC1000PAN) installation procedure

Caution: During installation, do not apply excessive force to the balance.


- When used as a mass comparator, the balance can achieve even more precise weighing, by using the auto-centering pan (sold separately), which reduces eccentric loading errors.
- After completing Steps 1 and 2 of the glass breeze break installation procedure in the previous section, secure the optional AX-MC1000PAN breeze break using the provided securing screws. At this time make sure to place the side of the breeze break with the screws at the back. Attach the pan support receptors to the shock absorbers.

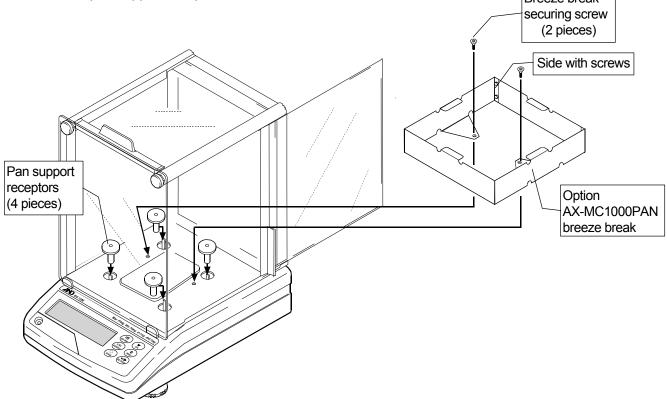

2. Place the auto-centering pan guide on the pan support receptors, and then place the auto-centering pan on the guide.

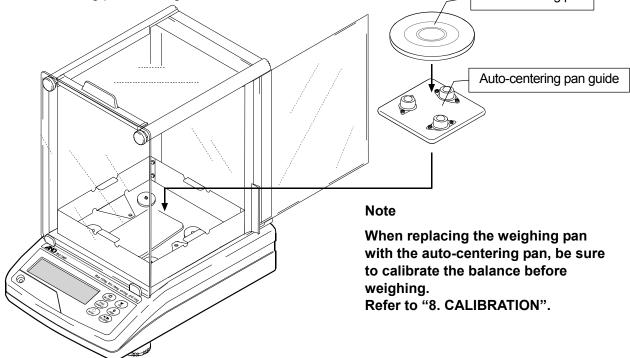


## 2-2 MC-6100

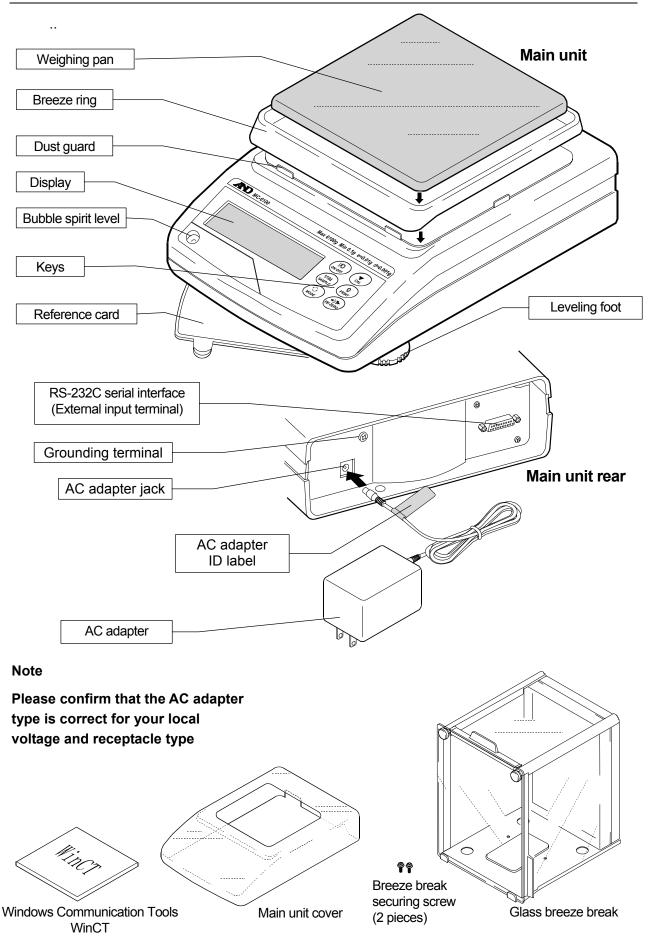


#### Glass breeze break installation procedure

#### Caution: During installation, do not apply excessive force to the balance.


 Remove the weighing pan, breeze ring, pan support receptors, and dust guard from your balance. To remove the dust guard, remove the two screws that secure it to the balance.

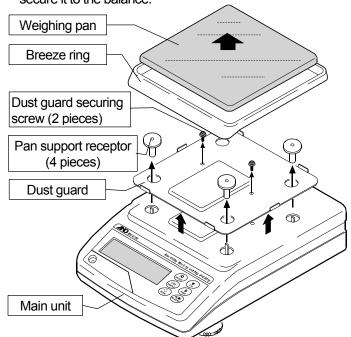

 Secure the breeze break to the main unit using two screws provided with the breeze break.
 Replace the pan support receptors and weighing pan.

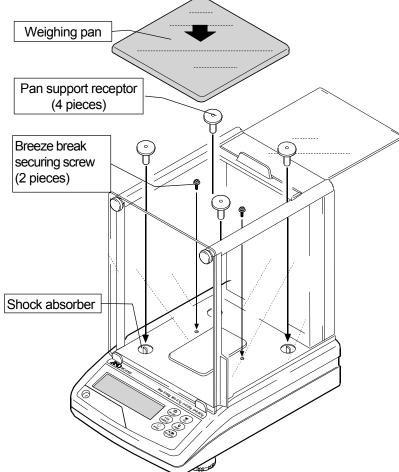

2. Place the breeze break on the main unit

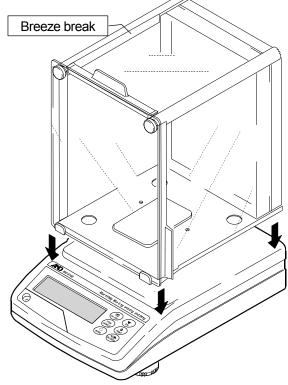

• When installing the pan support receptors:


If the pan support receptors are hard to push in, hold the head of the receptor as shown below and push it in while rotating the head to the right and left.




- Keep the removed dust guard, dust guard securing screws (2 pieces) and breeze ring.
- To reinstall the dust guard after removing the breeze break, use the two screws to secure the dust guard.
- The dust guard securing screws and breeze break securing screws are different in length.

#### Auto-centering pan (AX-MC6100PAN) installation procedure

- Caution: During installation, do not apply excessive force to the balance. • When used as a mass comparator, the balance can achieve even more precise weighing, by using the auto-centering pan (sold separately), which reduces eccentric loading errors.
- 1. After completing Steps 1 and 2 of the glass breeze break installation procedure in the previous section, secure the optional AX-MC6100PAN breeze break using the provided securing screws. At this time make sure to place the side of the breeze break with the screws at the back. In addition, place the breeze break toward the back of the balance as much as possible to prevent the breeze break from touching the auto-centering pan guide.

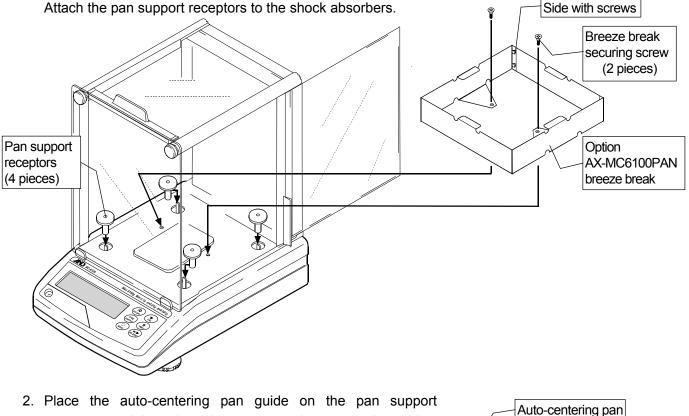

receptors, and then place the auto-centering pan on the guide.

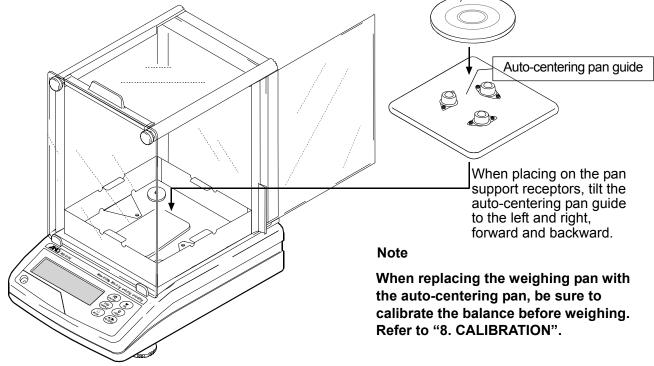

## 2-3 Installing the Balance

Install the balance as follows:


- 1. Refer to "3-1 Before Use" for installing the balance. Place the balance on a solid weighing table.
- 2. Assemble the balance as shown in "2. UNPACKING THE BALANCE".
- 3. Adjust the leveling feet to level the balance. Confirm it using the bubble spirit level.
- 4. Confirm that the adapter type is correct for the local voltage and power receptacle type.
- 5. Connect the AC adapter to the balance. Warm up the balance for at least 30 minutes with nothing on the weighing pan.

## 3. PRECAUTIONS

To get the optimum performance from the balance and acquire accurate weighing data, note the following:

### 3-1 Before Use

- The MC series is a high-precision balance with a resolution of 1/10000000. It is prone to changes in temperature and pressure, static electricity, drafts and vibration of the installation site. During calibration using an internal mass, it resolves up to one digit lower than the minimum weighing value. Therefore, be sure to place the balance in a stable environment.
- The best operating temperature is about 20°C±2°C at 45-60% relative humidity.
- Install the balance where it is not exposed to direct sunlight and it is not affected by heaters or air conditioners.
- Install the balance where it is free of dust.
- Install the balance away from equipment which produces magnetic fields.
- Install the balance in a stable place avoiding vibration and shock. Corners of rooms on the first floor are best, as they are less prone to vibration.
- The weighing table should be solid and free from vibration, drafts and as level as possible.
- Level the balance by adjusting the leveling feet and confirm it using the bubble spirit level.
- If static electricity is a problem at the installation site, use the electrostatic field meter and the static eliminator.
- Ensure a stable power source when using the AC adapter.

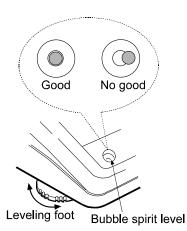


- Warm up the balance for at least 30 minutes. Plug in the AC adapter as usual.
- Calibrate the balance periodically for accurate weighing.
- Before using the balance for the first time or after having moved it to another location, warm up the balance for six hours or more to allow the balance to reach equilibrium with the room temperature, and then calibrate the balance.


#### Caution

#### Do not install the balance where flammable or corrosive gas is present.

### 3-2 During Use

- Static charge may cause weighing errors. When the ambient humidity is less than 45%RH, insulators such as plastic or glass are prone to static electricity. Discharge static electricity from the material to be weighed (hereinafter referred to as sample). Ground the balance and try the following:
  - Eliminate the static electricity by using an optional static eliminator, AD-1683,
  - Or try to keep the ambient humidity above 45%RH.
  - Or use a metal shield case.
  - Or wipe a charged plastic sample with the wet cloth.
- This balance uses a strong magnet as part of the balance assembly, so please use caution when weighing magnetic materials such as iron. If there is a problem, use the underhook on the bottom of the balance to suspend the material away from the influence of the magnet.

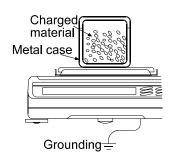

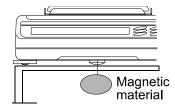


- Cancel the temperature difference between a sample, tare and the environment. When a sample is warmer (cooler) than the ambient temperature, the sample will be lighter (heavier) than the true weight. This error is due to a rising (falling) draft around the sample.
- Make each weighing gently and quickly to avoid errors due to changes in the environmental conditions.
- Do not drop things upon the weighing pan, or place a sample on the pan that is beyond the balance weighing capacity. Place a sample in the center of the weighing pan.
- Do not use a sharp instrument such as a pencil to press the keys. Use your finger only.
- Press the RE-ZERO key before each weighing to prevent possible errors.
- Take into consideration the affect of air buoyancy on a sample when more accuracy is required.
- Keep the balance interior free of dust and foreign materials.

## 3-3 After Use

- Avoid mechanical shock to the balance.
- Do not disassemble the balance. Contact the local A&D dealer if the balance needs service or repair.
- Do not use organic solvents to clean the balance. Clean the balance with a lint free cloth that is moistened with warm water and a mild detergent.
- Protect the internal parts of the balance from liquid spills and excessive dust.

## 3-4 Power Supply


- Do not remove the AC adapter while the internal mass is in motion, for example, right after the AC adapter is connected, or during calibration using the internal mass.
   If the AC adapter is removed under the conditions described above, the internal mass will be left unsecured, that may cause mechanical damage when the balance is moved.
   Before removing the AC adapter, press the ON:OFF key and confirm that zero is displayed.
- When the AC adapter is connected, the balance is in the standby mode if the standby indicator is on (refer to "4. DISPLAY SYMBOLS AND KEY OPERATION"). This is a normal state and does not harm the balance. For accurate weighing, warm up the balance for at least 30 minutes before use.

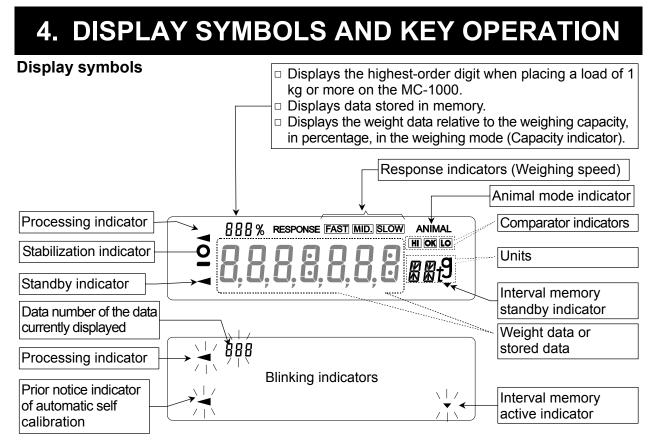

#### Key operation

Key operation affects how the balance functions. The basic key operations are:

"Press and release the key immediately" or "Press the key"
 = normal key operation during measurement

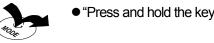




Key	When pressed	When pressed and held		
J/C ON:OFF	off. The weighing mode is enabled when the	indicator is displayed when the display is turned he display is turned on. e key during operation will interrupt the operation		
	and turn the display off.			
1/10d	In the weighing mode, turns the minimum weighing value on and off.	Enters the function table mode. Refer to "10. FUNCTION TABLE".		
SAMPLE	In the counting or percent mode, enters the sample storing mode.			
	No function at the factory setting	Performs response adjustment and self check.		
MODE	Switches the weighing units when units other than "g" are stored in the function setting.			
CAL	Performs calibration of the balance using the internal mass.	Displays calibration-related menu.		
	Stores the weighing data in memory or	No function at the factory setting		
	outputs to a printer or personal computer using the RS-232C interface (Factory setting), depending on the function settings.	By changing the function setting: Outputs "Title block" and "End block" for GLP report. (Refer to "11-2 GLP Report".) Displays the data memory menu.		
+0/T+ RE-ZERO	Sets the display to zero.			

## 5. WEIGHING UNITS

With the balance, only the unit "g" (gram) was set at the factory.

The following weighing units and weighing modes are available for selection:

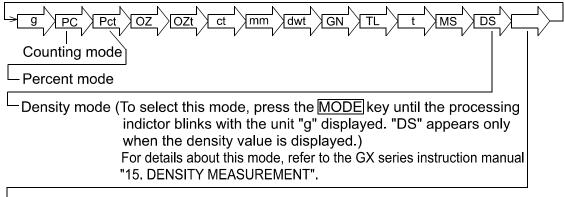

Programmable-unit (No unit displayed. For details, refer to the GX series instruction manual "14. PROGRAMMABLE-UNIT".)

A unit or mode can be selected and stored in the function table. For details, refer to the GX series instruction manual "5-2 Storing Units".

If a weighing mode (or unit of weight) has been turned off, that mode or unit will be missing in the sequence. Tael has four varieties, one of which can be selected and installed at the factory.

Name (unit, mode)	Abbrev.	Display	Function table (Storing mode)	Conversion factor 1 g =
Gram	g	g	g	1 g
Counting mode	PC	ΡĘ	PE	
Percent mode	Pct	Pct	$P_{\subseteq t}$	
Ounce (Avoir)	OZ	02	02	28.349523125 g
Troy Ounce	OZt	ΩZt	0 Z t	31.1034768 g
Metric Carat	ct	C t	c t	0.2 g
Momme	mm	ה הו	ריז ריז	3.75 g
Pennyweight	dwt	dint	dint	1.55517384 g
Grain (UK)	GN	БN	БN	0.06479891 g
Tael (HK general, Singapore)				37.7994 g
Tael (HK jewelry)	TL	ΤL	ΤL	37.429 g
Tael (Taiwan)		· <u>·</u>	· <u>-</u>	37.5 g
Tael (China)				31.25 g
Tola (India)	t	t	t	11.6638038 g
Messghal	MS	MS	M5	4.6875 g
Density mode	DS	jat g	D5	
(See note below)		∬ is used to		
		show the density.		
Programmable-unit (Multi-unit)	Mlt		ML t	

For details about the units and modes, see the table below:

Note: The blinking processing indicator with "g" indicates that the density mode is selected.

- The tables below indicate the weighing capacity and the minimum display for each unit, depending on the balance model.
- When a measurement unit other than gram is used, it is not possible to weight up to capacity
  of the balance because the display does not have sufficient digits.
  Use the MC series balance within the values shown in the tables below.

MC-1000 Unit Capacity Gram Minimum display Gram 1100 g 1100 g 0.0001 Ounce (Avoir) 19.7 OZ 560 g 0.000005 **Troy Ounce** 19.2 OZt 600 g 0.000005 Metric Carat 1950 ct 0.0005 390 g Momme 197 mom 740 g 0.00005 Pennyweight 707 dwt 1100 g 0.0001 16975 GN Grain (UK) 1100 g 0.002 Tael (HK general, Singapore) 19.8 TL 750 g 0.000005 Tael (HK jewelry) 19.7 TL 740 g 0.000005 Tael (Taiwan) 19.7 TL 740 g 0.000005 600 g Tael (China) 19.2 TL 0.000005 Tola (India) 94.3 t 1100 g 0.00001 Messghal 192 MS 900 g 0.00005

	MC-6100			
Unit	Capacity	Gram	Minimum display	
Gram	6100 g	6100 g	0.001	
Ounce (Avoir)	197 OZ	5600 g	0.00005	
Troy Ounce	196 OZt	6100 g	0.00005	
Metric Carat	19500 ct	3900 g	0.005	
Momme	1626 mom	6100 g	0.0005	
Pennyweight	3922 dwt	6100 g	0.001	
Grain (UK)	94137 GN	6100 g	0.02	
Tael (HK general, Singapore)	161 TL	6100 g	0.00005	
Tael (HK jewelry)	162 TL	6100 g	0.00005	
Tael (Taiwan)	162 TL	6100 g	0.00005	
Tael (China)	195 TL	6100 g	0.00005	
Tola (India)	522 t	6100 g	0.0001	
Messghal	1301 MS	6100 g	0.0005	

## 6. WEIGHING

## 6-1 Selecting a Weighing Unit (Mode)

Press the MODE key to select a unit or mode for weighing.

The unit "g" (gram) was set at the factory.

To use other units, select and store units and displaying order in the function setting of "Un L". For details on weighing unit storing procedure, refer to the GX series instruction manual "5-2 Storing Units".

## 6-2 Basic Weighing

#### For more stable weighing

To reduce the influence of drafts and vibration, set the following function settings as below.
 "Condition (*Land*)" of "Environment, Display (*bR5Fnc*)" to "Slow (*2*)"
 "Filter (*F IL*)" of ""Environment, Display (*bR5Fnc*)" to "Used (*1*)"

#### **Function settings**

Refer to "10. FUNCTION TABLE" on page 30 to check or change the function settings.

Class	Item and Parameter		eter	Description	
bASFnc	Cond	Condition	2	Slow response rate, stable value S	LOW
Environment Display	F IL	Filter	Ι	Used	

#### When using as a mass comparator

• To avoid eccentric loading errors, place the sample in the center of the weighing pan.

As an option, auto-centering pans are available: AX-MC1000PAN for the MC-1000 and AX-MC6100PAN for the MC-6100.

Using an AD-8922A remote controller, which is sold separately, the balance can be controlled remotely by the AD-8922A key operations in the same way as when the CAL or RE-ZERO key of the balance is pressed. For the connection procedure between the balance and the AD-8922A, refer to the AD-8922A instruction manual.

- Take measures against causes of weighing error at the installation site, such as changes in temperature, atmospheric pressure, drafts, vibration and static electricity. Perform weighing operations in a stable environment.
- The table below lists the weight class and recommended measuring range for each model of the MC series. The measuring range is determined so that the balance repeatability is to be less than one third of the maximum permissible error for each weight class.

	Model	N	1 C -	10(	0 0	Ν	ΛС-	61	00
	Class	F 1	F 2	M 1	M 2	F 1	F 2	M 1	M 2
	5 kg								
	2 kg					•			
	1 kg						•		
	500 g	▼							
	200 g							V	
	100 g								
	50 g		•						V
Weight	20 g								
(Displayed value)	10 g								
	5 g								
	2 g								
	1 g			•					
	500 mg								
	200 mg								
	100 mg				<b>V</b>				
			]F 1		F 2		M 1		M 2

Weight class and recommended measuring range

#### When building into a system

- When a special weighing pan is to be designed, the weight of the pan and the material to be weighed should not exceed the weighing capacity of the balance.
   To reduce influences of static electricity and magnetism, use materials other than resin and magnetic material such as iron.
- There is a function available to maintain the previous weight value in non-volatile memory, even if the AC adapter is removed.
   When "Zero upon power-on (*P-tr*)" of "Environment, Display (*bR5Fnc*)" is set to "*I*", the previous weight value is displayed upon power-on.
   For details, refer to "Zero upon power-on" on page 35.
- There is a function available to perform span calibration only, when performing calibration with a tare on the weighing pan.
   When "Span calibration (5Pn)" of "Environment, Display (bR5Fnc)" is set to "1", span calibration using the internal mass is performed, with a tare on the weighing pan.
   For details, refer to "Span calibration" on page 35.
- To set a higher response rate or to batch-weigh small amounts of material, such as a powdery material, refer to "15. EXTENDED FUNCTION".

## 7. RESPONSE ADJUSTMENT / SELF CHECK FUNCTION

This function detects the influence on weighing that is caused by drafts and/or vibration at the place where the balance is installed and sets the response characteristic automatically. When this function is selected, the balance self-checks the performance at the same time.

Two modes of response adjustment are available: automatic and manual.

The function has three rates as follows:

0.0000 9

0

Changing the response rate changes the display refresh rate.

0 0		0 1 3	
Indicator	Parameter	Response characteristic	Display refresh rate
FAST	Cond D	Fast response, Sensitive value	If the response rate is changed as follows:
MID.	Cond I		MID. or SLOW→ FAST =10 times/second
SLOW	[ond 2	Slow response, Stable value	FAST → MID. or SLOW = 5 times/second

#### Notes

- □ To set the refresh rate of 5 times/second when the response rate is FAST or 10 times/second when the response rate is MID. or SLOW, change the "Display refresh rate (5Pd)" parameter of "Environment, Display (b用5Fnc)" in the function table. For details, to "10. FUNCTION TABLE".
- □ If the weight value is not stabilized due to drafts or vibration when "[and 0 " is selected by the automatic response adjustment, change the parameter of "[and " manually in the function table.

## 7-1 Automatic Response Adjustment / Self Check Function

This function automatically updates the response adjustment by analyzing the influence of the environment on the weighing data and also self-checks the balance performance using the internal mass.

#### Operation


- 1 Press and hold the MODE key until RESPONSE is displayed, then release the key.
- 2 The balance automatically starts to check the balance performance and sets the response characteristic.

#### Caution

## Do not allow vibration or drafts to affect the balance during adjustment.

3 After automatic adjustment, the balance displays the updated response indicator and returns to the weighing mode. The response indicator remains displayed for about 30 seconds.

The example above indicates that the result of the self check is good and MID. is selected as the response rate.

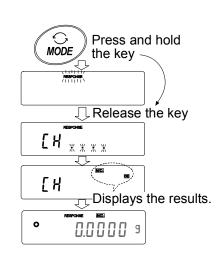

Notes

- □ If improper performance is found in the self check, the balance displays [H n□]. Contact the local A&D dealer for repair.
- If the automatic response adjustment fails, the balance displays <u>[H n[]</u>. Check the ambient conditions such as breeze and vibration, also check the weighing pan. Then, perform the adjustment again. To return to the weighing mode, press the <u>CAL</u> key.
- If the automatic response adjustment is not helpful, try the manual response adjustment.

## 7-2 Manual Response Adjustment

This function manually updates the response adjustment.

#### Operation


1 Press and hold the MODE key until RESPONSE is displayed, then release the key.

Press the MODE key again quickly.

- 2 Press the MODE key to select a rate of the response adjustment. Either FAST, MID. or SLOW can be selected.
- 3 After a few seconds of inactivity the balance displays *End*. Then, it returns to the weighing mode and displays the updated response indicator. The response indicator remains displayed for about 30 seconds.

#### Note

The response adjustment can be changed at "Condition (Land)" of "Environment, Display ( bR5Fnc )" in the function table. For details, refer to "10. FUNCTION TABLE".

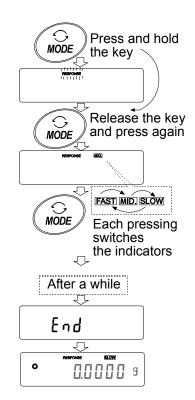

## 8. CALIBRATION

## 8-1 Calibration Group

<ul> <li>Automatic self calibration (calibration due to changes in temperature</li> </ul>	•	Automatic self calib	pration (calibratio	n due to change	s in temperature)
---	---	----------------------	---------------------	-----------------	-------------------

- Calibration using the internal mass (one-touch calibration)
- Calibration using an external weight

Calibration test

Calibration

• Calibration test using an external weight (Calibration test does not perform calibration.)

Correction of the internal mass value

#### Terms

The following terms are defined as follows:

· · · · · · · · · · · · · · · · · · ·		
Internal mass	=	Built-in calibration weight
External weight	=	A weight that you have. Referred to as a calibration weight when used for calibration.
Calibration weight	=	A weight used for calibration
Target weight	=	An external weight used for calibration test

#### Caution

- Calibration adjusts the balance for accurate weighing.
  - Besides periodic calibration and before each use, perform calibration when:
  - the balance is installed for the first time.
  - the balance has been moved.
  - the ambient environment has changed.
- Do not allow vibration or drafts to affect the balance during calibration.
- To output the data for GLP using the RS-232C interface, set "GLP output ( InFo )" of "Data output ( dout )". For details, refer to "10. FUNCTION TABLE". Time and date are added to the GLP report. If the time or date is not correct, adjust them. For details, refer to the GX series instruction manual, "10-9 Clock and Calendar Function".
- Calibration test is available only when "GLP output ( <code>\_\_\_\_\_\_b )</code>" of "Data output ( <code>doub</code> )" is set to " *I* " or " *2* ".
- The calibration and calibration test data can be stored in memory. To store them, set "Data memory ( dRLR )" to " 3". For details, refer to the GX series instruction manual, "12. DATA MEMORY".

#### Caution on using an external weight

• The accuracy of an external weight can influence the accuracy of weighing. Select an appropriate weight as listed below:

Model	Usable calibration weight	Adjustable range
MC-1000	<b>1000g</b> , 900g, 800 g, 700 g, 600 g, 500 g, 400 g, 300 g, 200 g	-0.0150 g to +0.0159 g
MC-6100	6000 g, <b>5000 g</b> , 4000 g, 3000 g, 2000 g	-0.150 g to +0.159 g

The calibration weight in bold type: factory setting

The calibration weight value can be adjusted within the range above.

#### Display



• This indicator means "the balance is measuring calibration data". Do not allow vibration or drafts to affect the balance while this indicator is displayed.

## 8-2 Automatic Self Calibration (Calibration due to changes in temperature)


This function automatically calibrates the balance when the balance detects an ambient temperature change. If GLP output is selected in the function table, the balance outputs the calibration report or stores the data in memory. Automatic self calibration functions even if the display is turned off (standby state).

#### Caution

If something is on the weighing pan, the balance judges that it is in use and does not perform automatic self calibration. To maintain the calibrated state, keep the weighing pan clear while not in use.

The displays shown below are related to the automatic self calibration.

Indicates that the balance detects a change in ambient temperature and automatic self calibration will start. If the balance is not used for a few minutes with this indicator blinking, the balance performs automatic self calibration. The blinking duration depends on the environment.

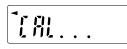

Indicates that the balance is measuring calibration data. Do not allow vibration or drafts to affect the balance while this indicator is displayed. After calibration, the balance returns to indicate the previous display.

#### Note

The balance can be used while the indicator blinks. But, it is recommended that to maintain the accuracy, stop using the balance and confirm that there is nothing on the pan and allow the balance to perform self calibration.

## 8-3 Calibration Using the Internal mass (One-touch calibration)

This function calibrates the balance using the internal mass. The only operation required is to press the CAL key

#### Operation

- 1 Connect the AC adapter and warm up the balance for at least 30 minutes with nothing on the weighing pan.
- 2 Press the CAL key.
- 3 The balance displays I and performs calibration using the internal mass. Do not allow vibration or drafts to affect the balance.
- 4 The balance displays <u>End</u> after calibration. If the "GLP output (*unFa*)" parameter of the function table is set to "*I*" or "*2*", the balance displays <u>*GLP*</u> and outputs the "Calibration Report" using the RS-232C interface or stores the data in memory. For details on the calibration report format, refer to "11-2 GLP Report".
- 5 The balance will automatically return to the weighing mode after calibration.

#### About the internal mass

The value of the internal mass may change due to corrosion or other damage caused by the operating environment, or due to aging. Check the internal mass periodically. Correct the internal mass value as necessary. For details, refer to "8-6 Correcting the Internal Mass Value".


To maintain the weighing accuracy, perform the calibration using an external weight periodically, as described below.

## 8-4 Calibration Using an External Weight

This function calibrates the balance using an external weight.

#### When the MC-1000 is used

- 1 Connect the AC adapter and warm up the balance for at least 30 minutes with nothing on the pan.
- 2 Press and hold the CAL key until [RLout] is displayed, then release the key.
- 3 The balance displays [RL 0].
  - If you want to change the calibration weight (a list of usable weights is shown on page 20), press the SAMPLE key and proceed to Step 4.
  - If you use the calibration weight value stored in the balance, proceed to Step 5.
- 4 Specify the calibration weight value as follows:
  - SAMPLEkeyTo switch between the<br/>calibration weight selection<br/>mode (All of the digits blinking)<br/>or the value adjustment mode<br/>(The selected digits blinking).<br/>With the MC-1000, the digits in<br/>the 2<sup>nd</sup>, 3<sup>rd</sup>, and 4<sup>th</sup> decimal<br/>places can be adjusted.<br/>With the MC-6100, the digits in<br/>the 2<sup>nd</sup> and 3<sup>rd</sup> decimal places<br/>can be adjusted.
  - RE-ZERO key To select the calibration weight (All of the digits blinking) or adjust the value (The selected digits blinking). Refer to page 20
  - PRINT key To store the new weight value. Even if the AC adapter is removed, the data is maintained in non-volatile memory.
  - CAL key To cancel the operation and return to [RL 0].

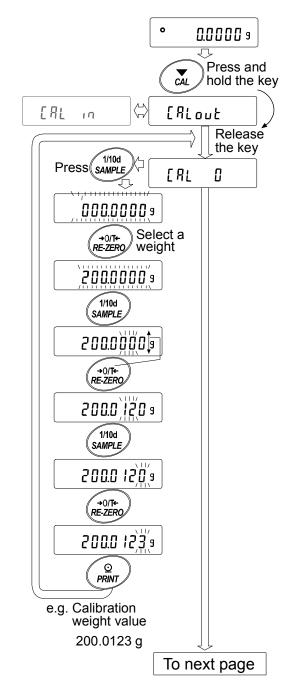


5 Confirm that there is nothing on the pan and press the PRINT key. The balance measures the zero point. Do not allow vibration or drafts to affect the balance.

The balance displays the calibration weight value.

- 6 Place the displayed calibration weight on the pan and press the PRINT key. The balance measures the calibration weight. Do not allow vibration or drafts to affect the balance.
- 7 The balance displays  $\boxed{End}$ . Remove the weight from the pan.
- 8 If the "GLP output (*inFa*)" parameter, of the function table, is set to "*I*" or "*2*", the balance displays <u>*GLP*</u> and outputs the "Calibration Report" using the RS-232C interface or stores the data in memory. For details on the calibration report format, refer to "11-2 GLP Report".
- 9 The balance will automatically return to the weighing mode.
- 10 Place the calibration weight on the pan and confirm that the value displayed is within  $\pm 2$  digits of the specified value. If it is not within the range, check the ambient conditions such as breeze and vibration, also check the weighing pan. Then, repeat Steps 1 to 10.



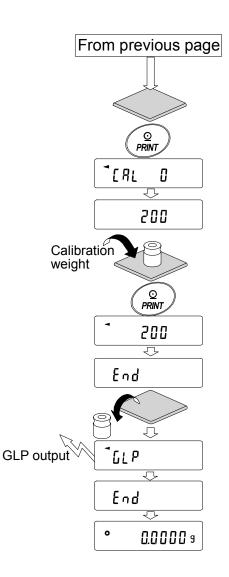


## 8-5 Calibration Test Using an External Weight

This function tests the balance weighing accuracy using an external mass and outputs the result. This is available only when the "GLP output (mFa)" parameter is set to "I" or "2". (Calibration test does not perform calibration.)

#### When the MC-1000 is used

- 1 Connect the AC adapter and warm up the balance for at least 30 minutes with nothing on the pan.
- 2 Press and hold the CAL key until [[out] is displayed, then release the key.
- 3 The balance displays [[[].
  - If you want to change the target weight (a list of usable weights is shown on page 20), press the SAMPLE key and proceed to Step 4.
  - If you use the target weight value stored in the balance, proceed to Step 5.
- 4 Specify the target weight value as follows:
  - SAMPLEkeyTo switch between the target<br/>weight selection mode (All of<br/>the digits blinking) or the<br/>value adjustment mode (The<br/>selected digits blinking).With the MC-1000, the digits in<br/>the 2<sup>nd</sup>, 3<sup>rd</sup>, and 4<sup>th</sup> decimal<br/>places can be adjusted.With the MC-6100, the digits in<br/>the 2<sup>nd</sup> and 3<sup>rd</sup> decimal places<br/>can be adjusted.
  - RE-ZERO key To select the target weight (All of the digits blinking) or adjust the value (The selected digits blinking). Refer to page 20
  - PRINT key To store the new weight value. Even if the AC adapter is removed, the data is maintained in non-volatile memory.
  - CAL keyTo cancel the operation and<br/>return to  $\begin{bmatrix} c & 0 \end{bmatrix}$ .

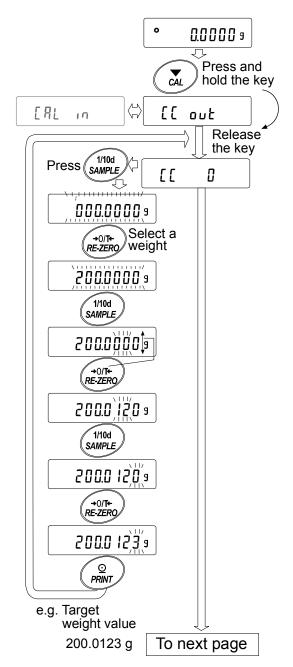

- From previous page 5 Confirm that there is nothing on the pan and press the PRINT key. The balance measures the zero point and displays the measured value. Do not allow vibration or drafts to affect the balance. The balance displays the target weight value. **1**] 0 • 0.0000 g 200 Target weight 6 Place the displayed target weight on the pan and press the PRINT key. The balance measures the target weight <u>o</u> and displays the measured value. Do not allow vibration PRINT or drafts to affect the balance. 200 6 E S1 0.005 End 7 The balance displays [End]. Remove the weight from the pan. ΓL Ρ GLP output 8 The balance displays  $\boxed{LP}$  and outputs the "Calibration" End Test Report" using the RS-232C interface or stores the  $\overline{\nabla}$ calibration test data in memory. For details on the 0 0.0000 g calibration test report format, refer to "11-2 GLP Report".
- 9 The balance will automatically return to the weighing mode.


## 8-6 Correcting the Internal Mass Value

The MC series balance can correct the internal mass value within the range shown below. This function corrects the internal mass value to conform to an external weight. The corrected mass value is maintained in non-volatile memory even if the AC adapter is removed.

Model	Correction reference value	Correction range
MC-1000	500.000 g	±0.020 g
MC-6100	2000.00 g	±0.20 g

The internal mass value is corrected as follows:

#### Example: Using the MC-6100

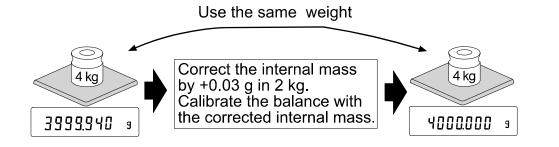


#### Operation


1 Calibrate the balance using the internal mass (one-touch calibration). Then, place an external weight and confirm the value to be corrected.

In the example, the value is off by -0.06 g and the correction reference value of the MC-6100 is 2 kg. So, the value is to be corrected by 0.03 g in 2 kg.


- 2 Press the ON:OFF key to turn off the display.
- 3 While pressing and holding the PRINT key and the SAMPLE key, press the ON:OFF key. The balance displays 75.

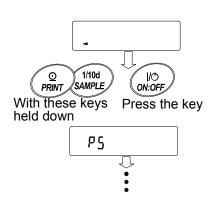

4 Press the **PRINT** key. Then the balance displays the function switches.

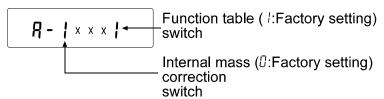

Set the function table switch and internal mass correction switch to "l" as shown above using the following keys.

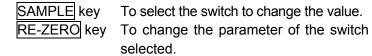


- 5 Press the PRINT key to store the new setting. The balance returns to the weighing mode.
- 6 Press and hold the <u>SAMPLE</u> key to enter the function table and release the key when <u>bR5Fnc</u> is displayed.
- 7 Press the SAMPLE key several times until [5 in] is displayed.
- 8 Press the PRINT key to enter the procedure for correcting the internal mass value.
- 9 Correct the internal mass value using the following keys.

RE-ZERO key	To select the value.					
	(-20 digits appear after +20 digits.)					
PRINT key	To store the new value and display the					
	next menu of the function table.					
CAL key	To cancel the correction and display					
	the next menu of the function table.					

- 10 Press the CAL key. The balance returns to the weighing mode.
- 11 Press the CAL key to calibrate the balance using the internal mass.
- 12 Place the external weight on the pan and confirm that the correction has been performed properly (Confirm that the value displayed is within the range that is described at "Accuracy right after calibration using the internal mass" of "18. SPECIFICATIONS".)

If the internal mass value has not been corrected properly, repeat the procedure, starting at Step 6 and adjust the correction value.

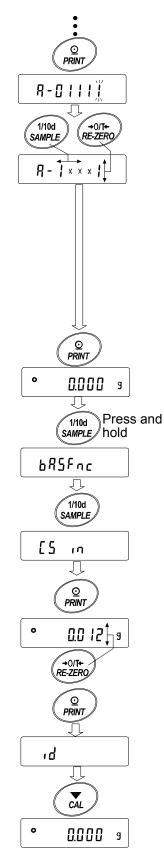


## 9. FUNCTION SWITCH AND INITIALIZATION

### 9-1 Permit or Inhibit

The balance stores parameters that must not be changed carelessly (e.g. Calibration data for accurate weighing, Data for adapting to the operating environment, Control data for the RS-232C interface). There are five switches for the purpose of protecting these parameters. Each switch can select either "permit" or "inhibit". "Inhibit" protects parameters against careless operations.

#### Switches

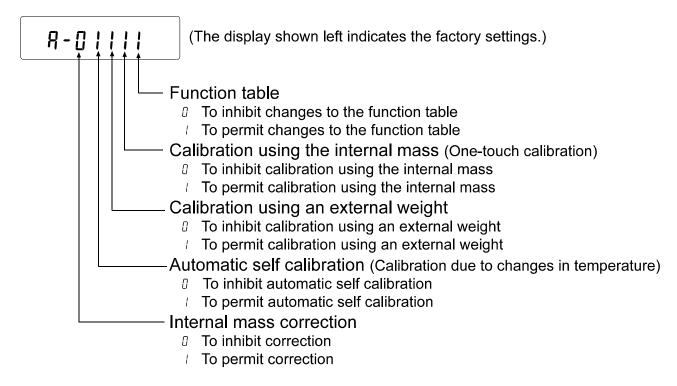

#### Operation

- 1 Press the ON:OFF key to turn off the display.
- 2 While pressing and holding the PRINT key and the SAMPLE key, press the ON:OFF key. The balance displays 2.
- 3 Press the PRINT key. Then the balance displays the function switches.
- 4 Set the switches using the following keys.

SAMPLE key	To select the switch to change the parameter.
RE-ZERO key	To change the parameter of the switch selected.

D: To inhibit changes. I: To permit changes

- PRINT keyTo store the new parameter and return to the weighing mode.
- CAL key To cancel the operation. ( [[] is displayed.)

To return to the weighing mode, press CAL key once again.

## 9-2 Initializing the Balance

This function returns the following parameters to factory settings.


- Calibration data
- Function table
- The sample unit mass value (counting mode), 100% reference mass value (percent mode)
- The data that is stored in the balance using the data memory function
- External calibration weight and target weight value
- Function switch settings

#### Note

Be sure to calibrate the balance after initialization.

#### Operation

- 1 Press the ON:OFF key to turn off the display.
- 2 While pressing and holding the PRINT key and the SAMPLE key, press the ON:OFF key. The balance displays P5.

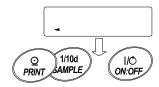

With these keys Press the key held down

PS

LI r (L r)))) (L r (L r (L r))) (L r (L r)) (L r (L r)) (L r) (

0.0000 s

- 3 Press the SAMPLE key to display [[1].
- 4 Press the PRINT key. To cancel this operation, press the CAL key.
- 5 Press the RE-ZERO key.
- 6 Press the PRINT key to initialize the balance.

The balance will automatically return to the weighing mode.

## **10. FUNCTION TABLE**


The function table reads or rewrites the parameters that are stored in the balance. These parameters are stored in non-volatile memory, and are maintained even if the AC adapter is removed.

## **10-1 Structure and Sequence of the Function Table**

The function table menu consists of two layers. The first layer is the "Class" and the second layer is the "Item". Each item stores a parameter.

#### Example

This example sets "Stores weighing data" for "Data memory" and "Every 1 minute" for "Interval time".

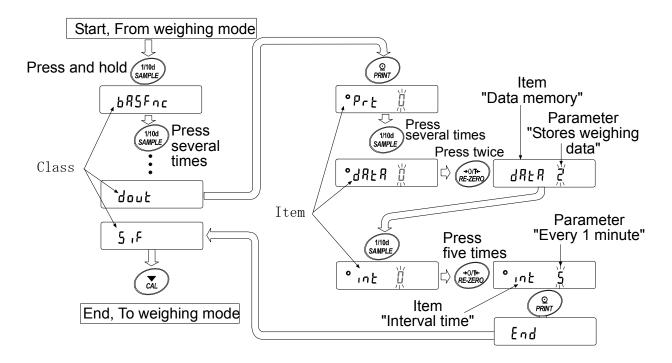

## 10-2 Display and Keys

Display/Key	Description
•	The symbol " <b>O</b> " indicates that the parameter displayed is in effect.
1/10d	When pressed and held in the weighing mode, enters the function table mode.
SAMPLE	Selects the class or item in the function table mode.
+0/JT+ RE-ZERO	Changes the parameter.
0	When a class is displayed, moves to an item in the class.
	When an item is displayed, stores the new parameter and displays the next class.
	When an item is displayed, cancels the new parameter and displays the next class.
( CAL	When a class is displayed, exits the function table mode and returns to the weighing
)	mode.

## **10-3 Details of the Function Table**

Class	Item and Paramete	er	Descri	ption			
	E and Condition	0 1 • 2	Fast response rate, sensitive value       FAST         MID.         Slow response rate, stable value       SLOW	Can be changed by response adjustment. With "hoLd I", sets the averaging time.			
	5と-6 Stability band width		Stable range is ±1 digit	The stabilization indicator illuminates when the display fluctuation is within the range. With "hoLd I",			
		2	Stable range is ±3 digits	sets the averaging time. Holds the display when stable			
	Hald Hold function	= 0 	OFF ON	in animal mode. With "Hold I", ANIMAL turns on.			
		0	OFF				
	trc Zana tanakina	<b>•</b> /	Normal	Keeps zero display by tracking			
	Zero tracking	2	Strong	zero drift.			
		3	Very strong				
	5Pd Diaplay refresh rate	• 0	5 times/second	Period to refresh the display			
	Display refresh rate	1	10 times/second				
bASFnc	PnE Decimal point	• 0	Point (.)	Decimal point format			
Environment Display	nt Decimal point	1	Comma (,)	Turns on the weighing mode			
Auto display-Of PoFF Auto display-Of		• 0	OFF	display when the AC adapter			
		1	ON	is plugged in.			
		• 0	OFF	Turns off the display after 10			
			ON (10 minutes)	minutes of inactivity.			
	65 , Capacity indicator	• 0	OFF	Capacity indicator Zero: 0%.			
			ON	Maximum capacity: 100%			
	FIL	• 0	Not used				
	Filter		Used (when the balance is used as a mass comparator)				
	P-tr	• 0	Sets the display to zero.				
	Zero upon power-on	1	Does not set the display to zero. Displays the previous value. With this setting, do not perform re-zero operations frequently. Refer to "Zero upon power-on" on page 35.				
		• 0	Performs zero and span calibration	on			
	5Pn Span calibration /		Performs span calibration only. Span calibration using the internal mass is possible with a tare on the weighing pan. Refer to "Span calibration" on page 35.				
EL AdJ Clock			to the GX series instruction al, "10-9 Clock and Calendar	Confirms and sets the time and date. The time and date are added to the output data.			
		- 0	No comparison	·			
	50	1	Comparison, excluding "near zero	o" when stable or overloaded			
	[P Comparator mode	2	Comparison, including "near zero	" when stable or overloaded			
		Э	Continuous comparison, excludin				
[P Fnc		Ч	Continuous comparison, including	g "near zero"			
Comparator		- 0	Digital input, upper/lower limits				
	EP in	1	Weighing input, upper/lower limits	[P H I, [P Lo can be selected.			
	Input method	2	Digital input, reference value	[P rEF, [P LnE can be			
		3	Weighing input, reference value	selected			
			torv setting				

Factory setting Note: "Digit" is a unit of minimum weighing value

Class	Item and Parameter		Description		
Upper limit				Displays when [P in ] or [P in ] is selected.	
<i>EP Lo</i> Lower limit			Refer to the GX series	Lr in TIS Selected.	
LOWER MIN			instruction manual, "10-10		
Reference val	lue		Comparator Function".	Displays when [P in 2 or	
[P Lit				[P in ] is selected.	
Tolerance					
		• 0	Key mode (when stable)	Accepts the PRINT key only when the display is stable.	
		1	Auto print mode A	Outputs data when the display is	
	Prt		(Reference = zero) Auto print mode B	stable and conditions of $RP-P$ , RP-b and the reference value are	
	Data output mode	2	(Reference = last stable value)	met.	
				With dALA D, outputs data	
		Э	Stream mode / Interval memory mode	continuously; with dALA 2, uses	
				interval memory.	
	RP-P	• 0	Plus only	Displayed value>Reference	
	Auto print polarity	1	Minus only	Displayed value <reference< td=""></reference<>	
		2	Both	Regardless of displayed value	
	<i>ЯР-Ь</i> Auto print difference	- 0	10 digits		
		1	100 digits	Difference between reference value and displayed value	
		2	1000 digits		
	dЯLЯ Data memory	• 0	Not used		
		1	Stores unit mass in counting mode	Related items:	
		2	Stores weighing data	Prt, int, d-no, 5-td, info	
		Э	Stores calibration data		
dout	וחנ Interval time	- 0	Every measurement		
Data output		1	2 seconds	-	
		2	5 seconds		
		Э	10 seconds	Interval time in the interval	
		Ч	30 seconds	memory mode when using Prt 3,	
		5	1 minute	AHFH S	
		6	2 minutes	-	
		٦	5 minutes		
		8	10 minutes		
	d-no	- 0	No output	Refer to the GX series	
	Data number output	1	Output	instruction manual, "12. DATA MEMORY".	
		• 0	No output	Selects whether or not the time	
	5-Ed		Time only	or date is added to the weighting	
	Time/Date output	2	Date only	data. For details, Refer to the GX series instruction manual,	
		Э	Time and date	"10-9 Clock and Calendar Function".	
	5- id	• 0	No output	Selects whether or not the ID	
	ID number output	1	Output	number is output.	
	PUSE	• 0	No pause		
	Data output pause		Pause (1.6 seconds)	Selects the data output interval.	
	1 .	Eac	tory setting		

Factory setting Note: "Digit" is a unit of minimum weighing value

Class	Item and Parame	eter	Description		
	RE-F	- 0	Not used	Selects whether or not automatic	
	Auto feed	- 1	Used	feed is performed.	
<b>לביט</b> Data output		- 0	No output	Selects GLP output method.	
	inFa	- 1	AD-8121 format	For how to set time and date to be added, refer to the GX series	
	GLP output	2	General data format	instruction manual, "10-9 Clock and Calendar Function".	
	Rr-d	- 0	Not used	Adjusts zero automatically after	
	Zero after output	1	Used	data is output.	
		0	600 bps		
	ьРБ	- 1	1200 bps		
	Baud rate	- 2	2400 bps		
	Bada rato	Э	4800 bps		
		Ч	9600 bps		
	<i>ЬЕРг</i> Data bit, parity bit	- 0	7 bits, even		
		- 1	7 bits, odd		
		2	8 bits, none		
<b>.</b> .	[rLF	- 0	CR LF	CR: ASCII code 0Dh	
	Terminator	- 1	CR	LF: ASCII code 0Ah	
5 ,F Serial	<i>Ł ЧРЕ</i> Data format	- 0	A&D standard format		
interface		- 1	DP format		
		2	KF format	Refer to the GX series instruction	
		3	MT format	manual, "10-6 Description of Item "Data Format".	
		Ч	NU format		
		5	CSV format		
	<i>Е - UP</i> Timeout	0	No limit	Selects the wait time to receive a command.	
		- /	1 second		
	Er[d	- 0	No output	AK: ASCII code 06h	
	AK, Error code	1	Output		
	[LS	- 0	Not used	Controls CTS and RTS.	
	CTS, RTS control	1	Used		
ป่า เŁ Unit		g	gram	Refer to "5. WEIGHING UNITS".	
d5 Fnc Density	Ldin	- 0	Water temperature	Available only when density mode is selected. Refer to the GX series	
function	Liquid density input	1	Liquid density	instruction manual, "15. DENSITY MEASUREMENT".	
תי 5 Internal mass value correction		Refer to "8. CALIBRATION".	Displayed only when the internal mass value correction switch is set to 1.		
ر <b>ا</b> ID number set	ting		Refer to "11. ID NUMBER AND	) GLP REPORT".	
<ul> <li>Factory setting</li> </ul>					

Factory setting

Class	Item and Parameter		Description
$E_r F_{nc}$ Extended function Only when $F_{l}L$ is $D$	F I-b Averaging range for the first moving average	0	Small
		1	
		2	
		Э	
		<b>-</b> 4	
		5	
		6	
		٦	▼ Large
	F I-Ł Averaging time for the first moving average	0	No averaging
		1	0.5 second
		• 2	1.0 second
		3	1.5 seconds
		Ч	2.0 seconds
		5	2.5 seconds
		6	3.2 seconds
		٦	4.8 seconds
		8	6.4 seconds
	F2-b Averaging range for the second moving average	0	Small
		- /	
		2	
		3	
		Ч	
		5	$\checkmark$
		6	Large
	F2-Ł Averaging time for the second moving average	0	No averaging
			0.5 second
		2	1.0 second
		Э	1.5 seconds
		Ч	2.0 seconds
		<b>■</b> 5	2.5 seconds
		6	3.2 seconds
<ul> <li>Factory setting</li> </ul>			

#### Caution

The balance may not transmit the data completely at the specified refresh rate, depending on the baud rate or data added to the weighing data such as time, date and ID number.

## 10-4 Description of the Class "Environment, Display"

Condition ( [and )	
Cond D	This parameter is for sensitive response to the fluctuation of a weight value.
$\bigcup$	Used for powder target weighing, weighing a very light sample or when quick response weighing is required. After setting, the balance displays FAST.
Cond 2	This parameter is for stable weighing with slow response. Used to prevent a weight value from drifting due to vibration or drafts. After setting, the balance displays SLOW.
	Notes
	In automatic response adjustment, the response rate is selected

automatically.
 With "Hold function (HoLd)" set to "ON (1)", this item is used to set the averaging time.

### Stability band width (5L-b)

This item controls the width to regard a weight value as a stable value. When the fluctuation per second is less than the parameter, the balance displays the stabilization indicator and outputs or stores the data. The parameter influences the "Auto print mode"



This parameter is for sensitive response of the stabilization indicator. Used for exact weighing.

This parameter ignores slight fluctuation of a weight value. Used to prevent a weight value from drifting due to vibration or drafts.

#### Note

With "Hold function (  $H_{oLd}$  )" set to "ON ( l )", this item is used to set the stabilization range.

### Zero upon power-on (P-tr)

When a hopper is attached to the weighing pan and loss-in weighing is performed, the remaining amount of the material will become unknown if tare is performed each time a weighing starts.

When "P-tr " is set to " l ", tare is not performed at weighing start. So, the remaining amount of the material can be monitored, when the power is turned on again after it was turned off.

### Span calibration (5Pn)

When a hopper is attached to the weighing pan and calibration is to be performed with the hopper attached, set "  $5P_n$  " to " 1". When the tare value (hopper and other devices attached) is within the value in the table below, calibration using the internal mass is possible.

Model	Tare value
MC-1000	500 g or less
MC-6100	5 kg or less

## **10-5** Description of the Item "Data output mode"

The parameter setting of the "Data output mode ( $P_r E$ )" applies to the performance when the "Data memory (dR ER)" parameter is set to "2" (to store the weighing data)and when the data is transmitted using the RS-232C interface.

### Key mode

When the **PRINT** key is pressed with the stabilization indictor turned on, the balance outputs or stores the weighing data and the display blinks one time.

Required setting dout Prt 0 Key mode

### Auto print modes A and B

When the displayed value is stable and the conditions of "Auto print polarity", "Auto print difference" and reference value are met, the balance outputs or stores the weighing data.

When the **PRINT** key is pressed with the stabilization indictor turned on, the balance outputs or stores the data and the display blinks one time.

Mode A: Example	•	"For weighing each time a sample is added (or removed), with "Ar - d" set to " /" (to adjust zero after the data is output)."					
Required setting	dout dout dout	РгЕ   АР-Р АР-Ь	Auto print mode A (reference = zero) Auto print polarity Auto print difference				
Mode B: Example	"For weigh	ning while a sar	nple is added."				
Required setting	dout	Prt 2	Auto print mode B (reference = last stable value)				
	dout dout	ЯР-Р ЯР-Ь	Auto print polarity Auto print difference				

### Stream mode

The balance outputs the weighing data continuously regardless of the display condition. The display does not blink in this mode. This mode is not available and the interval memory mode is used when the "Data memory (dRER)" parameter is set to "2" (to store the weighing data).

Example	"For monitoring data on a computer"						
Required setting	dout	Prt 3	Stream mode / Interval memory mode				
	dout	dAtA 0	Data memory function is not used.				
	bRSFnc	SPd	Display refresh rate				
	5 iF	bPS	Baud rate				

### Caution

The balance may not transmit the data completely at the specified refresh rate, depending on the baud rate or data added to the weighing data such as time, date and ID number.

### Interval memory mode

The weighing data is periodically stored in memory.

Example	"For perio	odical weighing	without a computer command and
	outputting	all of the data to	a computer at one time"
Required setting	dout dout	Prt 3 dAtA 2	Stream mode / Interval memory mode Data memory function is used. Stores weighing data.
Optional setting	dout dout	int 5-td I, 2, or 3	Interval time Adds the time and date.

## **10-6** Description of the Item "Data format"

### A&D standard format 5 IF LYPE 0

This format is used when the peripheral equipment can receive the A&D format. If an AD-8121B is used, set the printer to MODE 1 or 2.

- This format consists of fifteen or sixteen characters excluding the terminator. With the MC-1000, when the numerical data excluding the decimal point exceeds eight characters, the data format will be sixteen characters long.
- A header of two characters indicates the balance condition.
- The polarity sign is placed before the data with the leading zeros. If the data is zero, the plus sign is used.
- The unit, consisting of three characters, follows the data.

ST	, + 0 0 0 . 0	1 2 7 🗆 🗤 g	CR LF
Heade	er Data	Unit	Terminator
ST	, + 1 0 0 0 .		$g C_R L_F$ (Data example exceeding 8 characters)
Heade	er Data	Unit	Terminator
•			
ST	Stable header	Q T Stable hea	der (Counting mode)
US	Unstable header		
ΟL	Overload header		

### DP (Dump print) format 5 IF LYPE 1

This format is used when the peripheral equipment can not receive the A&D format. If an AD-8121B is used, set the printer to MODE 3.


- This format consists of sixteen characters excluding the terminator.
- A header of two characters indicates the balance condition. No overload header is used.
- The polarity sign is placed before the data, with spaces in place of leading zeros, if the data is not zero or overloaded.
- The unit, consisting of three characters, follows the data.

W T + 1 0 0	0.000		CRLF
Header	Data	Unit	Terminator
▼ W T Stable heade U S Unstable hea		table head	er (Counting mode)

### KF format 5 IF ESPE 2

This is the Karl-Fischer moisture meter format and is used when the peripheral equipment can only communicate using this format.

- This format consists of fourteen characters excluding the terminator.
- This format has no header characters.
- The polarity sign is placed before the data, with spaces in place of leading zeros, if the data is not zero or overloaded.
- This format outputs the unit only for a stable value.

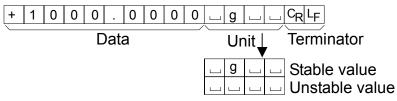

### MT format 5 ,F LYPE 3


- A header of two characters indicates the balance condition.
- The polarity sign is used only for negative data.
- The weighing data uses spaces in place of the leading zeros.
- The character length of this format changes dependent upon the unit

S 1 0	00.00	000gC <sub>R</sub> L <sub>F</sub>
Header	Data	Unit Terminator
	header le header ad header	

### NU (Numerical) format 5 .F LYPE 4

This format outputs only numerical data.

- This format consists of nine or ten characters excluding the terminator. With the MC-1000, when the numerical data excluding the decimal point exceeds eight characters, the data format will be ten characters long.
- The polarity sign is placed before the data with the leading zeros. If the data is zero, the plus sign is used.

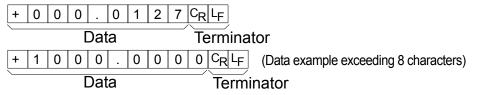

### CSV format 5 IF LYPE 5

- Separates the data of A&D standard format and the unit by a comma (,).
- Outputs the unit even when the data is overloaded.
- When the ID number, data number, time and date are added, outputs the ID number, data number, date, time and weighing data in this order and separates each item by a comma and treats all the items as one group of data.

LAB-123, No,012, 2009/12/31, 12:34:56, ST,+1000.0000, \_\_ g<CR><LF>

ID number Data number Date Time Weighing data

S	Т	,	+	1	0	0	0		0	0	0	0	,			g	$C_R$	LF	
0	L	,	+	9	9	9	9	9	9	9	Е	+	1	9	,			g	C <sub>R</sub> L <sub>F</sub>

### Note

To add the ID number, data number, time and date, the function settings must be changed.

## **10-7** Description of the Data Format Added to the Weighing Data

ID number dout 5-id 1

The number to identify a specific balance.

• This format consists of seven characters excluding the terminator.

L	А	В	-	1	2	3	$C_R$	$L_{F}$
---	---	---	---	---	---	---	-------	---------

### Data number dout dono l

This format outputs the data number just before the data is transmitted using the RS-232C interface.

- This format consists of six characters excluding the terminator.
- When CSV format (5, F LYPE 5) is selected, the period (.) is replaced with a comma (, ).

N 0 . 0 0 1 C<sub>R</sub> L<sub>F</sub>


Data number Terminator

### Note

The data number is added only to the weighing data that is stored in memory.

**Date** dout 5-td 2 or 3

• The date output order can be changed in "Clock ([L RdJ)". Outputs the year in four-digit format.

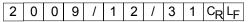


Time dout 5-td / or 3

- Outputs time in 24-hour format.
  - $\begin{array}{|c|c|c|c|c|c|c|c|}\hline 1 & 2 & \vdots & 3 & 4 & \vdots & 5 & 5 & C_R & L_F \\ \hline \end{array}$

### Note

When the data described above is added to the weighing data, the output is in the following order: ID number, Data number, Date, Time and Weighing data.

### **10-8 Data Format Examples**

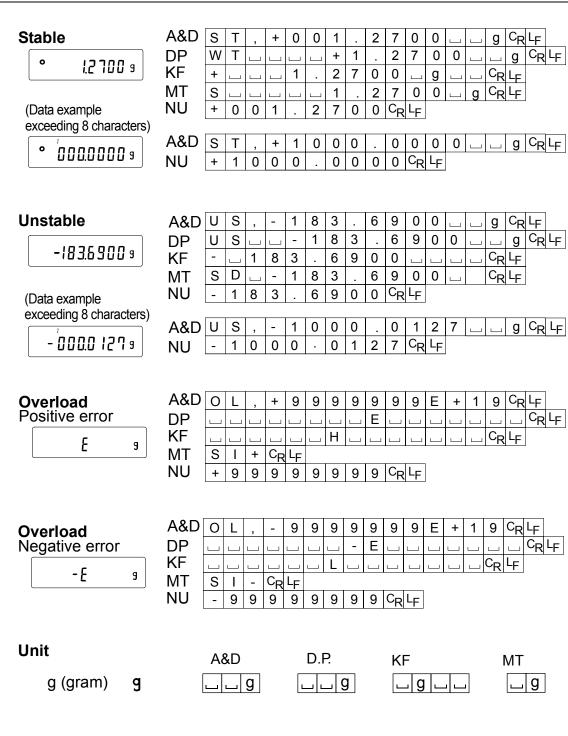

- □ Space, ASCII 20h
- <sup>C</sup><sub>R</sub> Carriage Return, ASCII 0Dh
- <sup>L</sup>F Line Feed, ASCII 0Ah

## **11. ID NUMBER AND GLP REPORT**

- The ID number is used to identify the balance when Good Laboratory Practice (GLP) is used.
- The ID number is maintained in non-volatile memory even if the AC adapter is removed.
- The GLP output format is selected at "GLP output ( *mFa*)" of the function table and can be output to a personal computer or printer using the RS-232C serial interface.
- The GLP output format includes the balance manufacturer, model, serial number, ID number, date, time and space for signature for weighing data, and the weight used and results for calibration or calibration test data.
- The balance can output the following reports for GLP.

"Calibration report" of the calibration, using the internal mass (Calibration due to changes in temperature and one-touch calibration.)

"Calibration report" of the calibration, using an external weight.

"Calibration test report" of the calibration test, using an external weight.

"Title block" and "End block" for the weighing data.

- Calibration and calibration test data can be stored in memory to output several reports at the same time. Refer to the GX series instruction manual, "12. DATA MEMORY" for details.
- For details on confirming and setting the time and date, refer to the GX series instruction manual, "10-9 Clock and Calendar Function".

## 11-1 Setting the ID Number

1	Press and hold the	SAMPLE key until	685Fnc	of the function table is displayed.
---	--------------------	------------------	--------	-------------------------------------

- 2 Press the SAMPLE key several times to display \_\_\_\_\_.
- 3 Press the PRINT key. Set the ID number using the following keys.

RE-ZERO key To set the character of the digit selected. Refer to the display character set shown below.

SAMPLE key To select the digit to change the value.

PRINT key To store the new ID number and display bRSFac

CAL key To cancel the new ID number and display bR5Fnc.

4 With bRSFnc displayed, press the CAL key to return to the weighing mode.

### **Display character set**



## 11-2 GLP Report

Set the following parameters to output the report.

- To print the report, set the "GLP output ( InFa )" parameter to "1", the "Data output pause (*PU5E*)" parameter to "1", and use MODE 3 of the AD-8121B. For details on using the printer, refer to "13-1 Connection to the AD-8121B Printer".
- To output the report to a personal computer using the RS-232C interface, set the "GLP output ( unFa )" parameter to " 2 ".
- If the time and date are not correct, set the correct time and date in "Clock (*LL* RdJ)" of the function table.

### Calibration report using the internal mass

When the setting is "inFo I":

When the setting is "InFo 2":

AD-8121 format

General data format

A & D MODEL MC-1000 S/N 01234567 ID ABCDEF6 DATE 2009/12/31 TIME 12:34:56 CALIBRATED(INT.) SIGNATURE	← Manufacturer → ← Model → ← Serial number → ← ID number → ← Date → ← Calibration type ← Signature	A.&.D <term> MODELMC-1000<term> S/N01234567<term> IDABCDEFG<term> DATE<term> 2009/12/31<term> TIME<term> CALIBRATED (INT.)<term> SIGNATURE<term> <term> <term></term></term></term></term></term></term></term></term></term></term></term>
Space, ASCII 2	<term> <term></term></term>	

Space, ASCII 20h
 <TERM> Terminator, C<sub>R</sub>, L<sub>F</sub> or C<sub>R</sub>
 <sup>C</sup>R Carriage return, ASCII 0Dh
 <sup>L</sup>F Line feed, ASCII 0Ah


### Calibration report using an external weight

When the setting is " I ":

### AD-8121 format

When the setting is " InFo 2 ":

General data format
→ \_\_\_\_\_A\_&\_D<TE

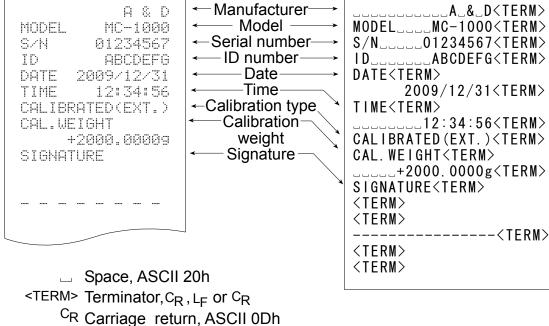

### Calibration test report using an external weight


LF Line feed, ASCII 0Ah

(Calibration test does not perform calibration.) When the setting is " a = l ":

AD-8121 format

When the setting is " In Fo 2 ":

General data format

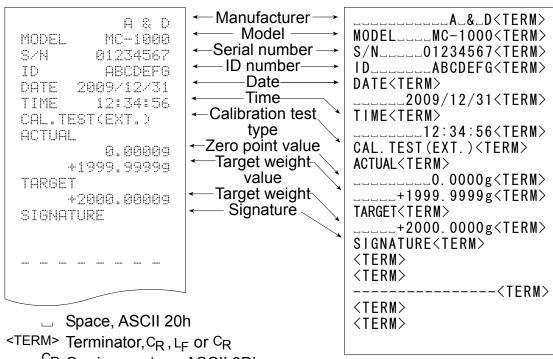

LF Line feed, ASCII 0Ah

### Title block and end block

When a weight value is recorded as the GLP data, "Title block" and "End block" are inserted at the beginning and at the end of a group of weight values, in the GLP report.

### Notes


- □ To output the report to an AD-8121B , use MODE 3 of the AD-8121B.
- □ If the data memory function is used, the "Title block" and "End block" can not be output.

### Operation

- 1 With the weighing data displayed, press and hold the PRINT key until 5tRrt is displayed. The "Title block" is output.
- 2 The weighing data is output according to the parameter setting of the data output mode.
- 3 Press and hold the PRINT key until  $\boxed{r \mathcal{E} \mathcal{E} \mathcal{E} \mathcal{I} d}$  is displayed. The "End block" is output.

When the setting is " I ":


When the setting is " In Fa 2 ":

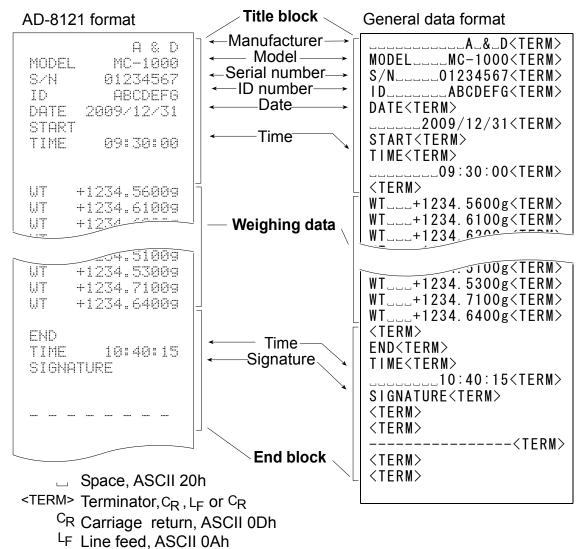

## 12. RS-232C SERIAL INTERFACE / EXTERNAL INPUT

### RS-232C

The balance is a DCE device. Connect the balance to a personal computer (DTE), using a straight through cable or a modem cable.

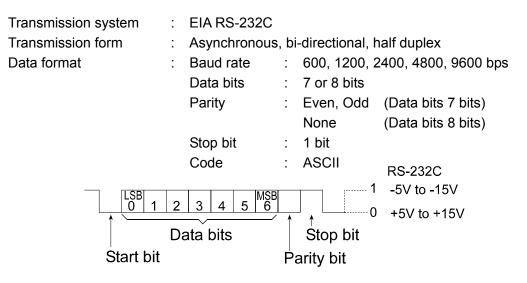


### **Pin connections**


Pin No.	Signal name	Direction	Description
1	FG	-	Frame ground
2	RXD	Input	Receive data
3	TXD	Output	Transmit data
4	RTS	Input	Ready to send
5	CTS	Output	Clear to send
6	DSR	Output	Data set ready
7	SG	-	Signal ground
18	PRINT	Input	Same as the PRINT key
19	RE-ZERO	Input	Same as the RE-ZERO key
8 – 25 (Excluding 18 and 19)	-	-	Not connected

### Note

The signal names for the RS-232C are the same as those for the DTE side except TXD and RXD.

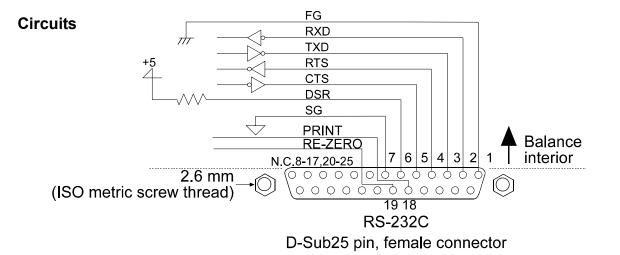


### **RS-232C** Terminals

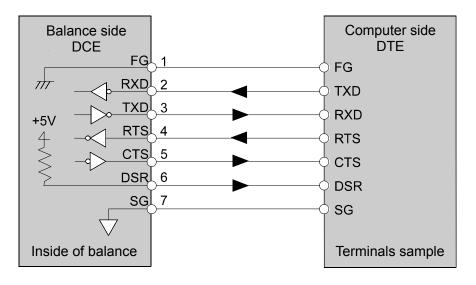

### **External contact input**


By connecting pin 18 (PRINT command) to pin 7, or pin 19 (RE-ZERO command) to pin 7 for 100 ms or more, the same operation as performed by pressing the PRINT key or the RE-ZERO key, will be performed.

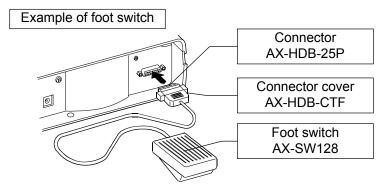

### Option

Connector

÷

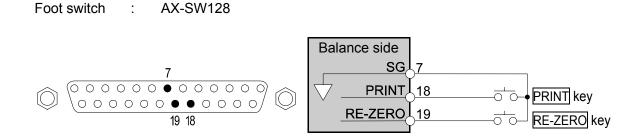

AX-HDB-25P/CTF

## **13. CONNECTION TO PERIPHERAL EQUIPMENT**

## 13-1 Connection to the AD-8121B Printer

Class Item		Factory			
Class		setting	Mode 1	Mode 2	Mode 3
	PrE Data output mode	0	0, 1,2,4,5	Э	0, <i>1</i> , 2, 4, 5
	<i>RP-P</i> Auto print polarity	0	*1	Not	*4
	<i>RP-b</i> Auto print difference	0	~1	applicable	*1
dout	d-חם Data number output	0	0	0	0, I
Data output	Time/date output	0	0	0	0, I, 2, 3
	5- ،d ID number output	0	0	۵	0, I
	PUSE Data output pause	0	0	0	0,1 * <b>2</b>
	<i>RŁ - F</i> Auto feed	0	0	0	0, I
	bР5 Baud rate	2	2	2	2
5 <i>،F</i> Serial interface	PEPr Data bit, parity bit	0	0	0	0
	[rLF Terminator	0	0	0	0
	<i>EYPE</i> Data format	0	0	۵	1
	[£5 CTS, RTS control	0	0	0	0

Set the following parameters to use the AD-8121B printer.

Set the parameter when the data output mode is set to the auto print mode (Prt I or Prt 2). \*1

\*2 When multiple lines are to be printed, set the parameter to 1.

### AD-8121B settings

Mode	AD-8121B DIP switches	Description
Mode 1		Prints upon data receipt Standard mode, statistical calculation mode
Mode 2		Prints using the AD-8121B DATA key or the AD-8121B built-in timer Standard mode, interval mode, chart mode
Mode 3		Prints upon data receipt Dump print mode

DIP switch 3: Handling unstable data

MODE 2 ON = To print unstable data

OFF = Not to print unstable data

DIP switch 4: Data input specification Q1534 ON = Use the current loop

OFF = Use the RS-232C



 $\square$  With dRLR D, the weight data can be output to the printer.

□ With dALA 2, the weight data stored in memory can be output to the printer.

 $\Box$  With dRLR 3, the calibration report stored in memory can be output to the printer.

□ Refer to "11. ID NUMBER AND GLP REPORT" for print samples.



## 13-2 Connection to a Computer

The MC series balance is of the DCE type (Data Communication Equipment), which can be connected to a personal computer using the RS-232C interface.

Before connection, read the personal computer manual thoroughly.

Use a standard DCE cable for connection (cable type: straight-through).

## 13-3 Using Windows Communication Tools (WinCT)

When Windows is used as an operating system in a personal computer, the provided WinCT can be used to transmit the weighing data to the personal computer.

The WinCT has two communication methods: "RsCom" and "RsKey". For details on WinCT, refer to the WinCT instruction manual.

The current version of the WinCT can be downloaded from the A&D website.

### RsCom

- Can transmit commands to control the balance.
- Can perform bi-directional communication between the balance and a personal computer using the RS-232C interface.
- Can display or store the data using a text file format. Can also print the data using a printer connected to the personal computer.
- When several ports of a personal computer have balances connected, the personal computer can communicate with each balance simultaneously.
- Can share a personal computer with other application software.
- Can receive the balance GLP report.

### RsKey

- Can transmit the weighing data output from the balance directly to other application software such as Microsoft Excel.
- Can be used with most application software.
- Can receive the balance GLP report.

### Note

### Windows and Excel are the registered trademarks of the Microsoft Corporation.

### Using the WinCT software, the balance can do the following:

1 Analyzing the weighing data and the statistics input by "RsKey"

The weighing data can be input directly into an Excel worksheet. Then, Excel can analyze the data to obtain total, average, standard deviation, maximum and minimum value, and display them in a graph.

2 Controlling the balance using commands from a personal computer By using "RsCom", the personal computer sends commands such as "re-zero" or "send weighing data" to the balance and controls the balance.

- Printing the balance GLP report using your printerThe balance GLP report can be printed using a printer connected to the personal computer.
- 4 Receiving weighing data at a certain interval The weighing data can be received at a certain interval and data characteristic with elapsed time can be obtained.
- 5 Using the MC series balance memory function The weighing data, instead of outputting it immediately to an external device, can be stored in the balance's memory. Later, all of the weighing data stored can be output to a personal computer at one time.
- 6 Using a personal computer as an external indicator With the "RsKey" test mode function, a personal computer can be used as an external weight indicator for the balance. (To do this, set the balance data output mode to stream mode.)

## 14. COMMANDS

### 14-1 Command List

### Commands to query weighing data

С	Cancels the <b>S</b> or <b>SIR</b> command.
Q	Requests the weighing data immediately.
S	Requests the weighing data when stabilized.
SI	Requests the weighing data immediately.
SIR	Requests the weighing data continuously.
<sup>E</sup> s <sub>C</sub> P	Requests the weighing data when stabilized.

Note: The "Q" and "SI" commands, the "S" and "<sup>E</sup>s<sub>c</sub>P" commands behave the same.

### Commands to control the balance

Note: The "R", "T", "Z" and " ${}^{E}s_{C}T$ " commands behave the same.

### Commands to control the comparator function

HI:***. ** ∟⊔ g	Sets the upper limit value. e.g.,HI:+2.34 uug (to set the upper limit	
_	value to 2.34 g:   represents a space.)	
LO:***. ** பப g	Sets the lower limit value. e.g.,LO:+1.23 g (to set the lower limit	
	value to 1.23 g: - represents a space.)	
?HI	Outputs the upper limit value.	
?L0	Outputs the lower limit value.	
lote: To use a command to control the comparator function, set the "Input meth		

Note: To use a command to control the comparator function, set the "Input method ([P in )" parameter to "[] " or " | ".

#### Commands to control the memory function

MCL	Deletes all data in memory.
MD:nnn	Deletes weighing data with the data number nnn.
?MA	Outputs all data in memory.
?MQnnn	Outputs weighing data with the data number nnn.
?MX	Outputs the number of weighing data in memory.

Note: "nnn" indicates a three-digit numerical value.

<sup>E</sup>s<sub>C</sub>: 1Bh in ASCII code

## 14-2 Acknowledge Code and Error Codes

When the "Serial interface function (5 F)" parameter is set to "ErEd I", the balance outputs <AK> code or error code to each command as follows:

<AK> (06h) Acknowledge in ASCII code.

- When the balance receives a command to request data and can not process it, the balance transmits an error code (EC, Exx). When the balance receives a command to request data and can process it, the balance outputs the data.
- When the balance receives a command to control the balance and can not process it, the balance transmits an error code (EC, Exx). When the balance receives a command to control the balance and can process it, the balance transmits the acknowledge code.

Among commands to control the balance, the following transmit the acknowledge code both when the balance receives the command and when the balance has accomplished the command. If the command can not be processed properly, the balance transmits an error code (EC, Exx). This error can be released using the CAL command.


CAL command (Calibration command)

ON command (Display ON command)

P command (Display ON/OFF command)

R command (RE-ZERO command)

### R command

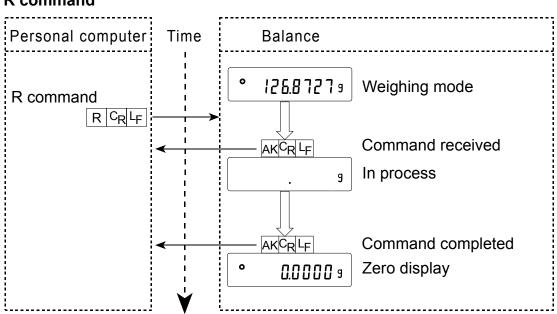

 When a communication error has occurred due to external noise, or a parity error has occurred due to transmission error, the balance transmits an error code. In this case, send the command again.

## 14-3 Control Using CTS and RTS

Depending on the "[L5" parameter of "Serial interface (5, F)", the balance performs as follows:

### CES 0

Regardless of whether the balance can receive a command or not, the balance keeps the CTS line HI. The balance outputs data regardless of condition of the RTS line.

### [25]

The CTS line is kept Hi normally. When the balance can not receive the next command (e.g. while the balance is processing last command), the balance sets the CTS line to Lo. The balance confirms the level of the RTS line before outputting a set of data. If the RTS level is Hi, the balance outputs data. If the RTS level is Lo, data is not output (The data is canceled).

## 14-4 Settings Related to RS-232C

Concerning the RS-232C, the balance has two functions: "Data output (  $d_{out}$  )" and "Serial interface (5 rF )". Set each function as necessary.

## **15. EXTENDED FUNCTION**

The MC series balance has several extended functions equipped for special applications or to troubleshoot when using the standard functions.

When the "Filter (F IL)" is set to "D", extra items are available as shown below.


Class	Item and Paramete	er	Description	
	F I-b Averaging range for the first moving average	0 1 2 3 • 4 5 6 7	Small Large	When the fluctuation of a weight value is within the averaging range, the averaging operation starts to stabilize the value. When the fluctuation is small, for example, when weighing or filling a small amount of sample, the averaging operation is always performed and the response rate may be slow. Under such a condition, change the parameter. Refer to "Averaging range ( $F I-b$ ) and averaging time ( $F I-b$ )"
F I - E Averaging time for the first moving average $Er Fnc$ Extended function $F2 - b$ Averaging range for the second moving average $F2 - E$ Averaging time for the second moving average	Averaging time for the	0 1 3 3 4 5 6 7 8	No averaging0.5 second1.0 second1.5 seconds2.0 seconds2.5 seconds3.2 seconds4.8 seconds6.4 seconds	When the fluctuation of a weight value is within the averaging range, the averaging operation starts. Once the duration of averaging operations reaches the averaging time, moving averaging starts. This parameter sets the time to start moving averaging. Refer to "Averaging range ( $FI-b$ ) and averaging time ( $FI-b$ )"
	Averaging range for the second moving	0 - 1 - 2 - 3 - 4 - 5 - 6 - 0 - 1	Small Large No averaging 0.5 second	Refer to "Filter depending on differences in the amount to deliver powdery and liquid material".
	Averaging time for the second moving average	2 3 4 • 5 6	1.0 second1.5 seconds2.0 seconds2.5 seconds3.2 seconds	

Factory setting

## 15-1 Description of "Averaging range" and "Averaging time"

### Averaging range (FI-b) and averaging time (FI-b)

- 1. When the fluctuation of a weight value is beyond the range that is selected in " *F I*-*b* ", the averaging operation is disabled and the display reflects the varying value.
- 2. Once the fluctuation becomes within the selected range, the averaging operation starts to stabilize the weight value.
- 3. The process of averaging increases. When the selected time is reached, moving averaging will be performed.

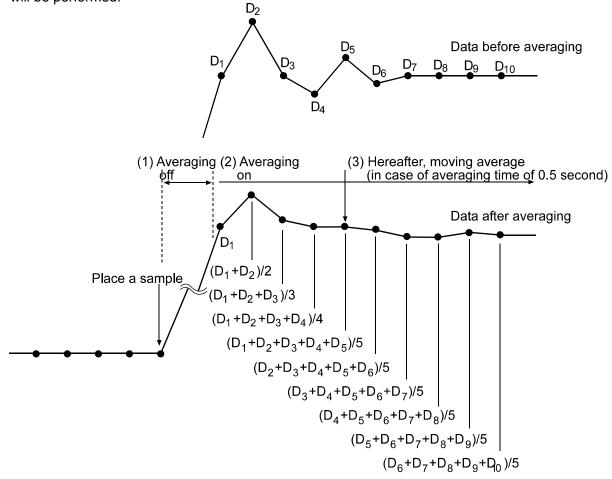

When a small amount of sample is weighed or is filled, the fluctuation of a weight value is too small to be beyond the selected range and the averaging operation is not disabled. Consequently moving averaging is always performed and it takes a longer time to reach the final weight value. Under such a situation, change the setting of "F I-b " to a smaller range. But please note that the smaller the range is, the more prone to external disturbance the value will become.

# Filter depending on differences in the amount to deliver powdery and liquid material

- When the weight value is not stable even without load fluctuations
  - $\Box$  Increase the averaging range for the first moving average parameter (*F I*-*b*).
  - $\Box$  Increase the averaging time for the first moving average parameter (*F I*-*L*).
  - Strengthen the digital filter. (Increase the function setting "Condition ( [and )" parameter.)
- When response is slow during full and medium flow
  - $\Box$  Decrease the averaging range for the first moving average parameter (*F I*-*b* ).
  - □ Weaken the digital filter. (Decrease the function setting "Condition (Land)" parameter.)
- When response is slow during dribble flow
  - $\Box$  Decrease the averaging time for the first moving average parameter (F I-L).
  - □ Weaken the digital filter. (Decrease the function setting "Condition ( Lond )" parameter.
- To increase stability without load fluctuations
  - $\Box$  Increase the averaging range for the second moving average parameter (*F2-b*).
  - $\Box$  Increase the averaging time for the second moving average parameter (*F2-L*).

## **16. MAINTENANCE**

- Do not use organic solvents to clean the balance. Clean the balance with a lint free cloth that is moistened with warm water and a mild detergent.
- Do not disassemble the balance. Contact the local A&D dealer if the balance needs service or repair.
- Use the original packing material for transportation.

## **17. TROUBLESHOOTING**

## 17-1 Checking the Balance Performance and Environment

The balance is a precision instrument. When the operating environment or the operating method is inadequate, correct weighing can not be performed. Place a sample on the pan and remove it, and repeat this several times. If the balance seems to have a problem with repeatability or to perform improperly, check as described below. If improper performance persists after checking, contact the local A&D dealer for repair.

### Checking that the balance performs properly

• Check the balance performance using the self-check function as described in "7-1 Automatic Response Adjustment / Self Check Function".

An error display appears when a malfunction is found.

- Check the balance repeatability using an external weight. Be sure to place the weight in the center of the weighing pan.
- Check the balance repeatability, linearity and calibrated value using external weights with a known value.

### Checking that the operating environment or weighing method is proper Operating environment

- Is the weighing table solid enough (especially for the balance with a minimum weighing value of 0.0001 g)?
- Is the balance level? Refer to "3-1 Before Use".
- Is the operating environment free from vibration and drafts? Has the breeze break been installed?
- Is there a strong electrical or magnetic noise source such as a motor near the balance?

### Weighing method

- Does the weighing pan touch the breeze ring or anything? Is the weighing pan installed correctly?
- Is the RE-ZERO key pressed before placing a sample on the weighing pan?
- Is the sample placed in the center of the weighing pan?
- Has the balance been calibrated using the internal mass (one-touch calibration)?
- Has the balance been warmed up for 30 minutes before weighing?

### Sample and container

- Has the sample absorbed or lost moisture due to the ambient conditions such as temperature and humidity?
- Has the temperature of the container been allowed to equalize to the ambient temperature? Refer to "3-2 During Use".

• Is the sample charged with static electricity? Refer to "3-2 During Use".

Static electricity is generated when the relative humidity is low.

• Is the sample of magnetic material such as iron? Caution is required for weighing magnetic materials. Refer to "3-2 During Use".

## 17-2 Error Codes

Display	Error code	Description
	EC, E11	Stability error
Errori		The balance can not stabilize due to an environmental problem. Prevent vibration, drafts, temperature changes, static electricity and magnetic fields.
		Refer to "3. PRECAUTIONS" for details on the operating environment and "7. RESPONSE ADJUSTMENT" about adapting the balance to the environment.
		To return to the weighing mode, press the CAL key.
	EC, E16	Internal mass error
Errorb		Applying the internal mass does not yield a change in the weighing value as specified.
		Confirm that there is nothing on the pan and perform the weighing operation from the beginning again.
	EC, E17	Internal mass error
Errari		The internal mass application mechanism does not function properly.
		Perform the weighing operation from the beginning again.
	EC, E20	Calibration weight error
		The calibration weight is too heavy. Confirm the calibration weight value.
		Press the CAL key to return to the weighing mode.
	EC, E21	Calibration weight error
		The calibration weight is too light. Confirm the calibration weight value.
		Press the CAL key to return to the weighing mode.

Display	Error code	Description
		Overload error
£		A sample beyond the balance weighing capacity has been placed on the pan.
		Remove the sample from the pan.
۲		Weighing pan Error
- [		The weight value is too light.
		Confirm that the weighing pan is properly installed and calibrate the balance.
		Sample mass error
Lū		The balance can not store the sample for the counting mode or for the percent mode because it is too light.
		Use a larger sample.
		Unit mass error
<u>25-</u> рс 50- рс		The sample unit mass for the counting mode is too light. Storing and using it for counting will cause a counting error.
		Add samples to reach the specified number and press the PRINT key.
	·	Pressing the PRINT key without adding samples will shift the balance to the counting mode. But, for accurate counting, be sure to add samples.
		Automatic response adjustment zero error
		The automatic response adjustment can not be performed because there is something on the pan.
		Clear the pan. Press the CAL key to return to the weighing mode.
		Automatic response adjustment unstable error
CHECK NG)		The automatic response adjustment can not be performed because the weight value is unstable.
		Check the ambient conditions such as breeze, vibration and magnetic fields, also check the weighing pan. Press the CAL key to return to the weighing mode.
		Internal error
<u> </u>		Indicates an internal error as the result of the self-check function.
(020		Repair is required. Contact the local A&D dealer.
		Clock battery error
rte Pf	•	The clock backup battery has been depleted.
		Press any key and set the time and date. The clock and calendar function works normally as long as the AC adapter is connected to the balance. If this error appears frequently, contact the local A&D dealer.

Display	Error code	Description
		Memory full
Ê	(Blinking)	The amount of weighing data in memory has reached the maximum capacity.
	(	Delete the data in memory to store new data. For details, refer to the GX series instruction manual, "12. DATA MEMORY".
<u> </u>		Memory full
FUL	(Illuminated)	The amount of calibration or calibration test data in memory has reached the maximum capacity (50 sets).
		The data in memory will be deleted automatically to store new data. For details, refer to the GX series instruction manual, "12. DATA MEMORY".
		Memory type error
		The type of memory set in the function table and the type of data stored are different.
		For details, refer to the GX series instruction manual, "12. DATA MEMORY".
	EC, E00	Communications error
		A protocol error occurred in communications.
		Confirm the format, baud rate and parity.
	EC, E01	Undefined command error
		An undefined command was received.
		Confirm the command.
	EC, E02	Not ready
		A received command can not be processed.
		e.g. The balance received a Q command, but not in the weighing mode.
		e.g. The balance received a Q command while processing a RE-ZERO command.
		Adjust the delay time to transmit a command.
	EC, E03	Timeout error
		If the timeout parameter is set to " $E$ - $UPI$ ", the balance did not receive the next character of a command within the time limit of one second.
		Confirm the communication.
	EC, E04	Excess characters error
		The balance received excessive characters in a command.
		Confirm the command.

Display	Error code	Description	
	EC, E06	Format error	
		A command includes incorrect data.	
		e.g. The data is numerically incorrect.	
		Confirm the command.	
	EC, E07	Parameter setting error	
		The received data exceeds the range that the balance can accept.	
		Confirm the parameter range of the command.	
Other errors		If the errors described above can not be released or other errors are displayed, contact the local A&D dealer.	
		Other symbol	
		When this indicator blinks, automatic self calibration is required. The indicator blinks when the balance detects changes in ambient temperature. If the balance is not used for several minutes with this indicator blinking, the balance performs automatic self calibration. The blinking duration depends on the operating environment.	

## 17-3 Asking For Repair

If the balance needs service or repair, contact the local A&D dealer.

The balance is a precision instrument. Use much care when handling the balance and observe the following when transporting the balance.

- Use the original packing material.
- Remove the weighing pan from the main unit.

## **18. SPECIFICATIONS**

	MC-1000	MC-6100
Weighing capacity	1100 g	6100 g
Maximum display	1100.0844 g	6100.844 g
Minimum weighing value	0.0001 g	0.001 g
Repeatability (Standard deviation)	0.0005 g / 1 kg to 500 g	0.004 g / 5 kg to 2 kg
See Note 1 below	0.0004 g / less than 500 g	0.0015 g / less than 2 kg
Linearity See Note 1 below	±0.003 g	±0.03 g
Sensitivity drift (10°C-30°C/50°F-86°F with)	±2 ppm/°C (When automatic self calibration is not used)	
Accuracy right after calibration using	±0.010 g	±0.15 g
the internal mass See Note 2 below	The value above is to	the weighing capacity.
Operating environment	5°C to 40°C (4	41°F to 104°F)
	85%RH or less (I	No condensation)
Internal mass	Yes	
Data memory function, weight data	Maxim	um 200
	(Maximum 100 when the	time and date are added)
Calendar function	Yes	
Display refresh rate	5 times/second or 10 times/second	
Display mode	g (g	ram)
Interface (Provided as standard)	RS-232C with Windows Co	ommunication Tools WinCT
External calibration weight	1000 g, 900 g	6000 g
	800 g, 700 g	5000 g
	600 g, 500 g	4000 g
	400 g, 300 g	3000 g
	200 g	2000 g
Weighing pan	128 x 128 mm	165 x 165 mm
Net weight	Approx. 4.6 kg	Approx. 5.1 kg
External dimensions	210 (W) x 317 (D) x 86 (H) mm	
AC adapter	Confirm that the adapter type is correct for the local voltage and power receptacle type	
Power consumption	Approx. 11VA (supplied to the AC adapter)	
Accuracy class F1	500 g, 1 kg	2 kg, 5 kg

Note 1: When the auto-centering pan is used or when loading and unloading are performed at the same place using the automatic loading machine under good ambient conditions.

Note 2: Accuracy right after calibration using the internal mass under good ambient conditions (within the temperature range of 10°C to 30°C (50°F to 86°F) with no abrupt changes in temperature or humidity, no drafts, no effect by magnetic fields or static electricity). The value of the internal mass may change due to corrosion or other damage caused by the operating environment, or due to aging. Check the internal mass using an external weight periodically.

## **19. OPTIONS**

### AX-MC1000PAN Auto-centering pan for the MC-1000

• Consists of an auto-centering pan, a breeze break, two breeze break securing screws and an auto-centering pan guide(128 x 128 mm).

### AX-MC6100PAN Auto-centering pan for the MC-6100

• Consists of an auto-centering pan, a breeze break, two breeze break securing screws and an auto-centering pan guide (165 x 165 mm).

### AX-HDB-25P/CTF Connector

### AX-KO1710-200 RS-232C Cable

• Length 2 m, straight type, D-sub 9pin - D-sub 25pin.

### AX-SW128 Foot switch

• Used to externally transmit a RE-ZERO or PRINT signal to the balance.

### AX-USB-25P-EX USB converter

• An RS-232C cable is provided to connect the USB converter to the balance.

### AD-8121B Printer

- Compact dot-matrix printer
- Statistical function, clock and calendar function, interval print function, graphic print function, dump print mode
- 5 x 7 dots, 16 characters per line
- Print paper (AX-PP143, 45 (W) x 50 (L) mm , ø65 mm)
- AC adapter or alkaline battery.

### AD-8524A/B Keyboard adapter

• Used to connect the balance to a personal computer with appropriate OS and applications.

#### AD-8920A Remote display

• Connected to the MC series using the RS-232C interface to display the weighing data away from the balance.

### AD-8922A Remote controller

• Connected to the MC series balance using the RS-232C interface to display the weighing data and to remotely control the balance.

#### AD-1682 Rechargeable battery

• Allows use of the balance in a place where AC power is not available.

#### AD-1683 DC static eliminator

• Used to minimize weighing errors due to static electricity on the material.



#### AD-1684 Electrostatic field meter

• Measures the amount of the static charge on the sample, tare or peripheral equipment and displays the result.

If those are found to be charged, discharge them using the AD-1683 DC static eliminator.

#### AD-1687 Weighing environment logger


- A data logger equipped with 4 sensors for temperature, humidity, barometric pressure and vibration that can measure and store environmental data.
- When connected to the RS-232C interface of the balance, the AD-1687 can store environmental data along with weighing data. Therefore, it is possible to store data in an environment where a computer can not be used.
- The stored data can be read to a personal computer using USB. As the AD-1687 is recognized as USB memory, special software is not required to read the data.

#### AD-1688 Data logger

- When connected to the RS-232C interface of the balance, the AD-1688 can store the data in an environment where a personal computer can not be used.
- The stored data can be read to a personal computer using USB. As the AD-1688 is recognized as USB memory, special software is not required to read the data.

#### AD-1689 Tweezers for calibration weight

• Used when calibrating the balance using an external weight.

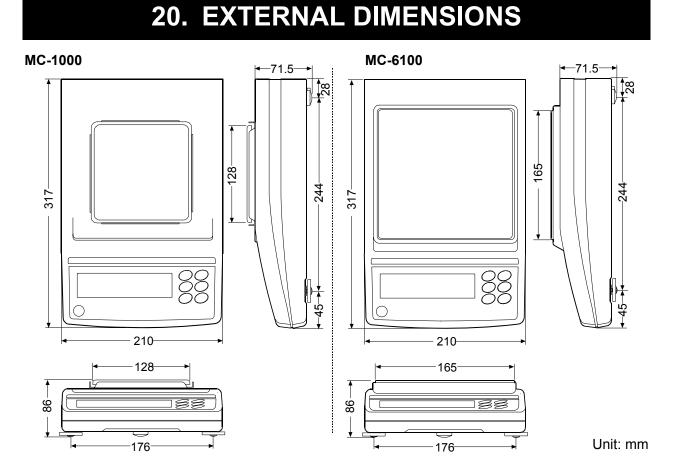

### 63

## 21. TERMS/INDEX

### Terms

Terms	
Stable value	The weight data when the stabilization indicator appears.
Environment	Ambient conditions such as vibration, drafts, temperature, static electricity and
	magnetic fields which affect the weighing operation.
Store	To save the weighing data, unit mass or calibration data using the data memory
	function.
Calibration	Adjustment of the balance so that it can weigh accurately.
Output	To output the weighing data using the RS-232C interface.
Zero point	A weighing reference point or the zero display. Usually refers to the value
	displayed when nothing is on the weighing pan.
Data number	Numbers assigned sequentially when weighing data or unit weight is stored.
Digit	Unit of digital resolution. Used for the balance, a unit of minimum weighing value.
Tare	To cancel the weight of a container which is not included in the weighing data.
Mode	Balance operational function.
Re-zero	To set the display to zero.
GLP	Good Laboratory Practice.
Repeatability	Variation in measured values obtained when the same weight is placed and
	removed repetitively. Usually expressed as a standard deviation.
	e.g. Standard deviation=1 digit: This means that measured values fall within $\pm 1$
	digit in the frequency of about 68%.
Stabilization time	Time required after a sample being placed, until the stabilization indicator
	illuminates and the weighing data is displayed.
Sensitivity drift	An affect that a change in temperature causes to the weighing data. Expressed
	as temperature coefficient.
	e.g. Temperature coefficient = 2 ppm/°C : If a load is 500 g and the temperature
	changes by 10°C, the value displayed changes by the following value.
	0.0002%/°C x 10°C x 500 g = 10 mg

## Index

## Keys and symbols

(CAL)	CAL key	13
	MODE key	13, 16
UC ON:OFF	ON/OFF key	13, 28
	PRINT key	13
+0/1+ RE-ZERO	RE-ZERO key	11, 13
1/10d SAMPLE	SAMPLE key	13
ANIMAL	Animal mode indicator	13
HI OK L	Comparator indicators	13
	Interval memory active indic	ator13
	Interval memory standby inc	dicator.13
*	Prior notice indication of aut	omatic
	self calibration	13
-	Processing indicator	13
FAST M	D. SLOW Response indicators	13
°	Stabilization indicator	13
	Standby indicator	13

### - A -

	37
4	, 7
4	, 7
4	, 7
	62
11,	62
	63
	62
,	
-	
53,	54
16,	62
	4 4 4 4  t 47,  32, 32,

AX-MC6100PAN	9, 16, 62
AX-SW128 Foot switch	46, 62
AX-USB-25P-EX USB converter	62

### - B -

ьрс	47
Breeze ring	
Bubble spirit level	4, 7
Building into a system	17

### - C -

Calibration Calibration test	
Capacity indicator	13, 31
EL AdJ	31
Cond	16
[Р	31
[P Fnc	31
СР НН	32
[P in	31
[P LL	32
[Р Lo	32
CSV format	39

### - D -

dRER	32
Data number	39, 64
Decimal point	31
Digit	31, 32, 64
Display refresh rate	31, 61
d-no	32
dout	32, 33, 47
DP (Dump print) format	37
Dust guard	4, 7

### - E -

EC, E00	59
EC, E01	59
EC, E02	59
EC, E03	59
EC, E04	59
EC, E06	60
EC, E07	60
EC, E11	57
EC, E16	57
EC, E17	57
EC, E20	57
EC, E21	57
Eccentric loading error	3, 6, 9

Environment	64
Extended function	34, 53
External dimensions	61, 63
External input	45

## - F -

F I-6	 3
FIL	 1
F I-E	 3
F2-B	 3
F2-E	 3
FAST	 3

## - G -

Glass breeze break	
GLP	
Grounding terminal4, 7	

## - H -

Hold function31
-----------------

## - / -

ıd	
ID number	41
Inhibit	
Initializing the Balance	
Insulators	11
Int	32
Interval memory mode	

## - K -

Key mode	
KF format	

## - L -

Leveling foot	4, 7
Linearity	. 56, 61

## - M -

Main unit cover	4, 7
Mass comparator	3, 16, 31
Memory full	59
MID	18
MT format	

### - N -

NU (Numerical) format38
-------------------------

- 0 -	
Output	64

## - P -

Pan sup	port receptor	5, 6, 8, 9
Permit		28
PoFF		31, 33
Prt		32
P-tr		31

## - R -

Reference card	4, 7
Repeatability	56, 61, 64
RESPONSE	
Response characteristic	
RS-232C serial interface	45
RsCom	
RsKey	48

## - S -

Self check function	18
Sensitivity drift	61, 64
SLOW	18
5Pn	31
5-Ed	32
Span calibration	17, 31, 35
Stability band width	
Stabilization time	64
Stable value	18, 35, 64
Static electricity	11, 57
Stream mode	36

## - U -

Underho	ook	11
טה וב		33
Units		14

## - W -

Warm up	10, 11, 12
Weighing capacity	11, 58, 61
Weighing pan	, 4, 7, 61
WinCT	4, 7, 48

### - Z -

Zero point	23, 25, 64
Zero tracking	31
Zero upon power-on	17, 31, 35

MEMO

MEMO

### 



### A&D Company, Limited

3-23-14 Higashi-Ikebukuro, Toshima-ku, Tokyo 170-0013 JAPAN Telephone: [81] (3) 5391-6132 Fax: [81] (3) 5391-6148

### A&D ENGINEERING, INC.

1756 Automation Parkway, San Jose, California 95131 U.S.A. Telephone: [1] (408) 263-5333 Fax: [1] (408) 263-0119

### A&D INSTRUMENTS LIMITED <UK Office>

Unit 24/26 Blacklands Way, Abingdon Business Park, Abingdon, Oxfordshire OX14 1DY United Kingdom Telephone: [44] (1235) 550420 Fax: [44] (1235) 550485

### A&D INSTRUMENTS LIMITED <German Sales Office>

Große Straße 13 b 22926 Ahrensburg Deutschland Telefon: [49] (0) 4102 459230 Telefax: [49] (0) 4102 459231

### A&D Australasia Pty Ltd.

32 Dew Street, Thebarton, South Australia 5031 AUSTRALIA Telephone: [61] (8) 8301-8100 Fax: [61] (8) 8352-7409

### A&D KOREA Limited 한국에이.엔.디(주)

대한민국 서울시 영등포구 여의도동 36-2 맨하탄 빌딩 8층 (8th Floor, Manhattan Bldg. 36-2 Yoido-dong, Youngdeungpo-ku, Seoul, KOREA) 전화: [82] (2) 780-4101 팩스: [82] (2) 782-4280

A&D RUS CO., LTD.Компания ЭЙ энд ДИ РУС121357, Российская Федерация, г.Москва, ул. Верейская, дом 17( Bldg. 17, Vereyskaya st., Moscow, 121357тел.: [7] (495) 937-33-44факс: [7] (495) 937-55-66

 A&D Instruments India Private Limited
 ऐ&डी इन्स्ट्रयूमेन्ट्स इण्डिया प्रा0 लिमिटेड

 509, उद्योग विहार , फेस -5, गुड़गांव - 122016, हरियाणा , भारत
 (509, Udyog Vihar, Phase-V, Gurgaon - 122 016, Haryana, India )

 फोन : 91-124-4715555
 फैक्स : 91-124-4715599